



DOC022.53.00654

**DR 5000**

USER MANUAL

January 2008, Edition 2



<b>Section 1 Specifications</b> .....	7
<b>Section 2 General Information</b> .....	9
2.1 Safety information .....	9
2.1.1 Use of hazard information .....	9
2.1.2 Precautionary labels .....	9
2.1.3 Class 1 LASER .....	10
2.1.4 Chemical and Biological Safety .....	10
2.1.5 Source Lamp Safety .....	10
2.2 Overview of product .....	11
<b>Section 3 Installation</b> .....	13
3.1 Unpack the instrument .....	13
3.2 Operating environment .....	13
3.3 Connection .....	14
3.4 Cell compartment, Multi Cell Holder, 100 mm Cell Adapter (optional) .....	15
3.4.1 100 mm adapter (optional) .....	16
3.4.2 Installation Multi Cell Holder .....	17
3.5 Beam path .....	18
<b>Section 4 Start Up</b> .....	19
4.1 Power the instrument on and off .....	19
4.2 Language selection .....	19
4.3 Self-Check .....	19
<b>Section 5 Standard Operations</b> .....	21
5.1 Overview .....	21
5.1.1 Tips for the use of the touch screen .....	21
5.1.2 Use of the alphanumeric keypad .....	21
5.1.3 Main Menu .....	22
5.2 Instrument Setup mode .....	23
5.2.1 Operator ID .....	23
5.2.2 Sample ID .....	24
5.2.3 Date and time .....	25
5.2.4 Display and sound preferences .....	25
5.2.5 Lamp Control .....	26
5.2.5.1 VIS-Lamp and UV-Lamp .....	26
5.2.5.2 Effect caused by the lamp settings on the Measuring Mode .....	27
5.2.6 PC and printer .....	29
5.2.6.1 Printer setup .....	29
5.2.6.2 Print data .....	31
5.2.6.3 HACH Data Trans .....	31
5.2.7 Password .....	32
5.2.7.1 Password deactivation .....	33
5.3 Store, recall, send and delete data .....	35
5.3.1 The Data Log .....	35
5.3.1.1 Auto/manual data storage .....	35
5.3.1.2 Recall stored data from the data log .....	35
5.3.1.3 Send data from the data log .....	36
5.3.1.4 Delete stored data from the data log .....	37
5.3.2 Store, recall, send and delete data from wavelength scan and time course .....	38
5.3.2.1 Data storage from wavelength scan or time course .....	38
5.3.2.2 Recall stored data from wavelength scan or time course .....	38
5.3.2.3 Send data from wavelength scan or time course .....	39
5.3.2.4 Delete stored data from wavelength scan or time course .....	41

## Table of Contents

---

5.4 Stored Programs .....	42
5.4.1 Select a saved test/method; enter user-specific basic data .....	42
5.4.2 Stored program options .....	42
5.4.3 Use of program timers .....	44
5.4.4 Set the reading mode .....	44
5.4.4.1 Take single wavelength measurements (single reading) .....	45
5.4.4.2 Take single wavelength measurements (continuous readings) .....	45
5.4.5 Set the dilution factor .....	45
5.4.6 Run a standard adjust .....	46
5.4.7 Set the chemical form .....	47
5.4.7.1 Change of the default setting of the chemical form .....	47
5.4.8 Run a reagent blank .....	48
5.4.9 Analysis of samples .....	49
5.4.10 Add stored programs to the favorite programs list .....	49
5.5 Barcode Programs .....	50
5.5.1 Complete a barcode 13 mm test/vial .....	51
5.5.2 Select the measuring range .....	51
5.5.3 Select the chemical evaluation form .....	51
5.5.3.1 Change of the default setting of the chemical form .....	52
5.5.4 Basic test-specific and sample-specific data settings .....	52
5.5.5 Sample blank .....	53
5.5.6 Update/edit barcode tests .....	54
5.5.6.1 Manual update of a barcode test .....	54
5.5.6.2 Update an existing barcode test .....	57
5.5.6.3 Program a new test .....	58
5.5.7 Upgrade of the instrument software .....	58
<b>Section 6 Advanced Operations .....</b>	<b>61</b>
6.1 User Programs .....	61
6.1.1 Program a user method .....	61
6.1.1.1 Single wavelength settings .....	63
6.1.1.2 Multi wavelength settings .....	64
6.1.1.3 Calibration settings for single and multi wavelength mode .....	66
6.1.1.4 Store a user program .....	70
6.1.1.5 Additional user-defined parameters and functions .....	70
6.1.2 Free programming program type .....	72
6.1.2.1 Measurement process .....	72
6.1.2.2 Enter a new element of a measuring sequence .....	73
6.1.2.3 Enter the calibration formula (evaluation formula) .....	75
6.1.2.4 Enter variables .....	78
6.1.2.5 Save a free programming user program .....	78
6.1.3 Select a user program .....	79
6.1.4 Add, edit and delete user programs from the favorites list .....	79
6.1.4.1 Add to Favorites .....	80
6.1.4.2 Edit .....	80
6.1.4.3 Delete .....	80
6.2 Favorite Programs .....	81
6.2.1 Recall a favorite program .....	81
6.2.2 Delete a favorite program .....	81
6.3 Standard Addition – monitoring/checking results .....	82
6.3.1 Complete a standard addition .....	83

6.4 Single Wavelength (absorbance, concentration and transmittance measurements)	87
6.4.1 Set up single wavelength mode	87
6.4.2 Take single wavelength measurements (single reading)	89
6.4.3 Take single wavelength measurements (continuous readings)	89
6.5 Multi-Wavelength mode – measurements at more than one wavelength	90
6.5.1 Set the reading mode at different wavelengths	90
6.5.2 Complete a measurement in the multi wavelength mode	93
6.6 Wavelength Scan mode – recording of absorbance and transmittance spectrums	94
6.6.1 Set up the wavelength scan	94
6.6.2 Wavelength scan reading	97
6.6.2.1 Navigation of the wavelength scan graph or a wavelength scan analysis	98
6.6.3 Work with reference scans	99
6.7 Time course of absorbance/transmittance	101
6.7.1 Time course setup parameters	101
6.7.2 Time course scan reading	102
6.7.3 Analysis of time course data	103
6.7.3.1 Navigation of a time scan or a time scan analysis	104
6.8 System checks	105
6.8.1 Instrument information	105
6.8.2 Upgrade of the instrument software	106
6.8.3 Optical checks	106
6.8.3.1 Wavelength check	107
6.8.3.2 Noise check	108
6.8.3.3 Stray light check	110
6.8.3.4 Absorbance check	111
6.8.3.5 Drift check	112
6.8.3.6 Verification kit	113
6.8.4 Output checks	115
6.8.5 Lamp history	115
6.8.6 Factory service	116
6.8.7 Service time	116
6.8.8 Instrument Backup	117
<b>Section 7 Maintenance</b>	121
7.1 Cleaning requirements	121
7.1.1 Spectrophotometer	121
7.1.2 Display	121
7.1.3 Cuvettes/sample cells	121
7.2 Lamp replacement	122
7.2.1 Tungsten lamp replacement	122
7.2.2 Deuterium lamp (UV) replacement	123
7.3 Fuse replacement	124
7.4 Filter pad maintenance	125
7.4.1 Filter pad replacement	125
<b>Section 8 Troubleshooting</b>	127
<b>Section 9 Replacement Parts</b>	129
<b>Section 10 Contact Information</b>	131
<b>Section 11 Limited Warranty</b>	133
<b>Index</b>	135



# Section 1 Specifications

Specifications are subject to change without notice.

Performance specifications	
Operating mode	Transmittance (%), Absorbance and Concentration
Source lamp	Gas-filled Tungsten (visible) and Deuterium (UV)
Wavelength range	190–1100 nm
Wavelength accuracy	± 1 nm in Wavelength Range 200–900 nm
Wavelength reproducibility	< 0.1 nm
Wavelength resolution	0.1 nm
Wavelength calibration	Automatic
Wavelength selection	Automatic, based on method selection
Scanning speed	900 nm/min 1 nm steps
Spectral bandwidth	2 nm
Photometric range	± 3.0 Abs in Wavelength Range 200–900 nm
Photometric accuracy	5 mAbs at 0.0–0.5 Abs 1% at 0.50–2.0 Abs
Photometric linearity	< 0.5%–2 Abs < = 1% at > 2 Abs with neutral glass at 546 nm
Stray light	KI-solution at 220 nm > 3.3 Abs / < 0.05%
Data storage	1000 measured values (result, date, time, sample ID, user ID)
User programs	50
Physical and environmental specifications	
Width	450 mm (17.7 in)
Height	200 mm (7.9 in)
Depth	500 mm (19.7 in)
Weight	15.5 kg (34.2 lbs)
Operating requirements	10–40 °C (50–104 °F), max. 80% relative humidity (non-condensing)
Storage requirements	–25–60 °C (–13–140 °F) max. 80% relative humidity (non-condensing)
Additional technical data	
Mains connection	100–240 V~; 50/60 Hz; automatic changeover
Interfaces	Use only shielded cable with max. length of 3 m. 1 x USB type B 2 x USB type A
Water-resistant to Enclosure	IP 31
Safety class	Safety class I





# Section 2 General Information

## 2.1 Safety information

Please read this entire manual before unpacking, setting up or operating this equipment. Pay attention to all danger, warning and caution statements. Failure to do so could result in serious injury to the operator or damage to the equipment.

Make sure that the protection provided by this equipment is not impaired, do not use or install this equipment in any manner other than that specified in this manual.

### 2.1.1 Use of hazard information

#### **DANGER**

*Indicates a potentially or imminently hazardous situation which, if not avoided, will result in death or serious injury.*

#### **WARNING**

*Indicates a potentially or imminently hazardous situation which, if not avoided, could result in death or serious injury.*

#### **CAUTION**




*Indicates a potentially hazardous situation that may result in minor or moderate injury.*

**Important Note:** *Indicates a situation which, if not avoided, may cause damage to the instrument. Information that requires special emphasis.*

**Note:** *Information that supplements points in the main text.*

### 2.1.2 Precautionary labels

Read all labels and tags attached to the instrument. Personal injury or damage to the instrument could occur if not observed. A symbol, if noted on the instrument, will be included with a danger or caution statement in the manual.

	This symbol, if noted on the instrument, references the instruction manual for operation and/or safety information.
	This symbol, if noted on the instrument, indicates a hot surface.
	Electrical equipment marked with this symbol may not be disposed of in European public disposal systems after 12 August of 2005. In conformity with European local and national regulations (EU Directive 2002/96/EC), European electrical equipment users must now return old or end-of life equipment to the Producer for disposal at no charge to the user. <b>Note:</b> <i>For return for recycling, please contact the equipment producer or supplier for instructions on how to return end-of-life equipment, producer-supplied electrical accessories and all auxiliary items for proper disposal.</i>
<b>LASER CLASS 1</b>	This symbol indicates that the instrument contains a Class 1 LASER device. Data: 0.3 mW; $\lambda = 650 \text{ nm}$

### 2.1.3 Class 1 LASER

A Class 1 LASER is installed in this instrument. Class 1 LASERS are products where the radiant power of the LASER beam accessible (the accessible emission) is always below the Maximum Permissible Exposure value. Therefore, for Class 1 LASERS the output power is below the level at which it is believed eye damage will occur. Exposure to the beam of a Class 1 LASER will not result in eye injury. Class 1 LASERS may therefore be considered safe. However, Class 1 LASER products may contain LASER systems of a higher Class but there are adequate engineering control measures to ensure that access to the beam is not reasonably likely. Examples of such products include LASER printers and compact disc players. CDRH assessment number 0510555-02.

Data: 0.3 mW; wavelength = 650 nm

### 2.1.4 Chemical and Biological Safety

#### **DANGER**

***Potential Chemical/ Biological Exposure Hazards. Handling chemical samples, standards and reagents can be dangerous. Users of this product are advised to familiarize themselves with safety procedures and the correct use of chemicals, and to carefully read all relevant Material Safety Data Sheets.***

Normal operation of this instrument may involve the use of hazardous chemicals or biologically harmful samples.

- The user must observe all cautionary information printed on the original solution containers and safety data sheet prior to their use.
- All waste solutions must be disposed in accordance with local and national law.
- The type of protective equipment must be selected according to the concentration and amount of the dangerous substance at the specific workplace.

### 2.1.5 Source Lamp Safety

The source lamp operate at high temperatures.

- To avoid a possible electric shock, disconnect the instrument from the power source before servicing the lamp. Make sure that the instrument cannot be accidentally switched on again.

#### **CAUTION**

***The UV lamp generates UV light.***

***Do not look directly at an operating lamp without wearing UV protective eye glasses.***

***Important Note: To prevent a possible burn, allow the lamp to cool at least 30 minutes before handling.***

## 2.2 Overview of product

The DR 5000 Spectrophotometer is a complete scanning UV/VIS spectrophotometer with a wavelength range of 190 to 1100 nm. The instrument comes with a complete set of application programs and multi-language support.

The DR 5000 Spectrophotometer contains the following application modes: Stored Programs, User Programs, Favorite Programs, Single Wavelength Mode, Multi-Wavelength Mode, Wavelength Scanning Mode and Time Course Mode.

The DR 5000 is used for testing in visible and ultraviolet wavelengths. A gas-filled tungsten lamp produces light in the visible spectrum (320 to 1100 nm), and a deuterium lamp available produces light in the ultraviolet spectrum (190 to 360 nm).

The DR 5000 Spectrophotometer provides digital readouts in direct concentration units, absorbance, or percent transmittance.

When a user-generated or programmed method is selected, the on-screen menus and prompts direct the user through the test.

This menu system also can be used to generate reports, statistical evaluations of generated calibration curves, and to report instrument diagnostic checks.



### **WARNING**

**Electrical and Fire Hazards. Use only the provided power supply. Only qualified personnel should conduct the tasks described in this section of the manual.**

### 3.1 Unpack the instrument

The DR 5000 Spectrophotometer comes packaged with the following items:

- DR 5000 Spectrophotometer
- Mains power cable
- Multi Cell Holder
- DR 5000 User Manual
- CD-ROM containing the HACH and HACH LANGE procedures

**Note:** *If any of these items are missing or damaged, contact the manufacturer or a sales representative immediately.*

### 3.2 Operating environment

The following conditions are necessary to ensure correct instrument operation and accurate results:

- Place the instrument firmly on an even surface. Do not push any objects under the instrument.
- Maintain an ambient temperature of 10 to 40 °C (50 to 104 °F) for proper instrument operation.
- The relative humidity should be less than 80%; moisture should not condense on the instrument.
- Leave at least a 15 cm (6 in.) clearance at the top and on all sides for air circulation to avoid overheating of electrical parts.
- Do not operate or store the instrument in extremely dusty, damp or wet locations.
- Keep the surface of the instrument, the cell compartment and all accessories clean and dry at all times. Splashes or spills on and in the instrument should be cleaned up immediately (see section 7.1 on page 121).

**Important Note:** *Protect the instrument from temperature extremes, including heaters, direct sunlight and other heat sources.*

### 3.3 Connection

The DR 5000 has three USB interfaces as a standard feature, located on the back of the instrument (Figure 1).

The USB Type A interfaces are used for communications with a printer, USB memory stick or keyboard. A USB memory stick is used to update instrument software.

The USB Type B interface is used for communications with a PC. The optional Hach Data Trans software (see section 5.2.6.3 on page 31) must be installed on the PC for this use.

A USB hub may be used to connect several accessories at a time.

**Note: USB cables must not be longer than 3 meters (10 feet).**

These USB interfaces enable data and graphics to be output to a Printer and a PC and upgrade of instrument software (see section 6.8.2 on page 106).

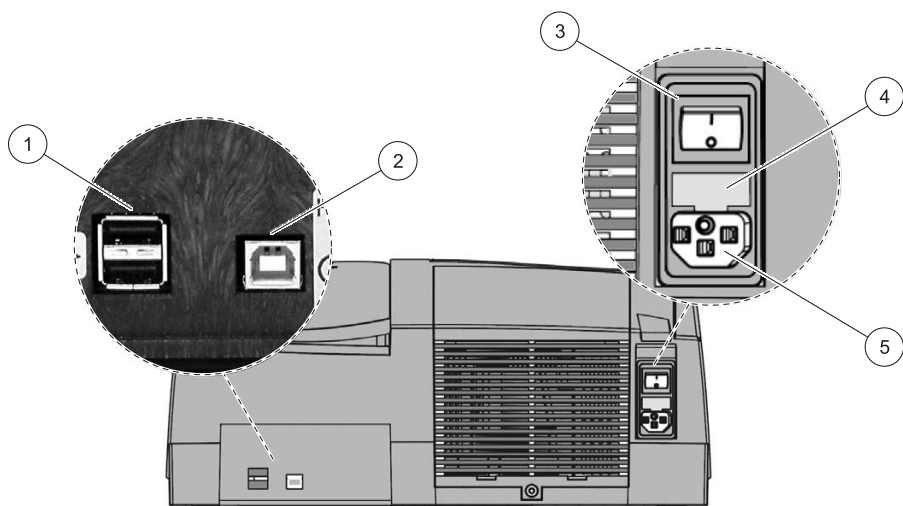


Figure 1 Interfaces

1	USB type A	4	Fuse
2	USB type B	5	Plug in power supply
3	On/Off switch		

### 3.4 Cell compartment, Multi Cell Holder, 100 mm Cell Adapter (optional)

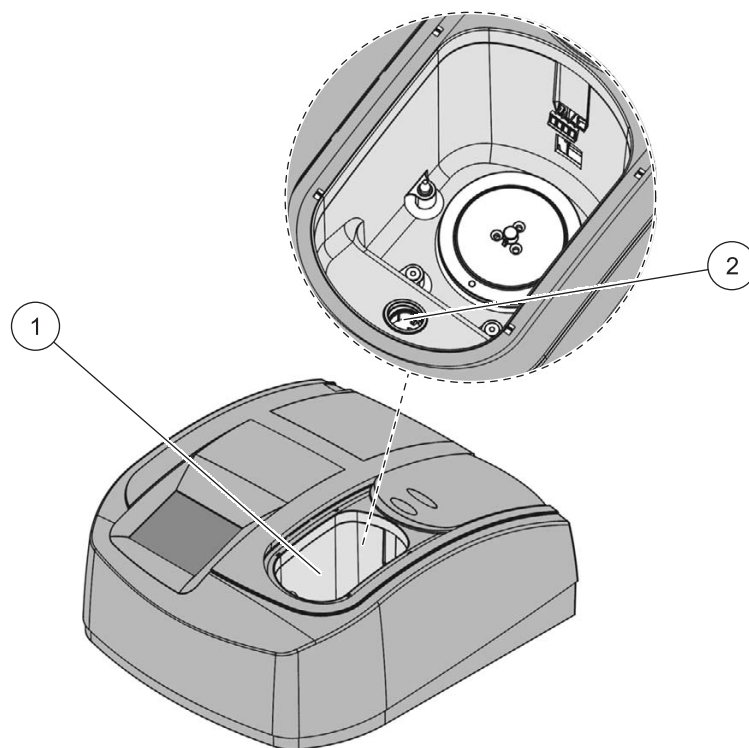


Figure 2 Cell compartment

1 Cell compartment	2 Cell compartment for 13 mm or 16 mm round cuvettes/vials only
--------------------	---

The DR 5000 Spectrophotometer comes equipped with a Multi Cell Holder (Figure 3 on page 16) which is the Standard Holder, provided with each instrument. The Multi Cell Holder can accommodate the following cell types:

- 10 mm square cells
- 20/50 mm rectangular cells
- 1" round cells
- 1" square cells

Only one cell type can be used for a measurement. On the top and bottom of the Multi Cell Holder are a variety of openings to accommodate different types of cells. Beside each opening is printed the type of cell for which it is intended.

Optionally there is a 100 mm adapter available.

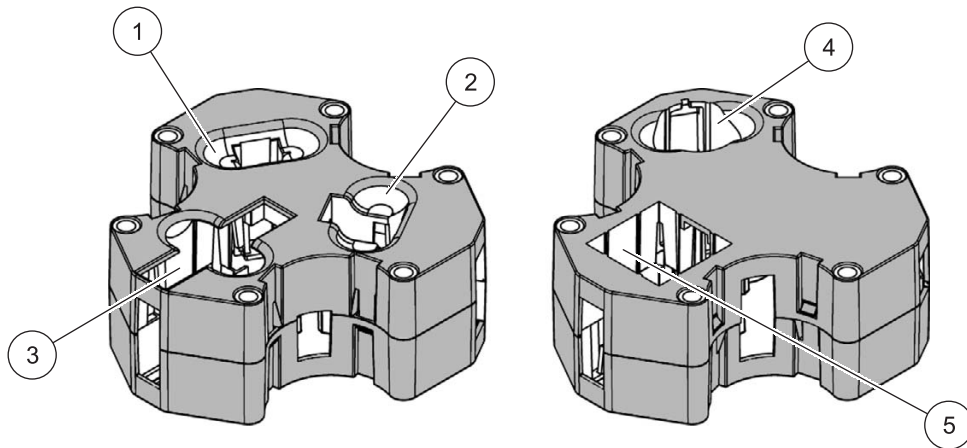


Figure 3 Multi Cell Holder (Top and Bottom)

<b>1</b> 10 mm square cell	<b>4</b> 1" round cell
<b>2</b> 20 mm rectangular cell	<b>5</b> 1" square cell
<b>3</b> 50 mm rectangular cell	

### 3.4.1 100 mm adapter (optional)

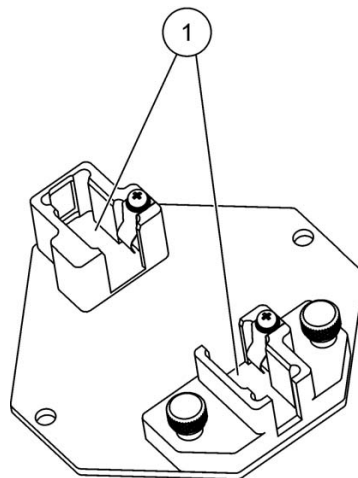


Figure 4 100 mm Cell Adapter

<b>1</b> 100 mm rectangular cell
----------------------------------



### 3.4.2 Installation Multi Cell Holder

Assembly of the Multi Cell Holder:

1. Open the cell compartment.
2. Identify the correct opening for the selected cell type in the Multi Cell Holder.
3. Insert the Multi Cell Holder in the cell compartment in such a way that the name of the selected cell type can be read directly and the cell opening is at the front. Secure the holder with two locking screws (Figure 5).

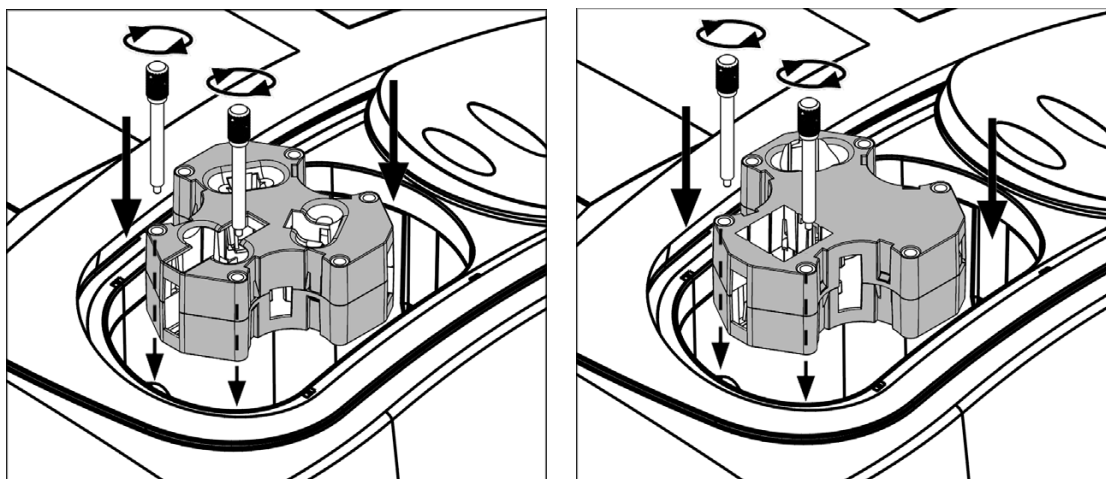


Figure 5 Installation Multi Cell Holder

### 3.5 Beam path

Figure 6 shows the beam path of the DR 5000:

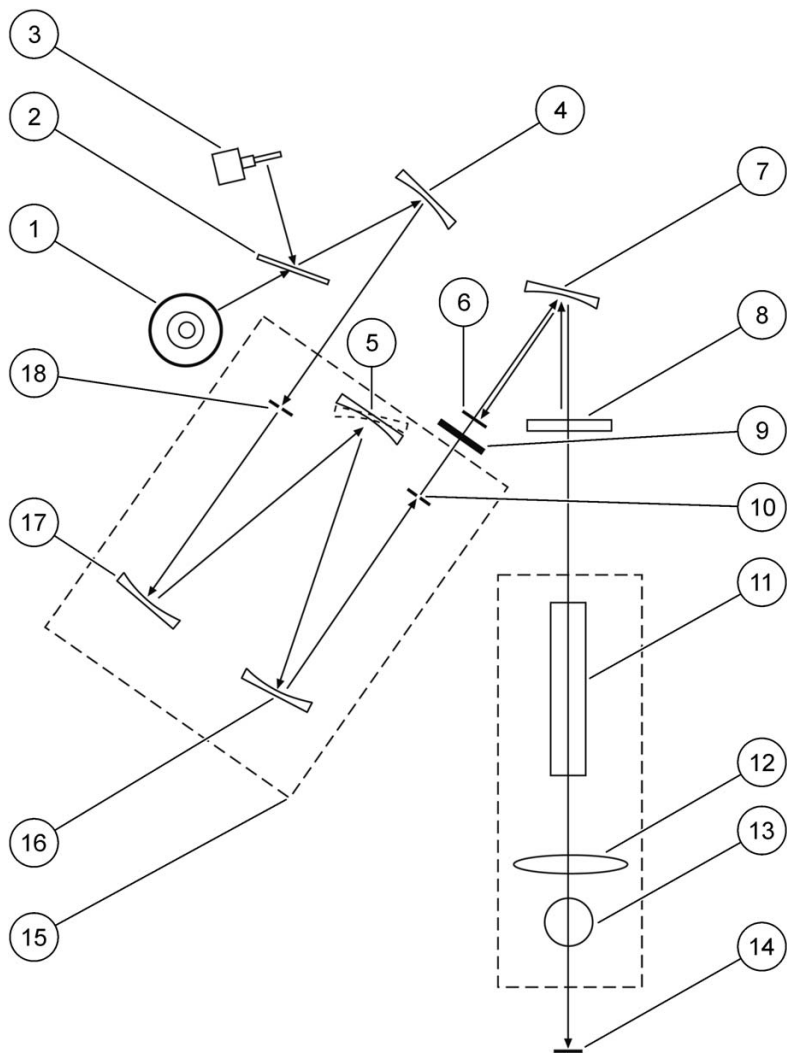


Figure 6 Beam path

1 Deuterium lamp	10 Exit slit
2 Lamp selection mirror	11 Cell compartment
3 Tungsten lamp	12 Lens
4 Mirror	13 Round cuvette/vial compartment
5 Grating	14 Measurement element
6 Reference element	15 Monochromator
7 Mirror	16 Mirror
8 Splitter mirror	17 Mirror
9 Filter	18 Entrance slit

## Section 4 Start Up

### 4.1 Power the instrument on and off

1. Plug in power supply.
2. Close the empty cell compartment.
3. Turn the instrument on by pressing the power switch on the back.

**Note:** Do not turn the instrument off and on in rapid succession. Always wait about **20 seconds** before turning the instrument on again, otherwise the electronic and mechanical systems will be damaged.

### 4.2 Language selection



The spectrophotometer software includes several language options. The first time the instrument is turned on, the language selection screen will appear.


1. Select the desired language.
2. Press **OK** to confirm the language selection. The self check will start automatically.

#### Changing the language setting

The instrument functions in the selected language until the option is changed.

1. While turning the instrument on, touch the screen at any point until the list for selecting a language appears (about 30 seconds).
2. Select the required language.
3. Press **OK** to confirm. The test program subsequently starts automatically.

### 4.3 Self-Check

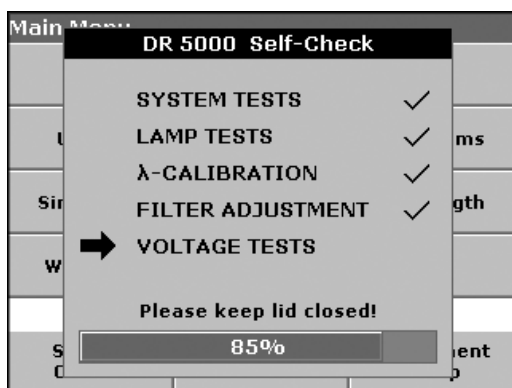

Each time the instrument is powered up, a series of diagnostic tests are performed automatically to ensure operation of major system components.

This procedure, which takes approximately two minutes, checks the system, lamp, filter adjustment, wavelength calibration and voltage. Each test which functions correctly is confirmed with a check mark.

The Main Menu is displayed when power up diagnostics are completed.

**Note:** Further error messages during self check, see Section 8 on page 127.



# Section 5 Standard Operations


## 5.1 Overview

### 5.1.1 Tips for the use of the touch screen

The entire screen is touch-activated. To make a selection, press the screen with a fingernail, fingertip, pencil eraser or a stylus. Do not press the screen with a sharp object, such as the tip of a ball point pen.

- Do not place anything on top of the screen, to prevent damage or scratching on the screen.
- Press keys, words or icons to select them.
- Use scroll bars to move up and down long lists very quickly. Press and hold the scroll bar, then move up or down to move through the list.
- Highlight an item from a list by pressing it once. When the item has been successfully selected, it will be displayed as reversed text (light text on a dark background).

### 5.1.2 Use of the alphanumeric keypad

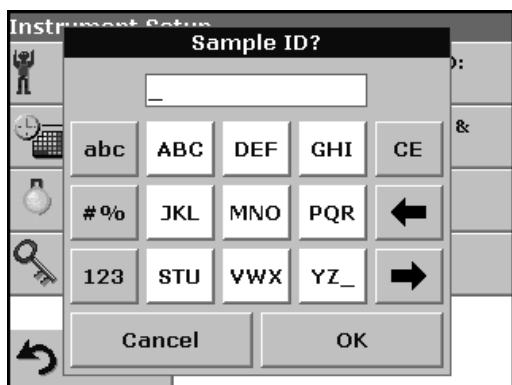

This display is used to enter letters, numbers and symbols as needed when programming the instrument. Unavailable options are disabled (grayed out). The icons on the right and left of the screen are described in Table 1.


The central keypad changes to reflect the chosen entry mode. Press a key repeatedly until the desired character appears on the screen. A space can be entered by using the underscore on the YZ\_ key.

**Note:** A USB keyboard (with US keyboard layout) or a USB Barcode handset scanner can be used for input (see Section 9 on page 129).

Table 1 Alphanumeric keypad

Icon / key	Description	Function
ABC/abc	Alphabetic	When entering alphabetic characters (ex. user-entered units), this key allows to toggle between upper and lower case letters.
# %	Symbols	Punctuation, symbols and numerical sub- and superscripts may be entered.
123	Numeric	For entering regular numbers.
CE	Clear Entry	Clear the entry.
Left Arrow	Backspace	Moves back one position. This deletes the character previously entered in the new position.
Right Arrow	Advance	Moves to the next space in an entry when two adjacent characters occur on the same key.

## 5.1.3 Main Menu


Main Menu			
Stored Programs			
User Programs		Favorite Programs	
Single Wavelength		Multi - Wavelength	
Wavelength Scan		Time Course	
System Checks		Recall Data	Instrument Setup

A variety of modes may be selected from the Main Menu. The following table briefly describes each menu option.

**Table 2 Main Menu options**

Option	Function
<b>Stored Programs (HACH LANGE Programs)</b>	Stored programs are pre-programmed methods that make use of HACH reagents and LANGE cuvette tests and pipette tests. The DR 5000 Procedures Manual contains illustrated, step-by-step procedures for analyses using HACH programs. The working procedures for LANGE tests are included in the test packs.
<b>User Programs</b>	User programs make "made to measure analysis" possible: –Users can program methods they have developed themselves –Existing HACH and LANGE methods can be stored as user programs. The LANGE tests can then be modified to suit the user's requirements.
<b>Favorite Programs</b>	List of methods/tests created by the user to suit his own requirements.
<b>Single Wavelength</b>	Single wavelength measurements are: <b>Absorbance measurements:</b> The light absorbed by the sample is measured in absorbance units. <b>Transmittance measurements (%):</b> The percentage of the light that passes through the sample and reaches the detector is measured. <b>Concentration measurements:</b> A concentration factor can be entered to enable the measured absorbance values to be converted into concentration values.
<b>Multi Wavelength</b>	In the multi-wavelength mode, absorbance (Abs) or percentage transmittance (%T) is measured at up to four wavelengths and absorbance differences and absorbance relationships are calculated. Simple conversions into concentrations can also be carried out.
<b>Wavelength Scan</b>	A wavelength scan shows how the light from a sample is absorbed over a defined wavelength spectrum. This function can be used to determine the wavelength at which the maximum absorbance value can be measured. The absorbance behavior is displayed graphically during the scan.
<b>Time Course</b>	The time scan records the absorbance or % transmittance at a wavelength over a defined time.
<b>System Checks</b>	The system checks menu offers a number of options, including optical checks, output checks, lamp history, instrument update, service time and instrument backup.
<b>Recall Data</b>	Stored data can be recalled, filtered, sent and deleted.
<b>Instrument Setup</b>	In this mode, user-specific or method-specific settings can be entered: Operator-ID, Sample-ID, Date & Time, Display & Sound, Lamp Control, PC & Printer, Password, and Module Options.

## 5.2 Instrument Setup mode

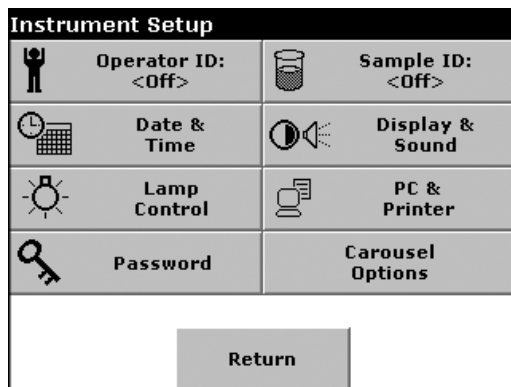

1. Select **Instrument Setup** in the Main Menu.

A selection of functions appears in order to configure the functions of the instrument.

This display appears when the multi-cell holder is installed or the 13 mm round cuvette/vial compartment is used. If the instrument recognizes a Sipper module or a Peltier module the **Carousel Options** key is replaced by **Sipper Options** or **Peltier Options**.


Optional modules for the DR5000:

- Carousel Holder (Sample Changer)
- Pour-Thru Cell module
- Sipper module
- Peltier module (Temperature control module)


Refer to the corresponding manuals/instruction sheets for more information.

### 5.2.1 Operator ID

Use this option to enter up to 30 sets of operator initials (up to 5 characters each) into the instrument. This feature helps record which operator measured each sample.

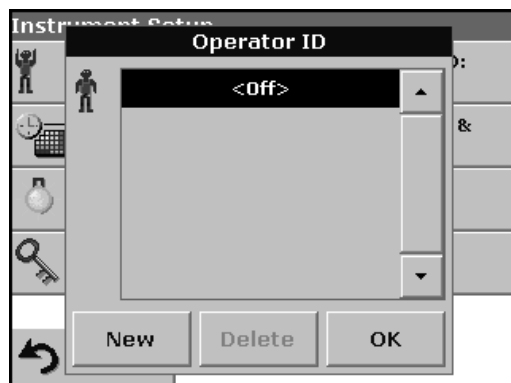

1. Press **Operator ID** in the Instrument Setup.
2. Press **New** to enter a new Operator ID.
3. Use the alphanumeric keypad to enter a new Operator ID.
4. Press **OK** to confirm.

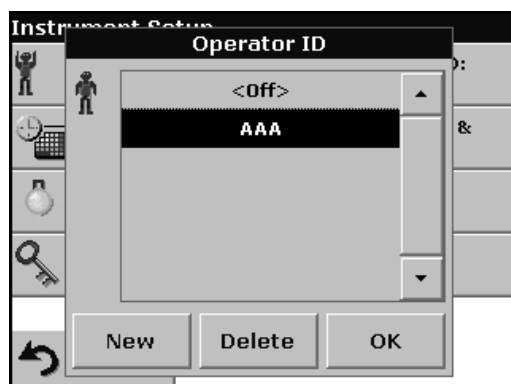


5. The display shows the chosen Operator ID.
6. Press **OK**. The instrument will return to the Instrument Setup screen and show the selected operator identifier.
7. The chosen Operator ID is activated.

**Note:** Press **Delete** to remove an Operator ID from the list.


**Note:** Alternatively, enter or change an Operator ID in measurement mode. In the results screen, press **Options>More>Instrument Setup** or if an Operator ID is already assigned, select the "Operator ID symbol immediately in the results screen.

### 5.2.2 Sample ID


Use this option to enter up to 30 Sample Identification tags (up to 13 characters each) into the instrument. Sample IDs can be used to specify the sample location or other sample specific information.

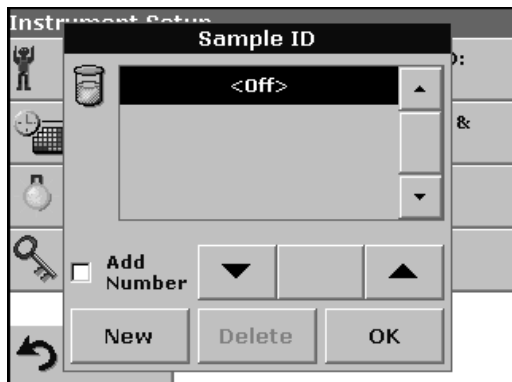

1. Press **Sample ID** in the Instrument Setup.
2. Press **New** to enter a new Sample ID.

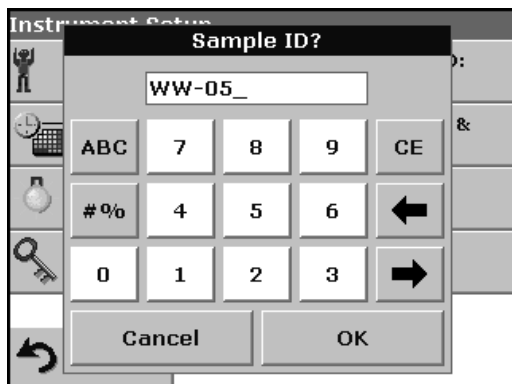

3. Use the alphanumeric keypad to enter a new Sample ID.  
*Note: If a USB Barcode handset scanner ( see Section 9 on page 129) is connected, Sample IDs can also be scanned.*
4. Press **OK** to confirm.

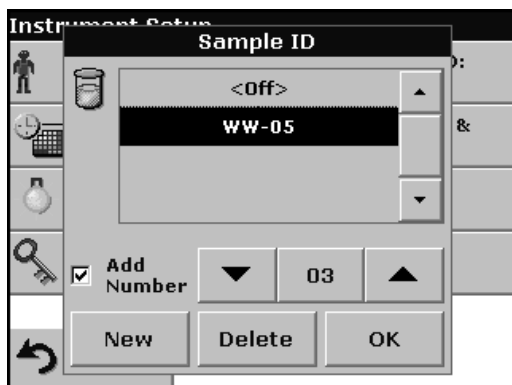

5. To number the Sample IDs sequentially (e.g. Inflow (01 etc.)), select **Add Number**.
  - Use the arrow keys to specify the first number of the sequence.
  - Use the key between the arrow keys to enter the first number of the sequence using the alphanumeric keypad.
6. Press **OK** to return to "Instrument Setup".
7. The Sample ID is activated. Each Sample ID is automatically numbered in ascending order after a measurement. The number is shown in parentheses behind the Sample ID.


To remove a Sample ID, highlight the ID and press **Delete**.



There are two ways to enter or change a Sample ID in the Measurement Mode:

**Option 1**


In the result display, press **Options>More>Instrument Setup**. If a Sample ID is already assigned, select the Sample ID symbol in the result display.

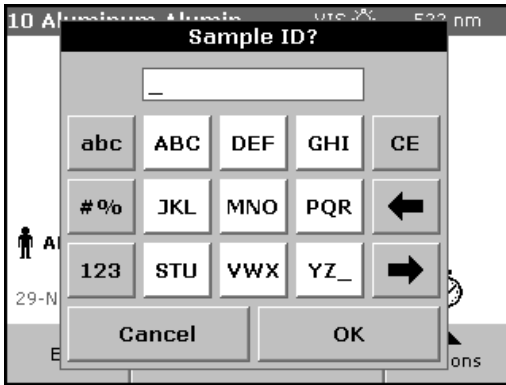

**Option 2**


With an USB keyboard (with US keyboard layout) or a scanner a Sample ID can be directly entered in the result display.

**5.2.3 Date and time**

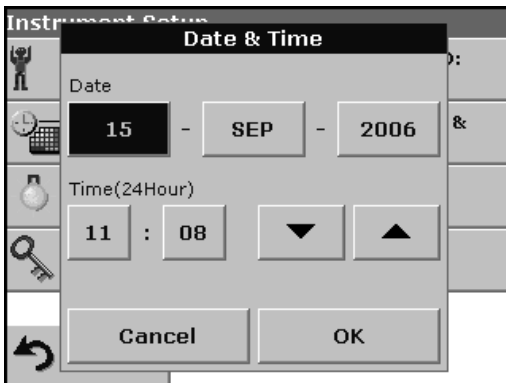

1. Press **Date & Time** in the Instrument Setup.
2. The date and time are subdivided over a number of fields. Press the appropriate field and use the arrow keys to change the value.
3. Press **OK** to confirm. The instrument will return to Instrument Setup.


**5.2.4 Display and sound preferences**

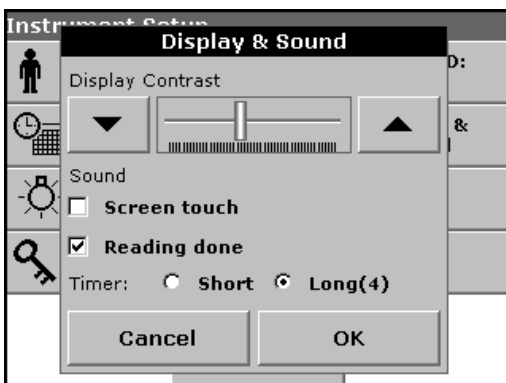

1. Press **Display & Sound** in the Instrument Setup.

The following options will be displayed:

- **Display Contrast**—Adjusts the display brightness to suit lighting conditions.
- **Screen touch**—Activates//Deactivates a short beep each time the screen is pressed (Default:off).
- **Reading done**—Activates/Deactivates a sound when a reading is complete (Default: short beep every time a reading is complete).
- **Timer**—Adjusts the length of the timer sound. Select Short or Long. Long beeps are recommended for noisy environments.

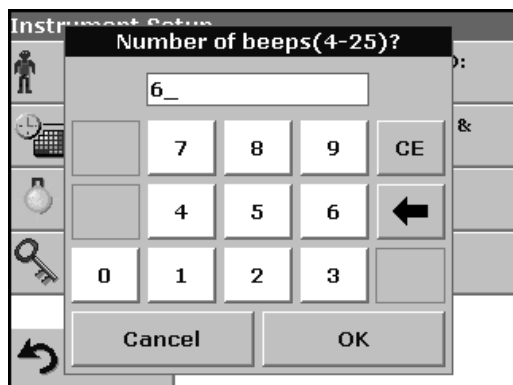

2. Select **Long** to change the number of audio signals.

Use the alphanumeric keypad to enter/specify the number of audio signals (4–25).

**Note:** A high number of audio signals increases the duration of the tones and a small number of audio signals reduces the duration of the tones.

3. Press **OK** to confirm. The selected number of the audio signals sounds as a corresponding acoustic signal.
4. Press **OK** to confirm. The instrument will return to Instrument Setup.

### 5.2.5 Lamp Control

The tungsten lamp produces light in the visible spectrum 320 to 1100 nm.

The deuterium lamp (UV-lamp) available produces light in the ultraviolet spectrum 190 to 360 nm.

In the overlap zone from 320 to 360 nm, either the deuterium lamp (UV-lamp) or the tungsten lamp can be used for measurements.

The lifetime of the lamps is influenced

- by on-off operation and the length of time in use.
- Typical use is to turn the instrument on for the entire 8–10 hour shift, and then off until the next day.

**Note:** In general, avoid on/off cycles of the lamp; this shortens the lamp's life span. For maximum life, turn the lamp off only if it will remain off for at least 4–5 hours.

The lamp switches on automatically

- if a lamp is needed for the selected program
- if the instrument is operating inside the lamp's spectrum

#### 5.2.5.1 VIS-Lamp and UV-Lamp




1. Press **Lamp Control** in the Instrument Setup.
2. Select **On** to switch on the VIS-Lamp and/or UV-Lamp.
3. Select **Save:** in order to define a time interval for the burning time of the VIS lamp and/or UV lamp.
4. Press the field below **Save** to select the lamp burning time.



5. Select the length of time the lamp will be switched on.

**Note:** After this period of time the lamp automatically turns off, after no measurement has been made outside the corresponding range (UV/VIS).

6. Touch **OK** to confirm.

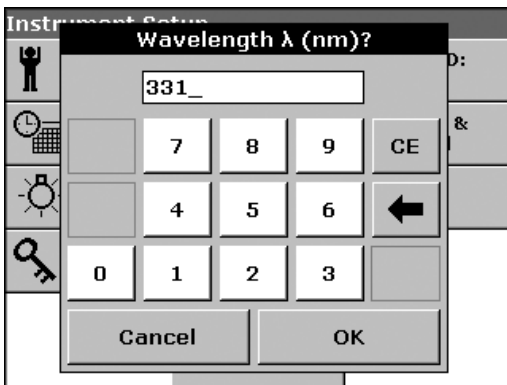

7. Touch **UV Switch** key to select the wavelength value between 320 nm and 360 nm, at which the instrument changes from the visible to UV source.

8. Use the alphanumeric keypad to enter the maximum wavelength for UV operation.

9. Touch **OK** to confirm and return to "Lamp Control".

10. Touch **OK** to confirm and return to "Instrument Setup".

**Note:** The selected program that requires a lamp has highest priority. If the lamp in the **Lamp Control** was turned off, the lamp will be turned on automatically if it is needed for the stored program.

### 5.2.5.2 Effect caused by the lamp settings on the Measuring Mode



UV-Lamp switches on manually. The icon "UV Lamp" appears in the display and flashes.

The instrument will display "Lamp Warmup...". This takes about 3 minutes.



When the lamp is warmed up and ready, the UV icon stops flashing.

**Note:** If both lamps are On, the icon UV-VIS is displayed in the selected Measurement Mode.

**Note:** Alternatively you can touch the UV- or VIS-Lamp icon on the measurement screen to change the actual setting.

## 5.2.6 PC and printer

The instrument is provided with 3 USB interfaces, which are located on the back of the instrument (see Figure 1 on page 14). These interfaces can be used for exporting data and graphics to a printer, updating data and for data communication to a personal computer. These interfaces can be used for the connection of a USB stick, an external USB keyboard or a USB Barcode handset scanner.

**Note:** A USB hub may be used to connect several accessories at a time.

A USB memory stick is used to upgrade data and software, see section 6.8.2 on page 106.

**Important Note:** A screened **USB cable** must not be longer than **3 m!**

**Table 3 USB connector**

USB Interfaces	Description
USB (Type B)	This USB interface is only intended to connect the instrument to a PC (with installation of the HACH Data Trans Software).
USB (Type A)	These USB ports can be used to connect a printer, a USB memory stick, a barcode scanner or a keyboard.



1. Press **PC & Printer** in the Instrument Setup.


A list with information about the connections opens.

### 5.2.6.1 Printer setup



For reasons of compatibility, the printer language must be **HP PCL 3**.

2. Press **Printer**.
3. Press **Setup** to display the Printer Setup screen.

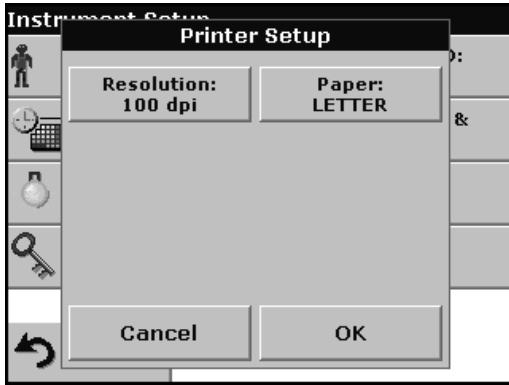

### Printer Setup:


- Resolution: Print quality
- Paper: Paper size

**Note:** If an optional Thermal Printer is connected, the function "Auto Send" on/off is available.



4. Select **Auto-Send: On** to send all measured data automatically to the Thermal printer.

**Note:** The option Auto-Send is **not** available for any other printer (e.g. ink jet printer).

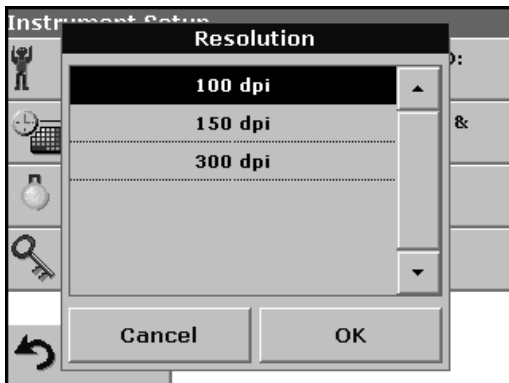


5. Press **Resolution** to select the print quality.

Select between

- 100 dpi
- 150 dpi and
- 300 dpi

6. Press **OK** to confirm.

**Note:** Press **OK** again to return to the **PC & Printer** menu.

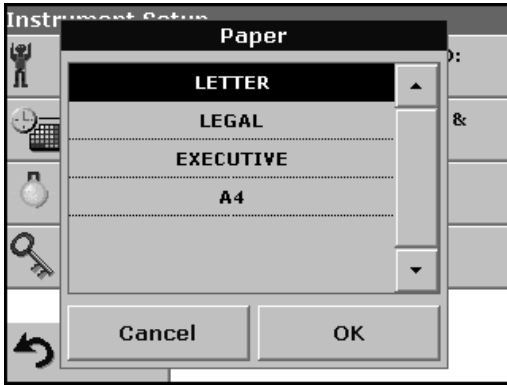

7. Press **Paper** to select the paper size.


Select between

- Letter
- Legal
- Executive
- A4

8. Press **OK** to confirm.

*Note: Press OK again to return to the PC & Printer menu.*

### 5.2.6.2 Print data

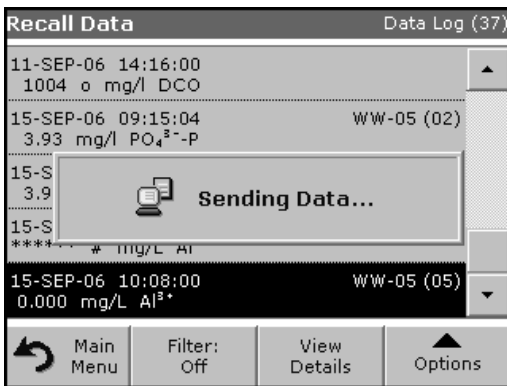

1. Press **Recall Data** in the Main Menu.

2. Select the data source, where the data to be printed are stored.

3. A list is displayed. Data can be filtered. For more information see section 5.3.1.2 on page 35.

4. Press the **PC & Printer** icon to send the data (table, curve) immediately to the printer.

5. Highlight **Single point** or **Filtered data** or **All data** and press **OK** to confirm.

**Sending Data...** is displayed until the data have been printed.

### 5.2.6.3 HACH Data Trans

The optional HACH Data Trans software must be installed on the PC for the subsequent to process for measurement data.




1. Press **PC & Printer** in the Instrument Setup.

2. Select **PC**.

3. Press **Setup** to display the PC Setup screen.

For further installation instructions, refer to the HACH Data Trans user manual.

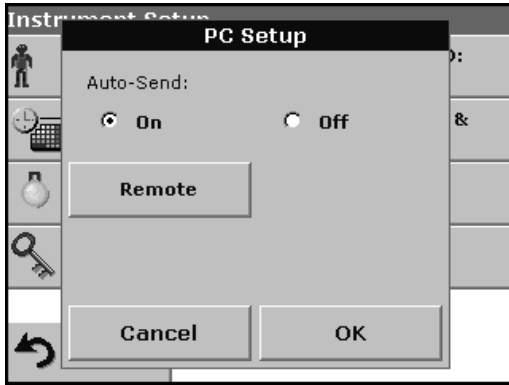

4. Select **Auto-Send: On** to send all measured data automatically to the PC.

*Note: If **Auto-Send: Off** is selected, the **PC & Printer** icon must be pressed, in order to send data to the PC.*


*Note: The remote function is only for monitoring the data transfer.*

### 5.2.7 Password

The Password menu contains a variety of security settings to control access to various functions. For example, prevent unauthorized changes to stored programs or instrument configurations.



1. Press **Password** in the Instrument Setup menu.
2. In order to highlight the **Security List** assign a password. Press **Set Password**.

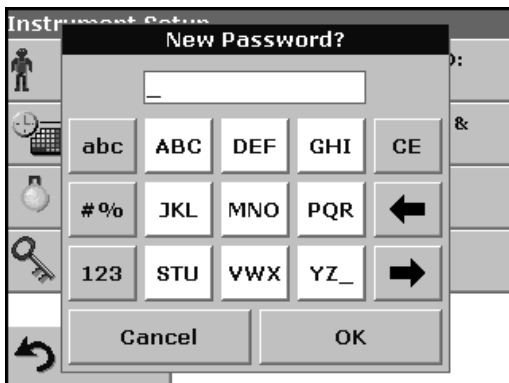

3. Use the alphanumeric keypad to enter a new Password (up to 10 characters each) and press **OK** to confirm.

The access to the Security List is activated.






4. Press **Security List** to lock various functions for unauthorized users.



5. Highlight the desired functions to control.
6. Confirm the **Security List** with **OK** to return to the Password menu.
7. Press **On** to highlight the new settings of the Security List.
8. Enter the new Password again to confirm.
9. Press **OK** to return to Instrument Setup.

*Note: The alphanumeric keypad to the Password inquiry appears when a user tries to reach a locked setting.*

### 5.2.7.1 Password deactivation

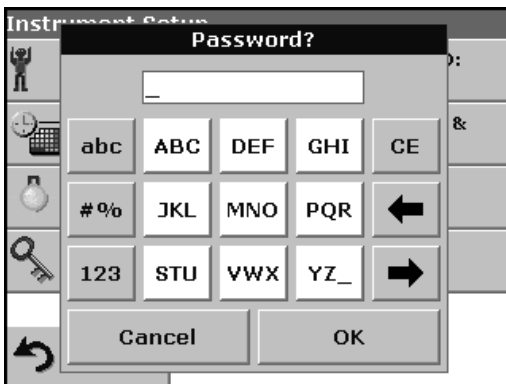

1. Press **Password** in the Instrument Setup.
2. Use the alphanumeric keypad to enter the former Password and press **OK** to confirm.



3. Press **Off** to deactivate the settings of the Security List.

4. Press **OK** to return to Instrument Setup.

**Note:** Use this function to delete the former Password or to enter a new one.


## 5.3 Store, recall, send and delete data

### 5.3.1 The Data Log

The Data Log will store up to 1000 readings taken in the modes: Stored Programs, User Programs, Favorite Programs, Single Wavelength and Multi Wavelength. A complete record of the analysis is stored, including the Date, Time, Results, Sample ID, and Operator ID.

#### 5.3.1.1 Auto/manual data storage

The data storage parameter indicates whether data are to be stored automatically or manually (in which case the user has to decide which data to store).

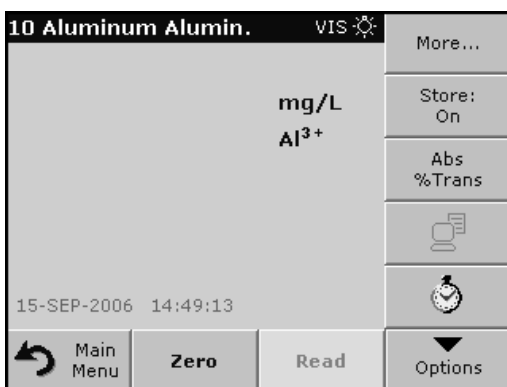


1. Press **Store: On/Off** in the Options menu.

- With the **Store On** setting, all measurement data are stored automatically.
- With the **Store Off** setting, no measurement data are stored. However, this setting can be changed to **Store On** in the result display through configuration. The reading currently shown in the display is then stored.

*Note: When the instrument's memory (data log) is full, the oldest data are automatically deleted allowing the new data to be stored.*

#### 5.3.1.2 Recall stored data from the data log

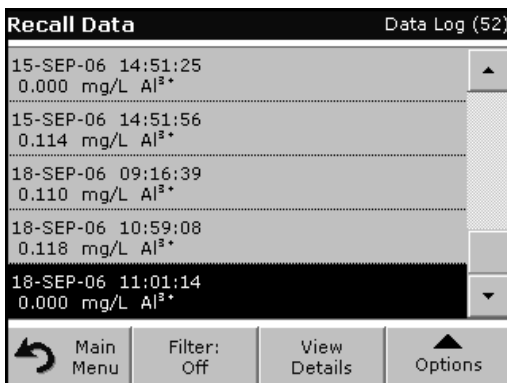

1. Press **Recall Data** in the Main Menu.

2. Press **Data Log**.

A listing of the stored data is displayed.

3. Press **Filter: On/Off**.




4. The function **Filter Settings** is used to search for specific items.

5. Highlight **On** to turn on the filters to select data by

- Sample ID
- Operator ID
- Start Date
- Parameter

or any combination of the four.

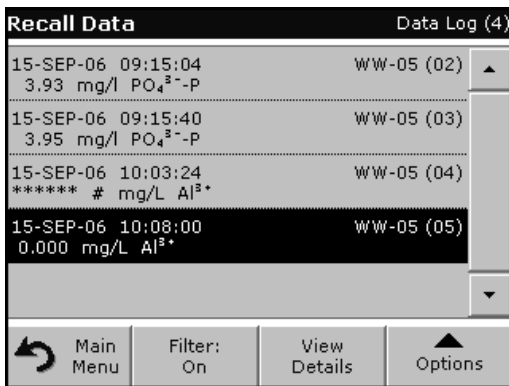

6. Press **OK** to confirm the selection. The chosen items are listed.


7. Press **View Details** to get more information.

### 5.3.1.3 Send data from the data log

Data is sent from the data log as CSV (Comma Separated Value) files through a USB memory stick to a file named DATALOG. The file can then be processed using a spreadsheet program. The file name will be formatted as:

DLYear\_Month\_Day\_Hour\_Minute\_Second. CSV.

To send data to a Printer, see section 5.2.6.2 on page 31.

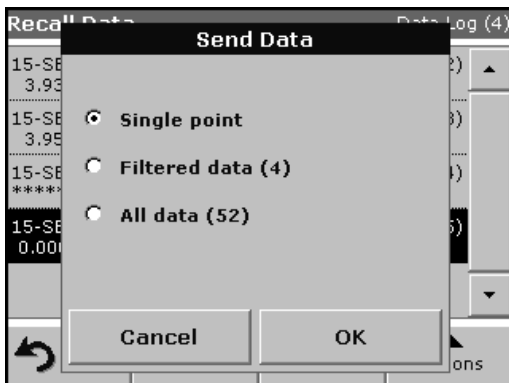

1. Plug in the USB device (Figure 1 on page 14).

2. Press **Recall Data** from the Main Menu. Press **Options** and then the **PC & Printer** icon


3. Select the data to send to the memory stick and press **OK**.

*Note: The number in parenthesis is the total number of data sets assigned to this selection.*

**To send measurement data to a PC:**

The optional HACH Data Trans software must be installed on the PC, see section 5.2.6.3 on page 31.

### 5.3.1.4 Delete stored data from the data log

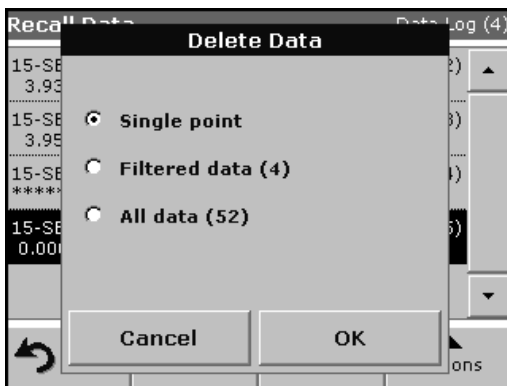

1. Press **Recall Data** in the Main Menu.
2. Press **Data Log>Options>Delete**.
3. Highlight **Single Point** or **Filtered data** or **All data** and press **OK** to confirm.


*Note: The number in parentheses is the total number of data sets assigned to this selection.*

## Standard Operations

### 5.3.2 Store, recall, send and delete data from wavelength scan and time course

The instrument can store 20 Wavelength Scans and 20 Time Course Data sets. The data can be stored manually at the user's discretion after viewing the data.

#### 5.3.2.1 Data storage from wavelength scan or time course

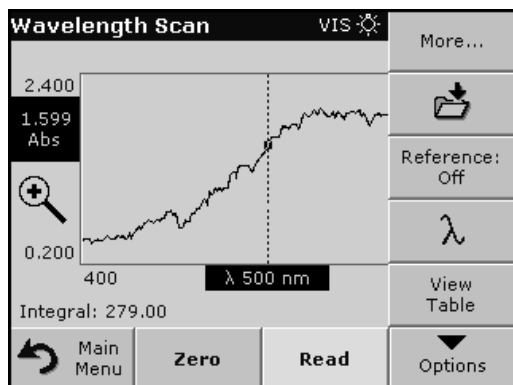

1. Press the **Store icon** in the Options menu after a reading is taken.

Date/Time	Scan	Wavelength Range
15-SEP-06 14:20:33	Scan 7	400 - 600 nm Δ 001 nm
18-SEP-06 09:19:40	Scan 8	400 - 600 nm Δ 001 nm
18-SEP-06 09:26:44	Scan 9	400 - 600 nm Δ 001 nm
18-SEP-06 09:29:14	Scan 10	400 - 600 nm Δ 001 nm
18-SEP-06 09:29:14	Scan 11	400 - 600 nm Δ 001 nm


The Store Data list will be displayed.

2. Press **Store** to save the current scan to the highlighted numbered - line. A scan can also be overwritten.

#### 5.3.2.2 Recall stored data from wavelength scan or time course

Date/Time	Scan	Wavelength Range
15-SEP-06 14:20:33	Scan 7	400 - 600 nm Δ 001 nm
18-SEP-06 09:19:40	Scan 8	400 - 600 nm Δ 001 nm
18-SEP-06 09:26:44	Scan 9	400 - 600 nm Δ 001 nm
18-SEP-06 09:29:14	Scan 10	400 - 600 nm Δ 001 nm
18-SEP-06 10:01:32	Scan 11	400 - 600 nm Δ 001 nm

1. Press **Recall Data** in the Main Menu.
  - a. Select **Wavelength Scan** or **Time Course** to recall data.
  - b. If a program is already in progress, press **Options>More>Recall Data**.

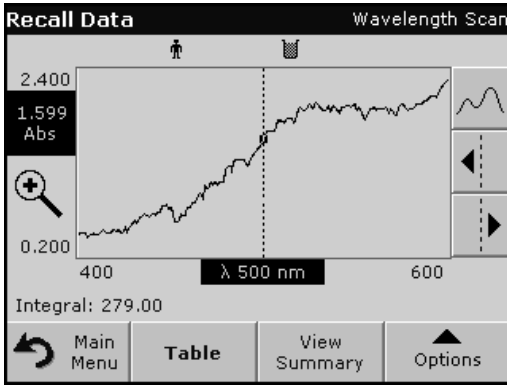

2. Press **Graph** to look at details.

*Note: Press **View Summary** to return to the Recall Data list.*

nm	Abs	Min/Max	nm	Abs	Min/Max
400	0.494		401	0.476	
402	0.504		403	0.500	Peak
404	0.500		405	0.448	
406	0.453	Valley	407	0.472	
408	0.475		409	0.503	Peak
410	0.481		411	0.481	Valley

3. Press **Table** to look at details.

*Note: Press **View Summary** to return to the Recall Data list.*


### 5.3.2.3 Send data from wavelength scan or time course

There are two ways to recall sent data to a USB memory stick, printer or PC with Hach Data Trans.

Date	Time	Scan	Wavelength Range	Action
15-SEP-06	14:20:33	Scan 7	400 - 600 nm Δ 001 nm	
18-SEP-06	09:19:40	Scan 8	400 - 600 nm Δ 001 nm	
18-SEP-06	09:26:44	Scan 9	400 - 600 nm Δ 001 nm	
18-SEP-06	09:29:14	Scan 10	400 - 600 nm Δ 001 nm	Delete
18-SEP-06	10:01:32	Scan 11	400 - 600 nm Δ 001 nm	PC & Printer

**Option 1:**

1. Press **Recall Data** in the Main Menu and then **Wavelength Scan** or **Time Course**.
2. Press **Options** and then the **PC & Printer** icon to send the data to a USB memory stick, to a printer or to a PC with Hach Data Trans.

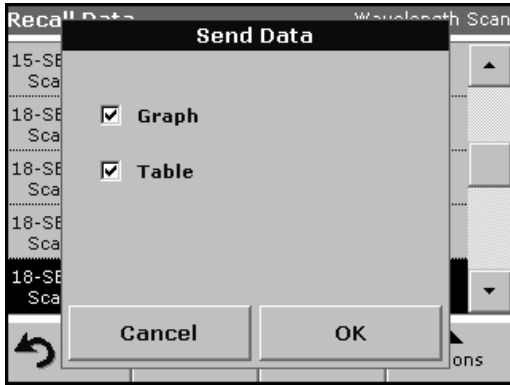


- When a printer is connected, select how to send the data to the printer (graph, table or both graph and table).
- When a USB memory stick is connected, the files will be automatically sent as CSV files (Comma Separated Value) to a file "WLData" (Wavelength Scan Data) or "TCData" (Time Course Data) to the USB memory stick.

The file name will be formatted as: "ScanData\_X.csv" (Wavelength Scan Data) or "TCData\_X.csv" (Time Course Data).

X = number of scans (1–20)

For further processing use a spreadsheet program.

**Note:** The advice "Data already exist. Overwrite?" appears when the files were already stored. Press **OK** to overwrite the stored data.

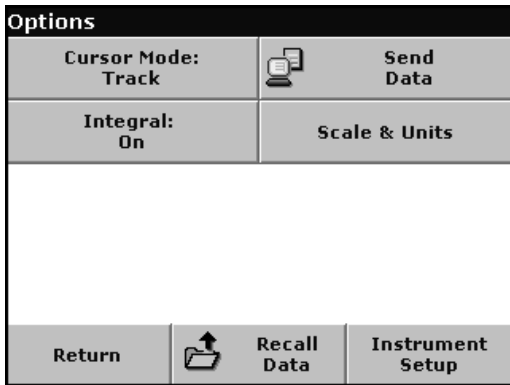

### Option 2:

1. Press **Wavelength Scan** or **Time Course** and then **Options>More>Send Data** to send the data to a USB memory stick or to a printer.

- When a printer is connected, select how to send the data to the printer (graph, table or both graph and table).
- When a USB memory stick is connected, the files will be automatically sent as CSV files (Comma Separated Value) to a file "WLData" (Wavelength Scan Data) or "TCData" (Time Course Data).

The file name will be formatted as:

"ScanData\_Year\_Month\_Day\_Hour\_Minute\_Second.CSV" (Wavelength Scan Data) or



The file name will be formatted as:

"TCYear\_Month\_Day\_Hour\_Minute\_Second.CSV" (Time Course Data).

For further processing use a spreadsheet program.



### 5.3.2.4 Delete stored data from wavelength scan or time course

Recall Data		Wavelength Scan	
15-SEP-06 14:20:33	Scan 7	400 - 600 nm $\Delta$ 001 nm	▲
18-SEP-06 09:19:40	Scan 8	400 - 600 nm $\Delta$ 001 nm	
18-SEP-06 09:26:44	Scan 9	400 - 600 nm $\Delta$ 001 nm	
18-SEP-06 09:29:14	Scan 10	400 - 600 nm $\Delta$ 001 nm	Delete
18-SEP-06 10:01:32	Scan 11	400 - 600 nm $\Delta$ 001 nm	
	Main Menu	Table	Graph
			Options

1. Press **Recall Data** from the Main Menu and then **Wavelength Scan** or **Time Course** or **Options>More>Recall Data**.

A listing of the stored data is displayed.

2. Highlight any data to delete.
3. Press **Delete** in the Options menu and press **OK** to confirm.

## 5.4 Stored Programs

The instrument contains more than 200 programmed procedures. They can be accessed through the **Stored Programs** menu.

### 5.4.1 Select a saved test/method; enter user-specific basic data

Stored Programs		
10	Aluminum Alumin.	0.800 mg/L
9	Aluminum ECR	0.250 mg/L
20	Barium	100 mg/L
771	Beer color	60.0 units
30	Benzotriazole	16.0 mg/L
40	Boron	14.0 mg/L
45	Boron LR	1.50 mg/L
50	Bromine	4.50 mg/L
55	Bromine AV	4.50 mg/L
395	CD 2	6.00 g/l

	Main Menu	Select by Number	Program Options	Start
--	-----------	------------------	-----------------	-------

4. Press **Stored Programs** in the Main Menu to view an alphabetical list of stored programs with program numbers.

The Stored Programs list will appear.

5. Highlight the required test.

**Note:** Select the program number by name or use the arrow keys to scroll through the list quickly and highlight the program or press **Select by number** to search for a specific program number. Use the alphanumeric keypad to enter the test number and press **OK**.

6. Press **Start** to run the program. After a program is selected, the screen for that parameter will appear.

**Note:** All corresponding data (wavelength, factors and constants) are already preset.

7. Follow the chemical procedures described in the corresponding Procedures Manual.

### 5.4.2 Stored program options

1. From the Main Menu, select **Stored Programs**. Select the necessary method and press **Start**.
2. Press **Options** for Parameter Setup. Refer to Table 4 for stored program descriptions.

10 Aluminum Alumin.	VIS	More...
mg/L		Store: On
Al <sup>3+</sup>		Abs %Trans
15-SEP-2006 14:49:13		
	Main Menu	Zero
		Read
		Options

Options	
Send Data	Reading Mode: Single
Dilution Factor: Off	Standard Additions
Standard Adjust: Off	Chemical Form: Al <sup>3+</sup>
Reagent Blank: Off	Save as User Program
Return	Recall Data
	Instrument Setup

Table 4 Stored programs options

Options	Description
More	For further Options


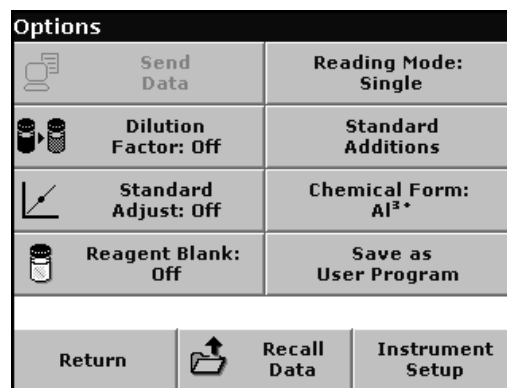


Table 4 Stored programs options (continued)

Options	Description
<b>Store Off/On</b>	With the <b>Store On</b> setting, all measurement data are stored automatically. With the <b>Store Off</b> setting, no measurement data are stored.
<b>% Trans/Conc/Abs</b>	To switch to % transmittance, concentration or absorbance readings
<b>Send Data icon / Send Data</b>	To send Data to a printer, computer or USB memory stick (Type A)
<b>Timer icon</b>	This functions as a stopwatch. It helps to ensure that the steps of an analysis are correctly timed (e.g. reaction times, wait times, etc., can be exactly specified). When the specified time has elapsed, an acoustic signal is emitted. The use of the timer has no influence on the measurement program.
<b>Reading Mode</b>	<b>Single Reading Mode:</b> A reading is only displayed after a measurement has been carried out (press Read; standard setting) (see section 5.4.4.1 on page 45). <b>Continuous Reading Mode:</b> After the zero measurement, all readings are displayed automatically and continuously (see section 5.4.4.2 on page 45).
<b>Dilution Factor Off/On</b>	A corrective dilution factor can be entered in order to take account of certain properties. The number entered at the dilution factor prompt will be multiplied by the result to compensate for the adjustment. For example, if the sample has been diluted by a factor of 2, enter 2. The default setting of the dilution factor is turned off. <b>Note:</b> <i>When a dilution is in effect, the dilution icon will appear on the display.</i>
<b>Standard Addition</b>	This enables the accuracy of the measurements to be checked. The (working) procedure for a test parameter contains a detailed explanation of how to use this function.
<b>Standard Adjust</b>	The (working) procedure for a test parameter indicates whether a standard adjustment is necessary and, if so, how to proceed.
<b>Chemical Form</b>	Some of the stored tests/methods allow to select the chemical form and the measuring range.
<b>Reagent Blank</b>	Some of the stored tests/methods include the "Reagent Blank" function. This enables the reagent blank value to be added to or subtracted from, the subsequent readings. The reagent blank value shifts the calibration curve along the y-axis, without changing the shape or gradient of the curve. The effect corresponds to a y-axis intercept of the calibration straight line. This is made clear by the following equation: Concentration = [(Conc. factor) * Abs] – (reagent blank value).
<b>Save as User Program</b>	To store the selected parameters as a User Program, see section 6.1 on page 61.
<b>Recall Data</b>	Call up saved measurement data, wavelength scans or time courses, see section 5.3 on page 35.
<b>Instrument Setup</b>	Basic data of the instrument, see section 5.2 on page 23.

### 5.4.3 Use of program timers

Some procedures do not require the use of timers. Other procedures require several timers. These timers are pre-programmed into each **Stored Program**, along with a description of the activity to be performed during the timed period.

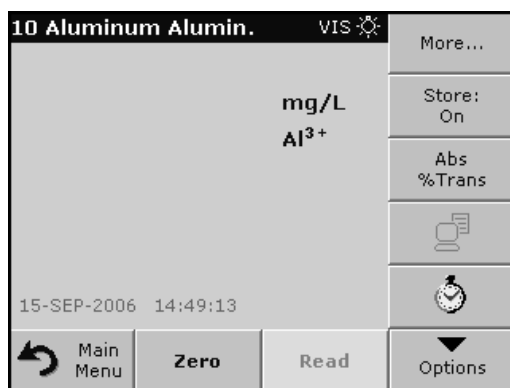

1. Press the **Timer** icon on the display.
2. Press **OK** to start the first timer.  
  
The timer will count down on the screen.
3. To start the next timed activity for the Stored Program, press the Timer icon and **OK**.

**Note:** Press **Close** to view the measurement screen while the timer is running. The time will be shown left side bottom instead of the date.


**Note:** Press **Cancel** to stop the timer midway through the countdown.

**Note:** The Timer will beep when the end of the running time is reached.




A general purpose timer is also available in many programs. When the timer icon is visible, press the icon and select **General Timer**. A new screen will appear. Enter the length of the timed interval and press **OK** to start the timer. The timer will beep when the timed interval ends.

### 5.4.4 Set the reading mode

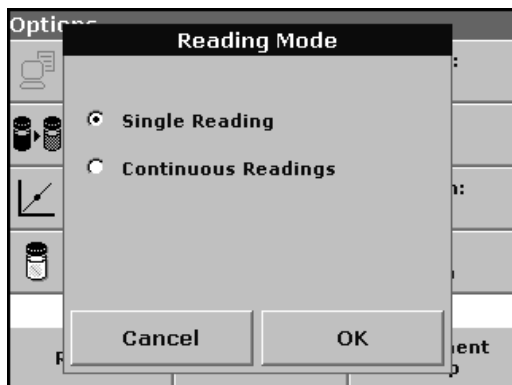

1. To highlight the required mode, press **Reading Mode**.
2. Select the required mode, then press **OK**, then **Return** to return to the result display.

### 5.4.4.1 Take single wavelength measurements (single reading)

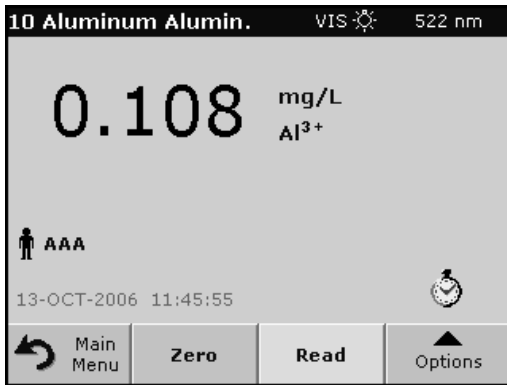


1. Insert the blank cuvette into the cuvette/sample cell holder. Press **Zero**.

*Note: The **Read** key is only active after the zero measurement has been carried out.*

2. Insert the sample cuvette into the cuvette/sample cell holder. Press **Read**.

*Note: For data storage, see section 5.3.2 on page 38.*

### 5.4.4.2 Take single wavelength measurements (continuous readings)

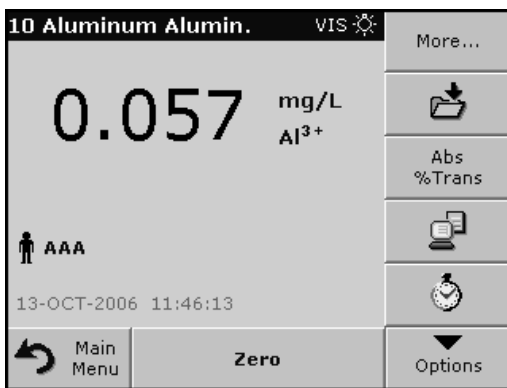

1. Insert the blank cuvette into the cuvette/sample cell holder. Press **Zero**.

*Note: In the reading mode "Continuous" only the Zero key is shown to start the reading. The reading sequence is started automatically.*

2. Insert the sample cuvette into the cuvette/sample cell holder.
3. Press **Options** and then the **Store** icon to store the displayed data in the Data Log.

*Note: For data storage, see section 5.3.2 on page 38.*

### 5.4.5 Set the dilution factor

The Dilution Factor function is used to calculate the original concentration of a sample that has been diluted by a known ratio. For example, dilution factor can bring the analyzed concentration within the test range.



1. Press **Options>More...>Dilution Factor**.

The number entered at the dilution factor prompt will be multiplied by the result to compensate for the adjustment.

For example, if the sample has been diluted by a factor of 2, enter 2. The default setting of the dilution factor is turned off.

2. Press **OK** to confirm.



3. Press **Return** to return to the result display.
4. Confirm 'Store present reading again?' to save the current measured value with the dilution factor.




**Note:** When a dilution is in effect, the dilution icon will appear on the display.

**Note:** If using undiluted samples, set the dilution factor off again.

### 5.4.6 Run a standard adjust

The Standard Adjust function allows the calibration curve for a stored program to be adjusted based on analysis of a known standard solution. The Accuracy Check section of written procedures often suggests a standard solution concentration for this purpose.


Read a standard before setting Standard Adjust to **On**.

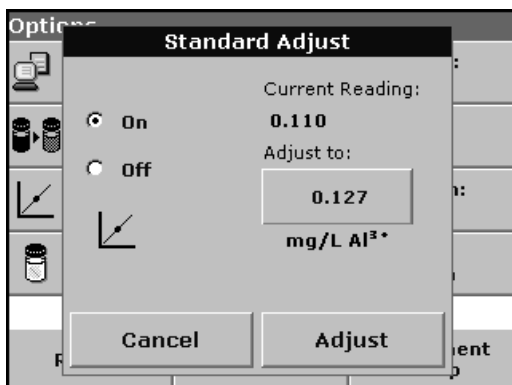

1. Follow the entire procedure, using a known standard for the sample.
2. After reading the concentration, press **Options>More>Standard Adjust**.
3. If Standard Adjust is set to **Off**, turn it **On**.

The Current Reading will show the concentration. The box on the right will show the default standard value for the test, as mentioned in the procedure.

4. If the measurement used a standard concentration that is different from the one displayed in the box, press the box on the right to enter a different standard value and enter the new value. Press **OK** to confirm.

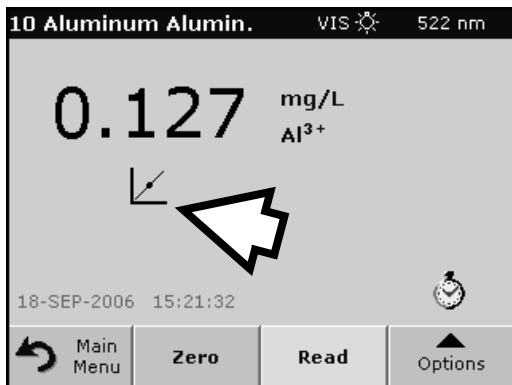

5. Press **Adjust** to enable the Standard Adjust. The Standard Adjust icon will appear.

*Note: The adjustment must be within certain limits, which vary with each program. The allowable percentage is shown after "Adjustment".*

*Note: When a Standard Adjust is in effect, the Standard Adjust icon will appear on the display.*


#### 5.4.7 Set the chemical form

Some Stored Programs allow a variety of chemical forms to be selected.

Press the unit (e.g. mg/L) or the chemical representation of the evaluation form (e.g. Al<sup>3+</sup>). A list of available evaluation forms is displayed. Select the required form by pressing the corresponding entry in the list.

*Note: When the program is exited, the evaluation form reverts to the standard setting.*

An alternative way of changing the standard setting:

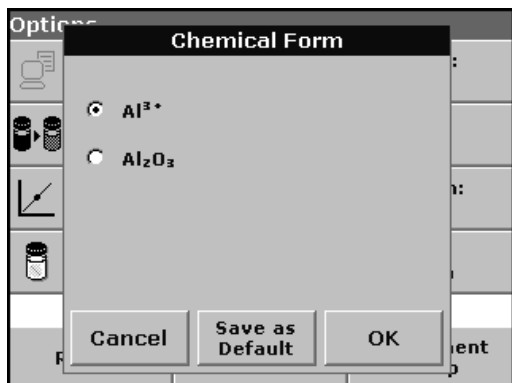

1. Press **Options>More>Chemical Form**.
2. Select the Chemical Form.

*Note: The stoichiometric conversion of the measurement result is carried out automatically.*

*Note: The selected Chemical Form will appear on the display. Test results will be calculated and displayed in this chemical form.*

##### 5.4.7.1 Change of the default setting of the chemical form

1. Insert the sample cuvette or blank (depending on the working procedure) into the cell compartment.
2. In the result display, press **Options>More>Chemical Form**.
3. A list of available evaluation forms appears. Select the new default setting.
4. Press **Save as Default**.


The current result and all further measurements will be displayed in the new chemical form.

### 5.4.8 Run a reagent blank

Some of the stored tests/methods include the "Reagent Blank" function. This enables the reagent blank value to be measured and then taken into account in calculating the measurement result.

#### Measurement/analysis of a reagent blank:

1. Prepare the test/method in accordance with the (working) procedure. Instead of a sample, deionized water is used to determine the reagent blank value.
2. Select the test. If required by the (working) procedure, insert the cell with deionized water into the cell compartment. Press **Zero**.
3. Insert the prepared sample cell into the cell compartment. Press **Read**. The result is displayed.

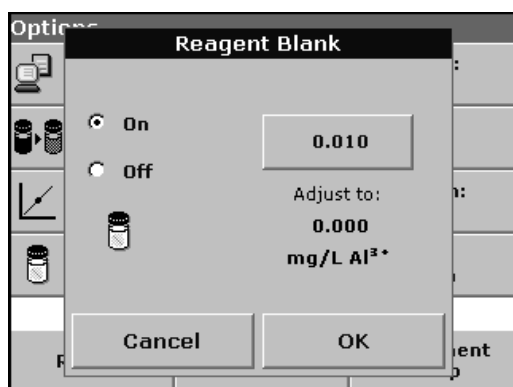


4. Press **Options>More>Reagent Blank**.
5. Press **On** to highlight the Reagent Blank function.
6. The concentration shown on the key is the measured value of the reagent blank. To use this value for more analyses of this parameter, press **OK**.

To enter a previously recorded reagent blank value, press the key and use the alphanumeric keypad.

7. Press **OK**.

**Note:** The Reagent Blank function is deactivated when the measurement program is left. To use the same blank value later for other tests using the same reagent lot, enter the value per step.


**Note:** The results calculated using the reagent blank value must lie within the limits of the measuring range of the test/method.

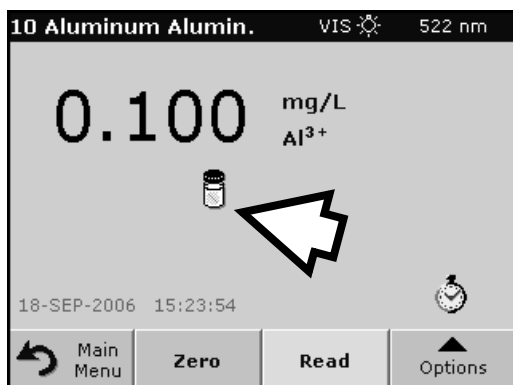


**Note:** The reagent blank icon is shown in the result display (see arrow) when the function is active.



### 5.4.9 Analysis of samples

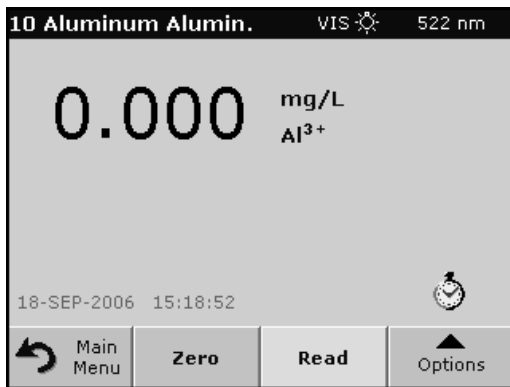

1. Press **Stored Programs** and select a program.
2. Insert the blank cuvette/cell into the cuvette/sample cell holder.
3. Press **Zero**.

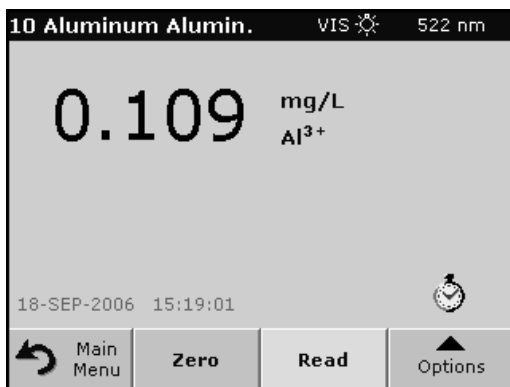

4. Remove blank cuvette/cell and insert sample cuvette/cell into the cell compartment.
5. Press **Read**. The result will be displayed.
6. For data storage, see section 5.3.2 on page 38.

### 5.4.10 Add stored programs to the favorite programs list

The Favorites menu simplifies test selection by creating a list of the most frequently used tests from the Stored Programs and User Programs.



1. Press **Stored Programs** in the Main Menu. The Stored Programs list will appear.
2. Highlight the selection by pressing it or **Select by Number** to search for the program by number.
3. Press **Add to Favorites** and press **OK** to confirm.

The program can now be selected from **Favorite Programs** menu in the Main Menu.

## 5.5 Barcode Programs

A special barcode reader in cell compartment #1 automatically reads the barcode on the 13 mm cuvette/vial as the cuvette/vial completes a single rotation. The instrument uses the barcode identification to automatically set the correct wavelength for the analysis and calculates the result immediately with the help of the stored factors.

In addition, measured values are recorded at 10 different positions during a rotation. A special outlier-elimination program is run and then the average of the measured values is calculated. Cuvette/vial errors and soiling are recognized and highly precise results are obtained.


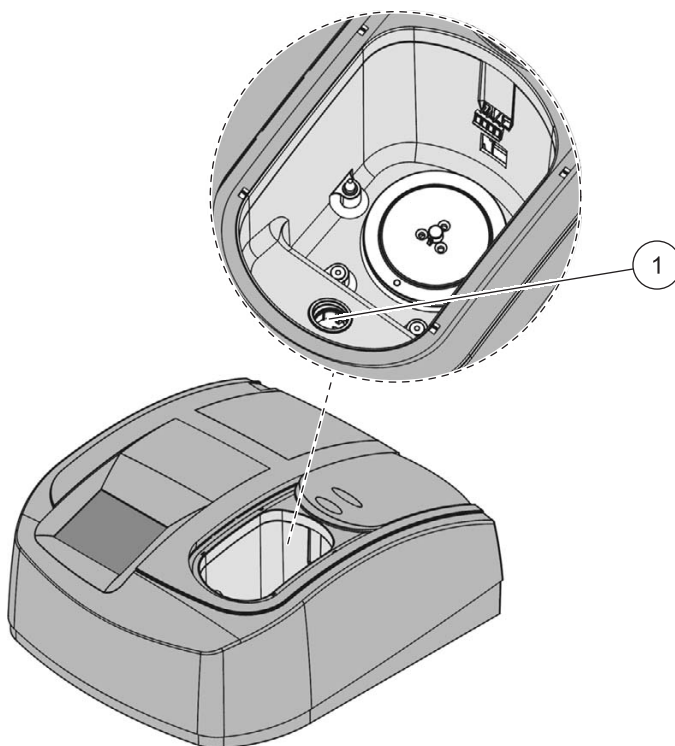


Figure 7 Cell compartment #1

1 Cell compartment #1 for 13 mm barcode cuvettes/or 16 mm vials


### 5.5.1 Complete a barcode 13 mm test/vial

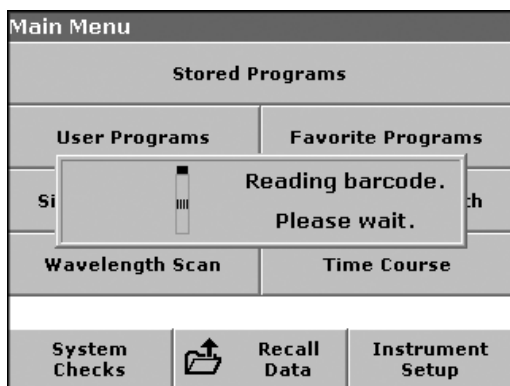

**Note:** The cell compartment must be closed before a measurement is made!

1. The instrument needs to be at the Main Menu.
2. Prepare the barcode 13 mm cuvette/vial in accordance with the working procedure and insert the cuvette/vial in cell compartment #1 and close the cell compartment.


When a coded cuvette/vial is placed in cell compartment #1 (Figure 7), the corresponding measurement program is automatically activated in the Main Menu.

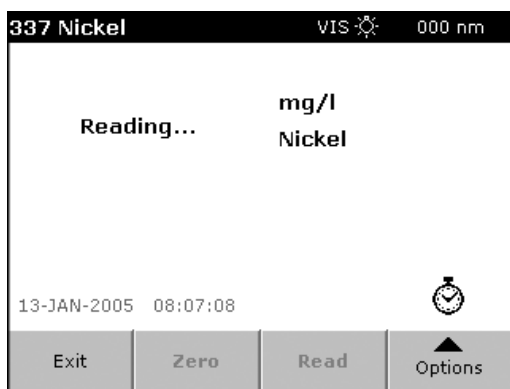

The measurement is started automatically and the results are displayed.

To evaluate other cuvette/vial tests and other parameters, insert the prepared cuvette/vial into the cell compartment and read the result.

### 5.5.2 Select the measuring range

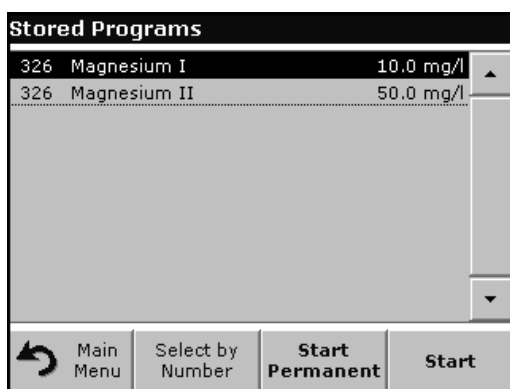

Some tests can be used for different measuring ranges. After the sample cuvette/vial has been inserted, a list of the different measuring ranges is displayed.

Select the required measuring range by highlighting the appropriate line.

Press **Start Permanent** if this measuring range is to apply to all subsequent measurements.

#### Changing the standard setting


In the result display, press **Options>More>Permanent: On**. The key will change to **Permanent: Off**.

### 5.5.3 Select the chemical evaluation form

The chemical form of the test result of a number of parameters can be selected individually.

In the result display, press the unit (e.g. mg/L) or the chemical representation of the evaluation form (e.g.  $\text{PO}_4^{3-}\text{-P}$ ). A list of possible evaluation forms is displayed, from which the required form can be selected. Press **OK** to confirm.

Another way of changing the standard setting is:

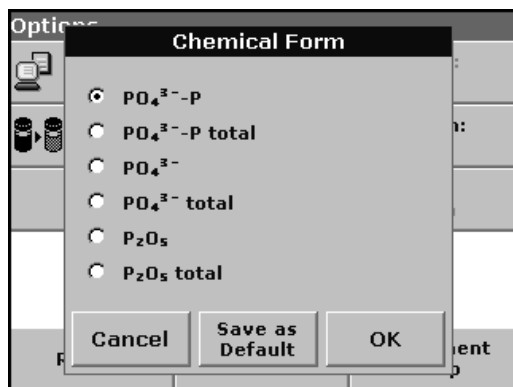

1. In the result display, press **Options>More>Chemical Form**.

A list of available evaluation forms appears.

2. Select the required chemical form and press **OK** to confirm.

*Note: The selected chemical form is displayed, but does not become the default. To change the default, see section 5.5.3.1.*



### 5.5.3.1 Change of the default setting of the chemical form

1. Insert the blank or sample cuvette/vial (depending on the working procedure) into the cell compartment and close the cell compartment.
2. In the result display, press **Options>More>Chemical Form**.
3. A list of available evaluation forms appear. Select the new default setting.
4. Press **Save as Default**.

The current result and all further measurements will be displayed in the new chemical form.

### 5.5.4 Basic test-specific and sample-specific data settings

Press **Options** to change test or sample specific settings.

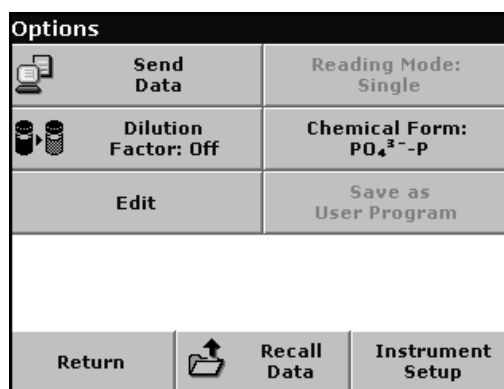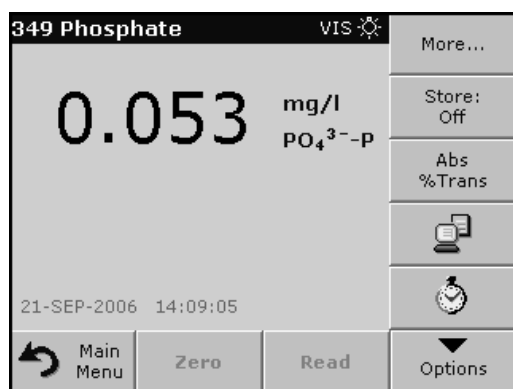

**Table 5 Barcode program options**

Options	Description
More	For further Options
Store Off/On	With the <b>Store On</b> setting, all measurement data are stored automatically. With the <b>Store Off</b> setting, no measurement data are stored.
Abs % Trans	To switch to % transmittance or absorbance readings



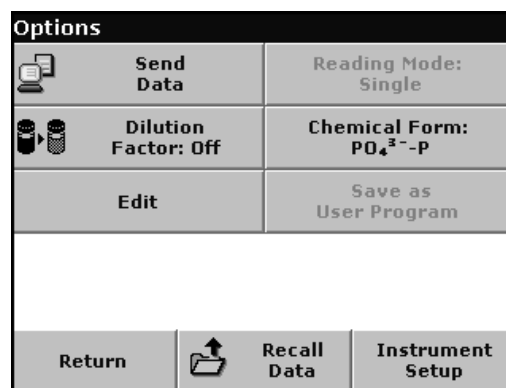


Table 5 Barcode program options (continued)

Options	Description
Send Data icon / Send Data	To send data to a printer, computer or USB memory stick (Type A)
Timer icon	This functions as a stopwatch. It helps to ensure that the steps of an analysis are correctly timed (e.g. reaction times, wait times, etc., can be exactly specified). When the specified time has elapsed, an acoustic signal is emitted. The use of the timer has no influence on the measurement program.
Dilution Factor Off/On	A corrective dilution factor can be entered in order to take account of certain properties. The number entered at the dilution factor prompt will be multiplied by the result to compensate for the adjustment. For example, if the sample has been diluted by a factor of 2, enter 2. The default setting of the dilution factor is turned off. <b>Note:</b> When a dilution is in effect, the dilution icon will appear on the display. <b>Note:</b> If undiluted samples are used, set the dilution factor off.
Chemical Form	Some of the stored tests/methods allow to select the chemical form and the measuring range.
Edit	To modify an existing program
Save as User Program	To store the selected parameters as a User Program, see section 6.1 on page 61.
Recall Data	Call up saved measurement data, wavelength scans or time courses, see section 5.3 on page 35.
Instrument Setup	Basic data of the instrument, see section 5.2 on page 23.

### 5.5.5 Sample blank

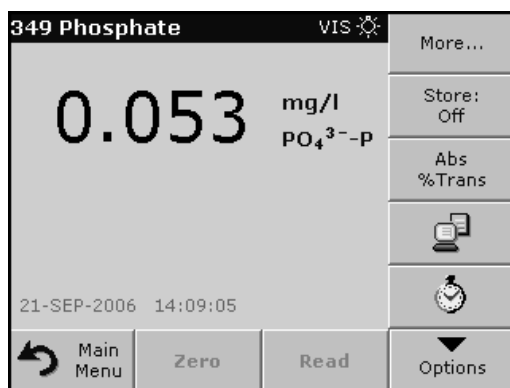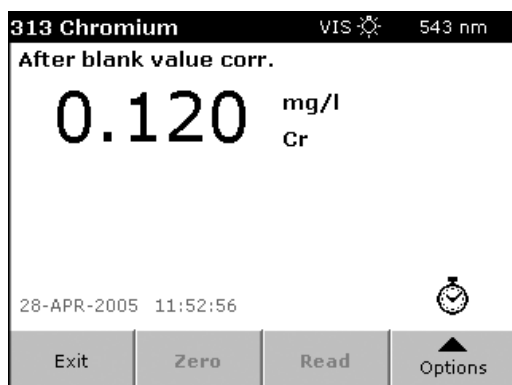

Turbidity and color in the sample matrix can falsify the results of a photometric analysis. The interference factors come from the sample or are created by reactions with the reagents.

The influence of turbidity and/or color can be eliminated or reduced by taking a sample blank reading.


In the barcode mode, a special cuvette/vial (TNT919) containing the sample blank is placed in cell compartment #1 after the sample reading has been taken and is automatically measured. The sample reading is then corrected by adding or subtracting the blank value. The final result is displayed, with the message "After blank value corr.".

Some barcode tests do not require a sample blank value to be determined, as turbidity and color are dealt with during the test

procedure. Other barcode tests use the reagent test vials to determine the sample blank. Please refer to the Hach procedure for method-specific instructions.

## 5.5.6 Update/edit barcode tests

### 5.5.6.1 Manual update of a barcode test

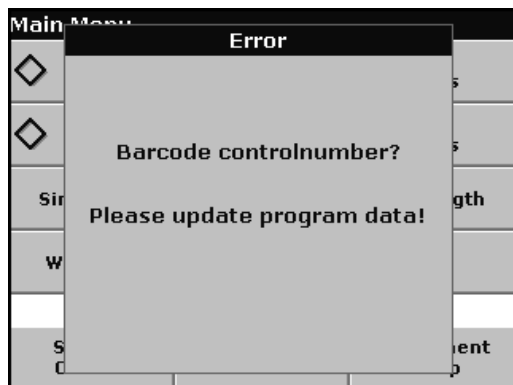


Using the data provided in the barcode, the instrument automatically sets the measurement wavelength and factors. If a discrepancy is detected between the barcode data and the stored data or a new test is identified, the instrument requests an update.

Data types are described in Table 6.

**Table 6 Program update options**


Program point	Description
<b>Name</b>	Name of the analysis parameter
<b>Version</b>	An abbreviation assigned by the user or the version number, etc. is entered here
<b>Measurement process</b>	Exact definition of the test: the number of wavelengths at which measurements are carried out, the number of absorbance measurements needed, the keys to be used, any waiting periods between measurements, etc.
<b>Formula</b>	Definition of the evaluation formulas with which the test result is calculated
<b>Variables</b>	The number of variables shown in the display depends on the definition of the measurement process and the formulas. Input of the numerical values of the wavelengths, factors, constants, etc.
<b>Timer 1, Timer 2, Timer 3, Timer 4</b>	This function can be used to enter abbreviations and defined times for up to four timers. Highlight the appropriate line and press <b>Edit</b> . The timers are activated or deactivated with the control boxes on the left of the display. In the next column, a selection can be made from a list of names that describe the corresponding work step. In the third column, the times for each active timer are entered.

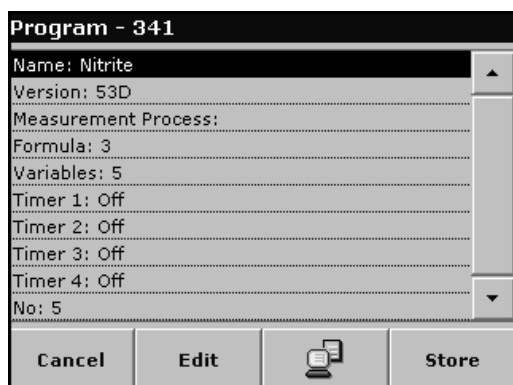

#### Enter the Name


Check the working procedure in advance to determine whether this point needs to be changed.

1. Highlight the line containing **Name** and press **Edit**. Use the alphanumeric keypad to enter the name indicated in the working procedure.
2. Press **OK** to confirm.

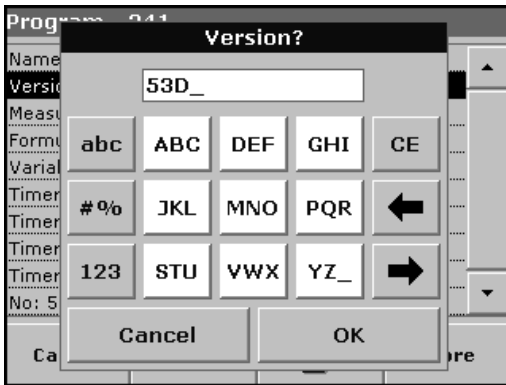

### Enter the Version

1. Highlight the line containing Version and press **Edit**. Use the alphanumeric keypad to enter the name indicated in the working procedure.
2. Press **OK** to confirm.

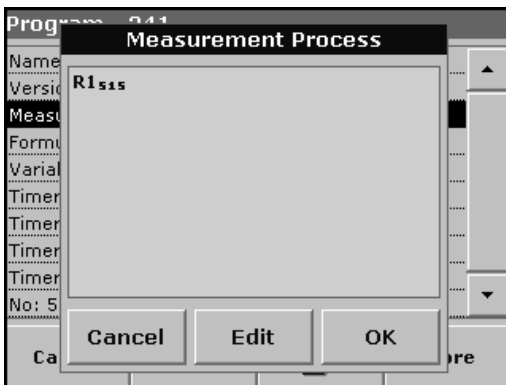


### Enter the Measurement Process

For detailed information on entering the measurement process, see section 6.1.2 on page 72.

Check the working procedure in advance to determine whether this point needs to be changed.

1. Highlight the **Measurement Process** line in the data overview and press **Edit**.
2. Press **Edit**, select the sequence to be edited and press **Delete**.
3. Press **New** and use the alphanumeric keypad to enter the process indicated in the working procedure.

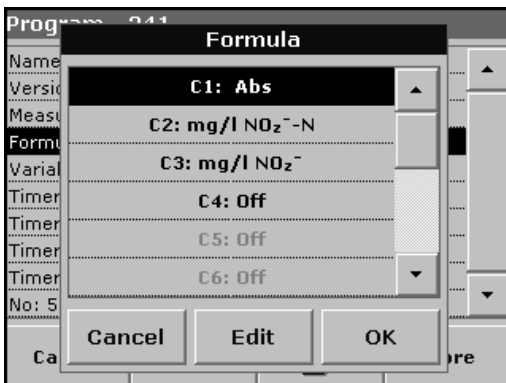


### Enter the Formula, Concentration Units, Designation, Measuring Ranges

For detailed information on entering the Formula, see section 6.1.2 on page 72.

Check the working procedure in advance to determine whether this information needs to be changed.

1. Highlight the **Formula** line in the data overview and press **Edit**.
2. Select the formula, press **Edit** and use the alphanumeric keypad to enter the data specified in the working procedure (for C1=, C2=, units, name, measuring range limits, etc.).

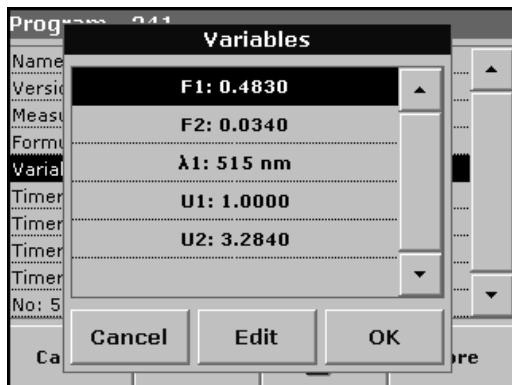

### Enter the Variables (Factors, Wavelength and Conversion Factors)

For detailed information on entering the Variables, see section 6.1.2 on page 72.

Check the working procedure in advance to determine whether this point needs to be changed.

1. Highlight the **Variables** line in the data overview and press **Edit**.
2. Select the variable, press **Edit** and use the alphanumeric keypad to enter the data specified in the working procedure (for F1, F2,  $\lambda$ 1, U1 etc.). Confirm each entry with **OK**.

Abbreviation of Variables:

F1: Factor 1

F2: Factor 2

$\lambda$ 1: Wavelength 1

U1: Conversion Factor 1 for the first chemical form

U2: Conversion Factor 2 for additional chemical forms etc.




### Timer 1, Timer 2, Timer 3 and Timer 4:

The timers are activated or deactivated with the control boxes on the left of the display. In the next column, a selection can be made from a list of names that describe the corresponding work step. In the third column, the times for each active timer are entered.



## 5.5.6.2 Update an existing barcode test

**Note:** Only in very few cases does the revision of a test require all test specifications to be updated.




Program - 341	
Name: Nitrite	
Version: 53D	
Measurement Process:	
Formula: 3	
Variables: 5	
Timer 1: Off	
Timer 2: Off	
Timer 3: Off	
Timer 4: Off	
No: 5	
Cancel	Edit
	
Store	

**Option 1: Manual update of test data**

1. The instrument needs to be at the Main Menu.
2. Insert the blank or sample cuvette/vial (depending on the working procedure) in cell compartment #1 and close the cell compartment.
3. After "Barcode controlnumber?", a data overview is displayed, including the specifications of the test that is to be revised.

**Note:** The working procedures of the corresponding test contains the new data (Wavelength, Factors, Measurement Ranges, Conversion Factor for chemical Form, etc.).

4. Highlight the corresponding line and press **Edit**.
5. When finished programming, press Store and then **Cancel** to return to the User Programs menu.
6. Press **Start** to complete the test with the new data.

Options	
 Send Data	Reading Mode: Single
 Dilution Factor: Off	Chemical Form: $\text{PO}_4^{3-}\text{-P}$
Edit	Save as User Program
Return	 Recall Data
Instrument Setup	

**Option 2: Manual check/revision of test data**


If an incorrect entry was made during the data update and is not recognized via the barcode of the cuvette/sample vial (e.g. an incorrect measuring range), another option is available to check and if necessary to update, the test data.

1. The instrument need to be at the Main Menu.
2. Insert the blank or sample cuvette/vial (depending on the working procedure) into the cell compartment #1 and close the cell compartment.
3. Press **Options>More>Edit**.
4. A data overview is displayed, including the specifications of the test that is to be revised. Compare the displayed data with the data in the working procedure and edit the displayed data in accordance with the working procedure.

### 5.5.6.3 Program a new test



1. The instrument needs to be at the Main Menu.
2. Insert the sample cuvette/vial or blank (depending on the working procedure) into the cell compartment #1 and close the cell compartment.

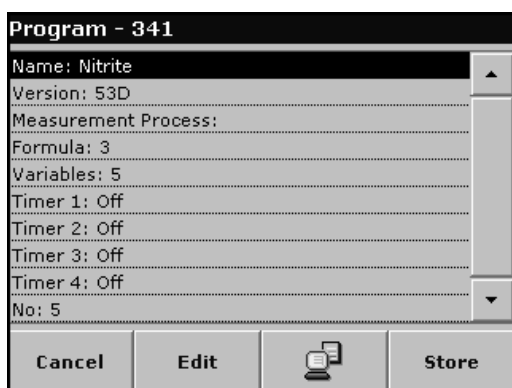

3. After "Program not available", a data overview is displayed, including the specifications of the test that is to be revised.  

The working procedures of the corresponding test contains the new data (Wavelength, Factors, Measurement Ranges, Conversion Factor for chemical Form etc.).
4. Highlight the corresponding line including the specifications of the test that is to be programmed and press **Edit**.
5. To finish programming, press **Store** and then **Cancel** to return to the Main Menu.
6. Press **Start** to run the test with the new data.

### 5.5.7 Upgrade of the instrument software

To obtain the software for the update from the Internet at [www.hach.com](http://www.hach.com):



1. Go to <http://www.hach.com>.
2. On the DR 5000 product page, click **Lab System Software/Software Update Downloads** under the heading **Downloads**.
3. Locate the appropriate download and follow the prompts for saving the file(s) to the USB memory stick.
4. On the DR 5000, press **Instrument Update** in the System Checks menu.
5. Connect the USB stick to the USB interface on the DR 5000 (section 3.3 on page 14). Press **OK**. The link is established automatically and the software is updated.

6. Press **OK** to return to the System Checks menu.

**Note:** *When the instrument software has been updated, a prompt to restart the instrument is displayed.*



# Section 6 Advanced Operations


## 6.1 User Programs

User programs provide the opportunity to complete "made to measure" analysis.

The User Programs database is empty when the instrument leaves the factory and is used to accommodate programs created by users specific needs. Here are a few examples of entries:

- Programming of user-created procedures. The analysis procedure must be developed first, before it can be programmed. The user must define or determine the program sequences, calculation formulas, measurement wavelengths, factors, measuring range limits, etc.
- Modified tests
- Assignment of user programs to the favorites menu for frequently used tests.
- Creation of a specific selection of methods and tests.

Press **User Programs** in the Main Menu and then **Program Options**. The **Program Options** menu contains several input and editing options (Table 7):

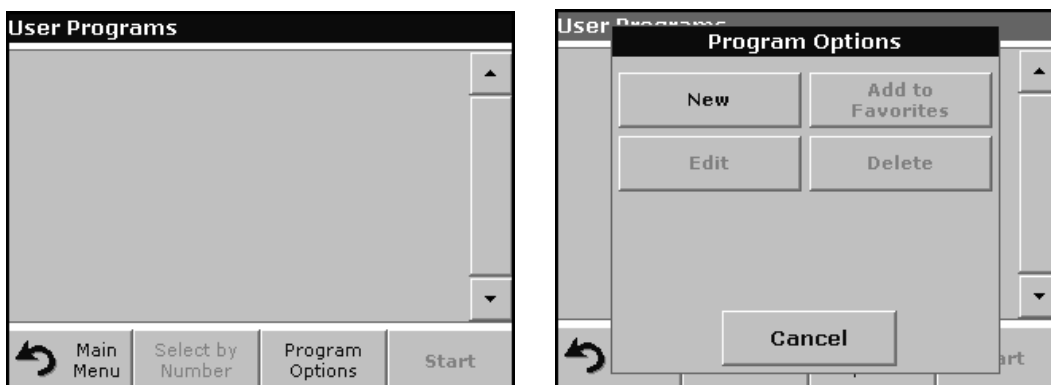


**Table 7 Program options user programs**

Options	Description
1 <b>New</b>	Select New to program a new user program. <i>Note: The first time <b>Program Options</b> is selected, only the <b>New</b> option is available. The other options remain inactive (gray) until the first program has been created.</i>
1 <b>Add to favorites</b>	Select <b>Add to favorites</b> to add an existing user program to the list of frequently used programs.
1 <b>Edit</b>	Select <b>Edit</b> to modify an existing program
1 <b>Delete</b>	Select <b>Delete</b> to remove a program from the list of user programs. The program will be simultaneously deleted from the Favorites list.

### 6.1.1 Program a user method

All input steps and their significance and options are explained in the following sections.

1. Select **New** in the Program Options menu.

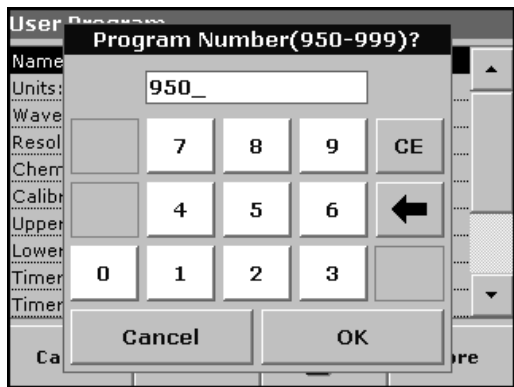


### Program Number:

Specific test number, with which the program can subsequently be called up from the selection list in the **User Programs** menu or the **Favorites** menu.


2. Use the alphanumeric keypad to enter a program number between 950 and 999. The lowest available number appears automatically.
3. Press **OK**.

*Note: If the program number is already assigned to another user program, a message appears, asking whether the existing program should be replaced. Press **OK** to overwrite the existing program.*

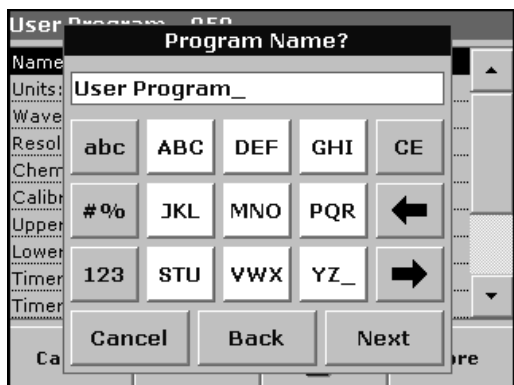

### Program Name:

4. Use the alphanumeric keypad to enter a program name. The name can be a maximum of 28 characters long.
5. Press **Back** to go back to the previous program point or press **Next** to continue with the input of the program data.

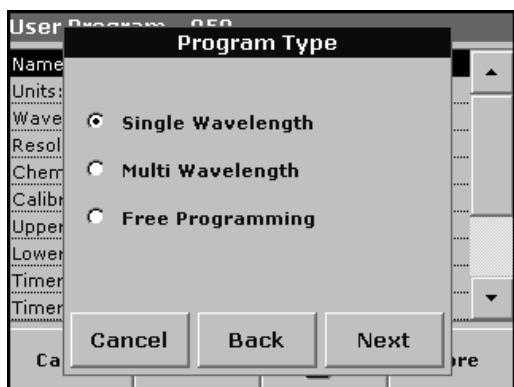

### Program Type:


6. Select the required option (Table 8) and press **Next**.
7. If the Single Wavelength (section 6.1.1.1 on page 63) or Multi Wavelength (section 6.1.1.2 on page 64) is selected, define the unit, wavelength, absorbance formula, wavelength  $\lambda$ , concentration factor Kx, concentration resolution, chemical form and calibration equation. More information on Free Programming parameters see section 6.1.2 on page 72.

**Table 8 Program descriptions**

Program Type	Description
<b>Single Wavelength</b>	Measurements at a defined wavelength
<b>Multi Wavelength</b>	In the Multi Wavelength mode, absorbance values can be measured at up to four wavelengths and the results can be mathematically processed to obtain sums, differences and relationships.
<b>Free Programming</b>	This is an advanced form of programming for original user-developed methods. In Free Programming, the user defines the measurement process, variables and calculations involved in obtaining a reading.

### 6.1.1.1 Single wavelength settings


If the Single Wavelength mode is selected, the following parameters can be defined:

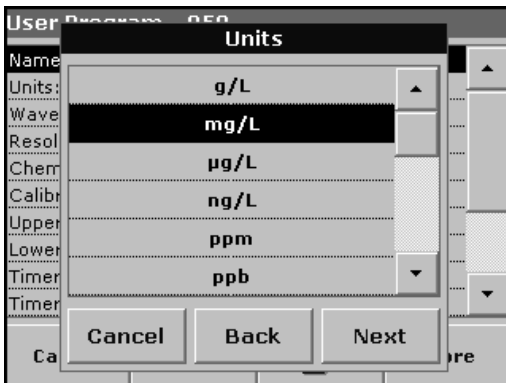

**Units:**

Select the required unit from the list and press **Next**.


*Note:* Units of measure not included in this list can be added in the edit program under **Program Options, Edit**. Select **Units, Edit** and then **New**.

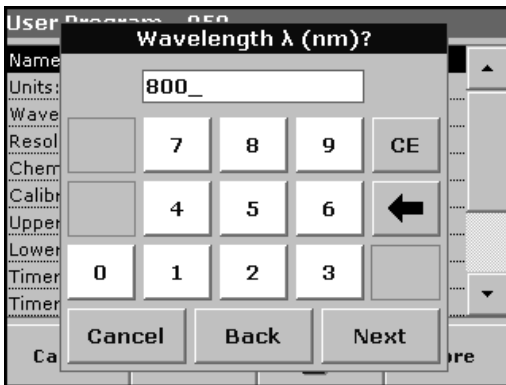

**Wavelength (single wavelength program type):**


Use the alphanumeric keypad to enter the measurement wavelength. The entered wavelength must be in the range from 190–1100 nm.

Press **Next** to proceed.

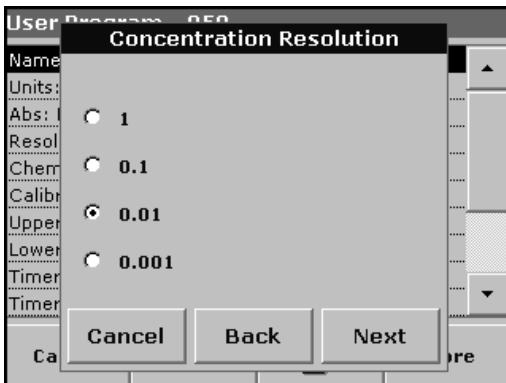

**Concentration resolution (number of decimal places)**

Select the required number of decimal places from the displayed list and press **Next**.

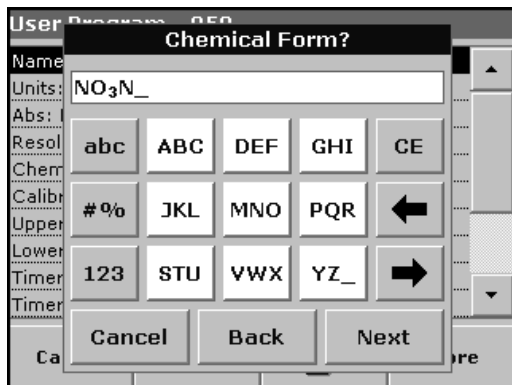


### Chemical form:

Enter the chemical formula used in the display to represent the analysis parameter.

Use the alphanumeric keypad to enter the chemical form and press **Next** to enter Calibration settings.

### 6.1.1.2 Multi wavelength settings

If the Multi Wavelength mode is selected, the following parameters can be defined:

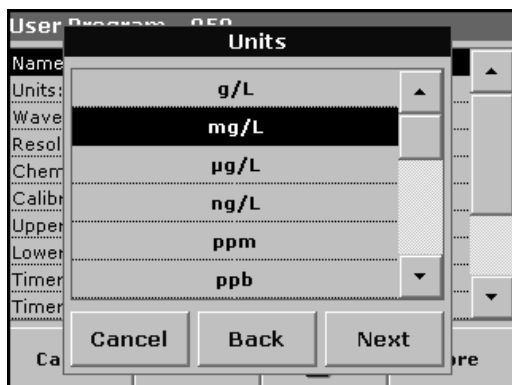

### Units:

Select the required unit from the list and press **Next**.

*Note: Units of measure not included in this list can be added in the edit program under **Program Options, Edit**. Select **Units, Edit** and then **New**.*



### Absorbance formula (multi-wavelength program type):


The Absorbance Formula menu is used to define the wavelengths and the coefficients used in the formula. The absorbance formula defines the calculation for the multi-wavelength measurement.

Press the appropriate key to edit the input.

Press the **Formula** key.

In the displayed list, select the formula for the program and press **OK** to confirm.



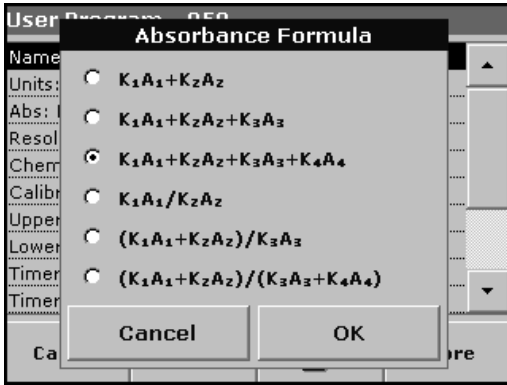

**List of available absorbance formulas**


A<sub>1</sub> is the absorbance at wavelength 1,

A<sub>2</sub> is the absorbance at wavelength 2 and so on

K<sub>1</sub> is the factor at wavelength 1,

K<sub>2</sub> is the factor at wavelength 2 and so on

If a subtraction has to be completed, the factors can be entered with a minus sign.

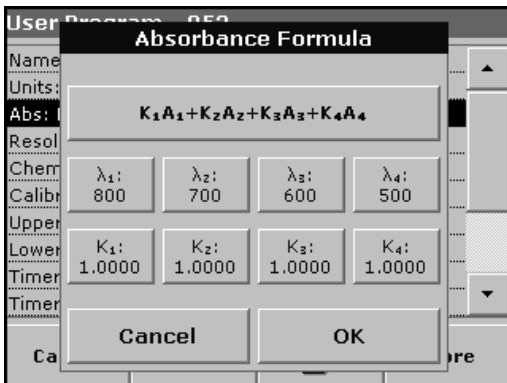

**Wavelength λ<sub>x</sub>:**


Press a λ<sub>x</sub> key and use the alphanumeric keypad to enter a wavelength. Press another λ<sub>x</sub> key and enter the next wavelength. If necessary, repeat until all the wavelengths for the formula have been entered. The wavelengths must be in the range from 190–1100 nm. Press **OK**.

**Concentration Factor K<sub>x</sub>**

Multiplication factor for converting absorbance values into concentration values.


Press a factor key and use the alphanumeric keypad to enter a factor. If the formula includes more than one factor, press another factor key and enter another factor and repeat until all the factors have been entered. Press **OK**. When all the relevant data have been entered, press **Next**.

*Note: Up to five digits can be entered, including a maximum of 4 decimal places to the right of the decimal.*

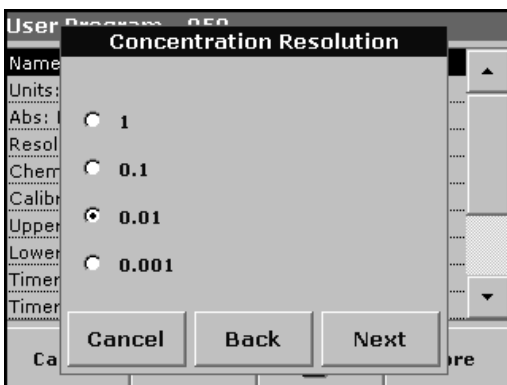

**Concentration resolution (number of decimal places)**

Select the required number of decimal places from the displayed list and press **Next**.

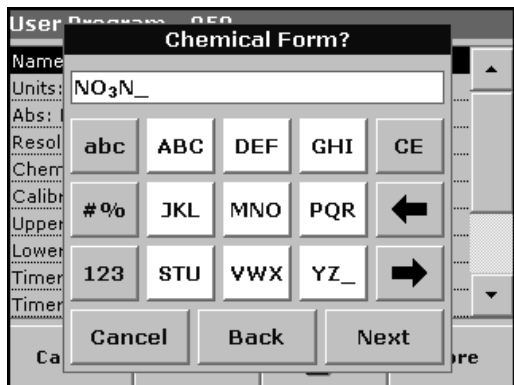

### Chemical form:

Enter the chemical formula used in the display to represent the analysis parameter.

Use the alphanumeric keypad to enter the chemical form and press **Next** to enter Calibration settings.

### 6.1.1.3 Calibration settings for single and multi wavelength mode

A method is calibrated by determining the absorbance values of several standard solutions of known concentration.


There are three ways to create and store a calibration curve. Instructions for each method follows (Table 9).

**Table 9 Calibration settings**

Mode	Descriptions
<b>Enter values</b>	A calibration table is created by entering the concentration values and the absorbance values of the analyte solution. The absorbance values are plotted versus standard concentrations and the calibration curve is displayed as a graph (page 66).
<b>Read Standards</b>	A calibration table is created by entering the concentration values of the standard solutions and then measuring the absorbance of the analyte solutions. The absorbance values are plotted versus standard concentrations and the calibration curve is displayed as a graph (page 68).
<b>Enter Formula</b>	If the calibration curve can be determined from the mathematical relationship between concentration and absorbance by linear regression, etc., the corresponding formula can be selected (linear, 2nd or 3rd order polynomial) from a list and the appropriate factors can be entered (page 69).

### Calibration by entering calibration values

Enter concentration/absorbance

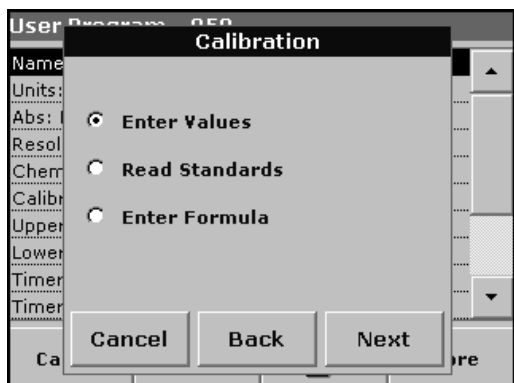

1. Select the **Enter Values** and press **Next**.

Calibration Values	
mg/L	Abs
0.0000	0.000
1.0000	0.200
2.0000	0.400
3.0000	0.600
4.0000	0.800


Buttons: Exit, mg/L, Abs, Graph

- To enter the standard concentrations and corresponding absorbance values in the displayed table, press the "+" symbol. Use the alphanumeric keypad to enter the values.

Press **OK** and enter the corresponding absorbance value. Press **OK**.


The entered data are displayed in the table. Repeat the sequence for each data point to enter it.

- To change a value in the table, highlight the appropriate line, press the unit key (e.g. **mg/L**) or Abs and enter the changed value via the alphanumeric keypad.

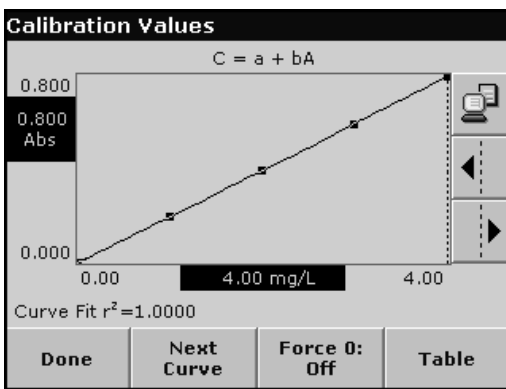

- When the data have all been entered, press **Graph**.

*Note: The correlation coefficient ( $r^2$ ) is shown on the left below the axes.*

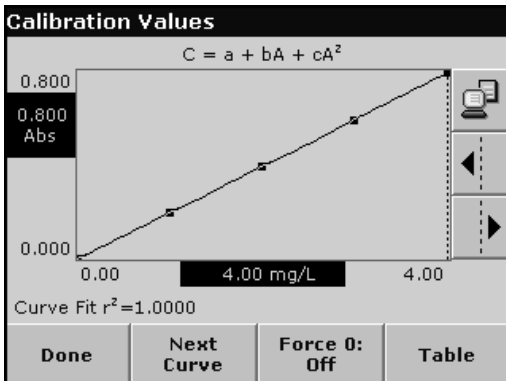

- The linear equation corresponds to the standard setting. Press **Next Curve** to display the polynomial 2nd order curve. Press **Next Curve** again, to display the polynomial 3rd order curve.


- Press **Force 0** to change the setting from **Off** to **On**. The curve now passes through the origin of the coordinate system.

*Note: This may have an adverse effect on the correlation coefficient ( $r^2$ ).*


- Press **Table** to display the table again.

- When the table has been completed and the curve type has been chosen, press **Done** when the graph is displayed or **Exit** when the table is displayed. Go to section 6.1.1.4 on page 70.

## Calibration by reading standards

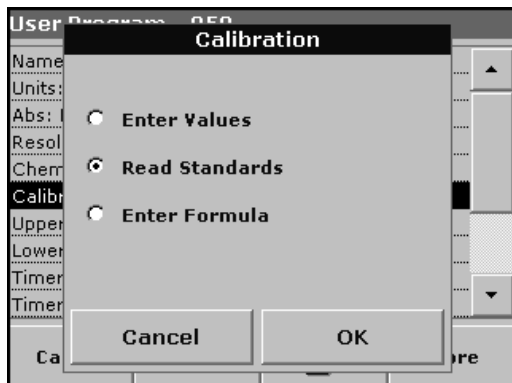

1. Press **Read Standards** and press **Next**.
2. To enter the standard concentrations in the displayed table, press the "+" symbol. Use the alphanumeric keypad to enter the standard concentration. Press **OK**.

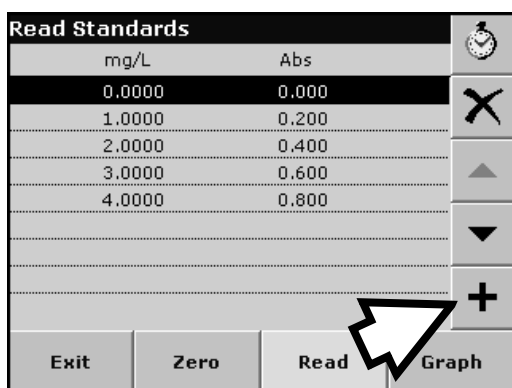

3. Press the "+" symbol again (see arrow) and enter the next standard concentration. Repeat this sequence until all standard concentrations (maximum of 24 solutions) have been entered.
4. Highlight the line with the appropriate concentration and insert the cuvette/cell with the corresponding standard solution.
5. Insert the zero solution into the cell compartment. Press **Zero**.
6. Insert the **first** standard solution into the cell compartment. Press **Read**.

Insert the **second** standard solution into the cell compartment. Press **Read**.

Repeat this sequence until all the standard solutions have been measured (maximum of 24 solutions).

The entered and measured data are displayed in the table.

**Note:** To delete a standard concentration, highlight the appropriate line and press the **Delete** icon.

The timer icon shown in the display helps to ensure, when necessary, that the steps of an analysis are correctly timed (e.g. reaction times, wait times, etc., can be exactly specified). When the specified time has elapsed, an acoustic signal is emitted. The use of the timer has no influence on the measurement program.


7. When the data have all been entered and the measurements have all been completed, press **Graph**.
8. The linear curve corresponds to the standard setting. Press **Next Curve** to display the polynomial 2nd order curve. Press **Next Curve** again to display the polynomial 3rd order curve.

9. Press **Force 0** to change the setting from **Off** to **On**. The curve then passes through the origin of the coordinate system.


*Note: This may have an adverse effect on the correlation coefficient ( $r^2$ ).*

10. Press **Table** to display the table again.
11. When the table has been completed and the curve type has been chosen, press **Done** when the graph is displayed or **Exit** when the table is displayed. Go to section 6.1.1.4 on page 70.

### Calibration by entering the formula

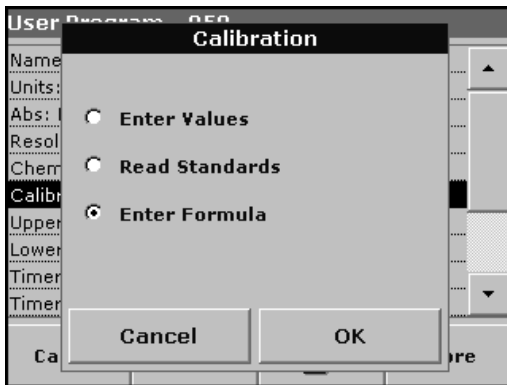

1. Press **Enter Formula** and press **Next**.

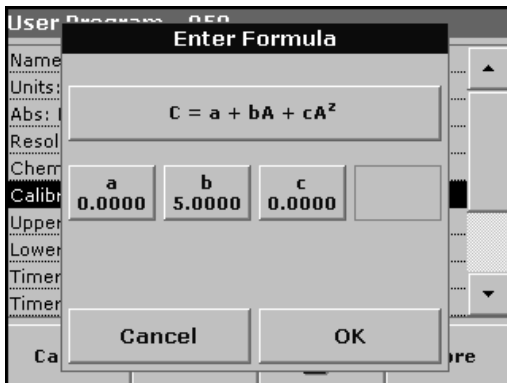

2. Press the formula key.


A list of available formulas (linear and 2nd and 3rd order polynomial) is displayed. Up to 4 coefficients can be entered, depending on the selected formula. Press the required formula.

3. Depending on the selected formula, the required coefficients (a, b, c...) are displayed. Press the coefficient keys and enter the corresponding values via the alphanumeric keypad. After each entry, press **OK** to confirm.

*Note: The coefficients can have 5 digits and can have a positive or a negative sign.*

## Advanced Operations

### 6.1.1.4 Store a user program

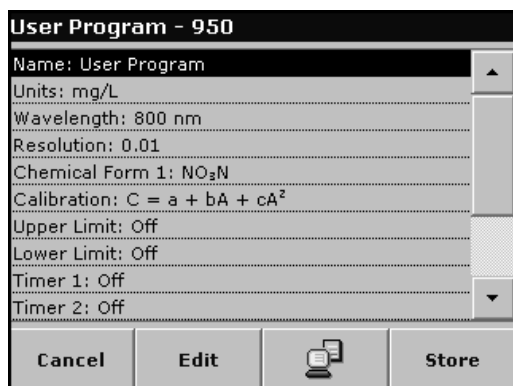

The input of the basic data is complete. An overview of the variable program data is displayed.

4. To enter more specifications or change existing ones, highlight the appropriate line and press **Edit**.
5. Select **Store** to save the user program.
6. Press the **PC & Printer** icon to send the program data to a printer or to a USB memory stick (connect the USB memory stick to the USB interface first).

*Note: The test data will be formatted in the preinstalled PrgData folder as a .csv and as a .lst file.*

7. To transfer the program data from one instrument to another create a new folder on the USB memory stick named **dbhl** and copy the **.lst** files into this folder. (see section 6.8.2 on page 106).

### 6.1.1.5 Additional user-defined parameters and functions

In addition to the previously defined basic data, additional parameters and functions can be defined for user programs:

- measuring range upper and lower limits
- timer functions
- chemical forms



#### Upper and lower limits of the measuring range


It is possible to enter a maximum (upper) and minimum (lower) measurement value. An error message is displayed if a reading is above the upper limit or below the lower limit.

1. Highlight the appropriate line in the overview of the program data and press **Edit**.
2. Set **On** and press the **0.000** key to enter the measuring range limit. Confirm the entry by pressing **OK**.

**Timer 1 / Timer 2 / Timer 3 / Timer 4:**

This function can define time intervals for up to four timers. Timer designations such as Shake, Wait and Swirl can be assigned.

1. Highlight the appropriate line in the overview of the program data and press **Edit**.
2. The timers are activated or deactivated with the check boxes in the left part of the display. In the next column, a selection can be made from a list of names that designate the corresponding work step. In the third column, the times for each timer are entered (in mm:ss).

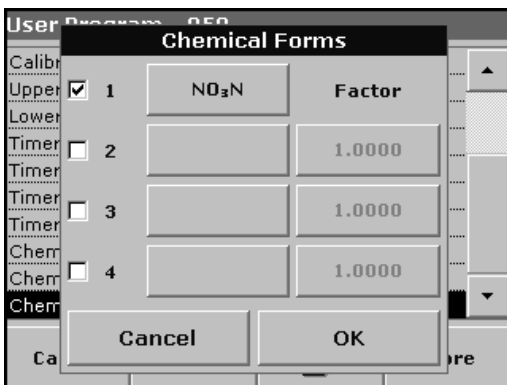**Chemical form 2 / chemical form 3 / chemical form 4:**


If a **Chemical Form 1** has been defined, up to three additional alternative forms can be entered here.

1. Highlight the appropriate line in the overview of the program data and press **Edit**.
2. The chemical forms are activated or deactivated with the check boxes in the left part of the display.
3. Press the left key to enter another chemical form with the alphanumeric keypad and press **OK** to confirm. Press the right key to enter the conversion factor to calculate the concentration of the additional chemical form from the concentration of Chemical Form 1 and press **OK** to confirm.
4. Press **Store** to save the program data. Press **Cancel** to return to the Main Menu.

## 6.1.2 Free programming program type

Free Programming is an advanced option for entering original user-developed methods. When the Free Programming option is selected, an overview of the specifications of the programmed test is displayed. Each input option can be modified to develop the user method. Refer to Table 10 for more information. To modify an input option, select the appropriate line and press **Edit**.

**Important Note:** Perform steps 1–6 of Programming a New User Method (section 6.1.1 on page 61) before proceeding with Free Programming.

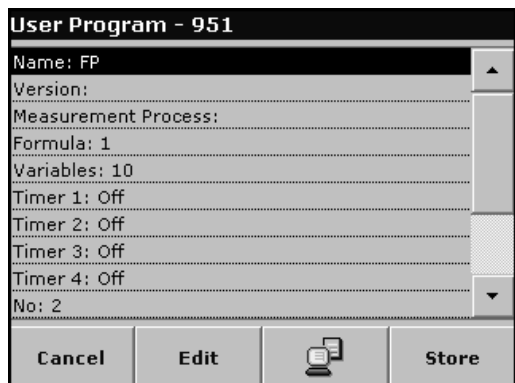

Highlight the line containing the program point that is to be edited or defined and press **Edit**.

**Table 10 Definitions of the program points**

Program point	Description
<b>Name</b>	Name of the analysis parameter
<b>Version</b>	An abbreviation or version number assigned by the user is entered here.
<b>Measurement Process</b>	Exact definition of the test: the number of wavelengths at which measurements are made, the number of absorbance measurements needed, the keys to be used, any waiting periods between measurements, etc.
<b>Formula</b>	Definition of the formulas with which the test result is calculated.
<b>Variables</b>	The number of variables shown in the display depends on the definition of the measurement process and the formulas. Input of the numerical values of the wavelengths, factors, constants, etc.
<b>Timer 1, Timer 2, Timer 3, Timer 4</b>	Used to enter abbreviations and defined times for up to four timers. Highlight the appropriate line and press <b>Edit</b> . The timers are activated or deactivated with the control boxes on the left of the display. In the next column, a selection can be made from a list of names that describe the corresponding work step. In the third column, the times for each active timer are entered.

### 6.1.2.1 Measurement process

The measurement process defines the handling and the measurements of the test:

- At which and how many wavelengths should measurements be completed
- How many absorbance measurements must be completed?
- When should the zero measurement and the sample measurement be completed?
- Are waiting times necessary between measurements?



- Should individual program sequences be repeated?

The elements of a measuring sequence, such as zero and sample measurements and the timer(s) (reaction times, waiting times, etc.) are individually defined.


### 6.1.2.2 Enter a new element of a measuring sequence

**Important Note:** Each component of the measurement process **must** be entered in the order in which it will be completed.



1. Highlight the **Measurement Process** line in the data overview and press **Edit**.
2. Press **Edit** again and then **New**.

#### Content and definition of the keys

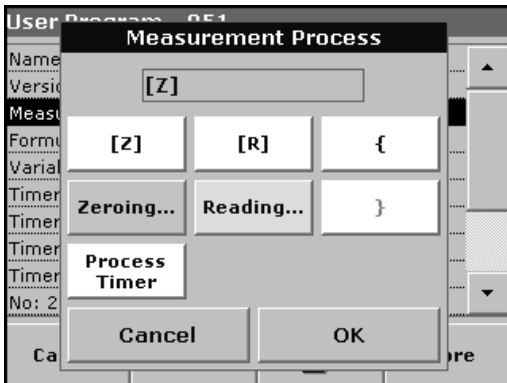

#### [Z] key / Zeroing

1. Press the **[Z]** key to program a zero measurement. Confirm with **OK**.
2. Press **New** and then **Zeroing...** and use the alphanumeric keypad to enter the wavelength at which the zero measurement is to be completed. Press **OK** and confirm the input by pressing **OK** again.
3. If zero measurements are to be carried out at a number of wavelengths, repeat the above two steps for each wavelength.

**Note:** The entered measurement sequence is displayed.




### Process Timer key

1. Press the **Process Timer** key to enter any waiting, reaction or handling times that have to be taken into account. Use the alphanumeric keypad to enter the time. Press **OK** and confirm the input by pressing **OK** again.

*Note: This time is integrated into the measurement process.*


*Note: The entered measurement sequence is displayed.*

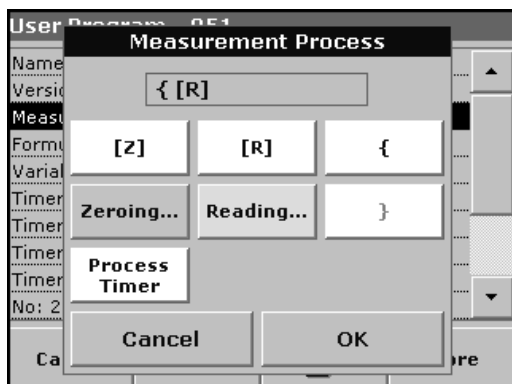

### [R] key / Reading...

1. Press the **[R]** key to program a measurement of the substance that is to be analyzed. Confirm with **OK**.
2. Press **New** and then **Reading...** and use the alphanumeric keypad to enter the wavelength at which the measurement is to be completed. Press **OK** and confirm the input by pressing **OK** again.
3. If measurements are to be completed at a number of wavelengths, repeat the above two steps for each wavelength.

*Note: The entered measurement sequence is displayed.*

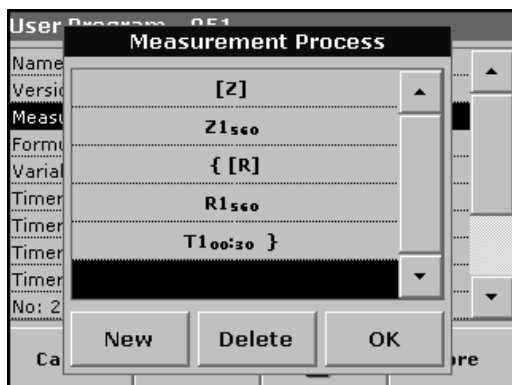

### { } key


Elements of the measurement sequence that are to be repeated are placed in brackets.

The left bracket "{" marks the start of the sequence that is to be repeated and the right bracket "}" marks the end.

*Note: The key showing the right bracket remains inactive until a left bracket is entered.*

1. Press {.
2. Press the key that defines the sequence that is to be repeated: **[Z]** or **[R]**. Confirm with **OK**.
3. Press **New** and then press **Zeroing...** or **Reading...** and use the alphanumeric keypad to enter the wavelength at which the measurement should be completed. Press **OK** and confirm the input by pressing **OK** again.
4. Press } to end the sequence.

*Note: If an action such as a zero measurement recurs at different stages of a measurement sequence, the series of actions is numbered sequentially (e.g. Z1, Z2, etc.).*

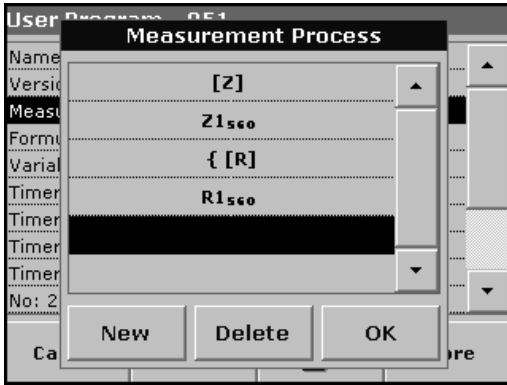

**Delete an element of a measuring sequence**

Select the appropriate line and press **Delete**. The element is deleted.

**Insert an element of a measuring sequence**


Select the line in the measuring sequence where the insertion is to be made and press **New**.

A new element can be entered at the selected position.

When the input is complete, press **OK** in the "Measurement Process" display. The data overview is then displayed.


**6.1.2.3 Enter the calibration formula (evaluation formula)**

The calibration formula (evaluation formula) defines the calculation and display of intermediate and final results. The previously defined elements of the measuring sequence are the basis for calculating the concentrations.

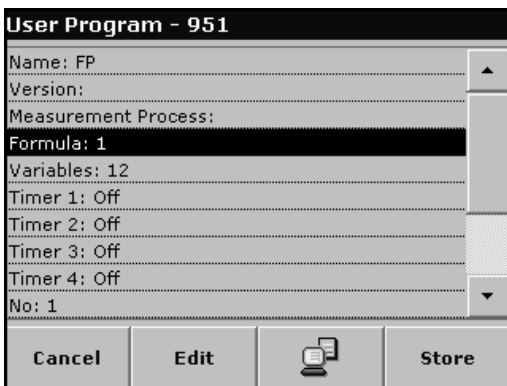


**Enter calibration formula C1**

1. Highlight the **Formula** line in the data overview and press **Edit**.

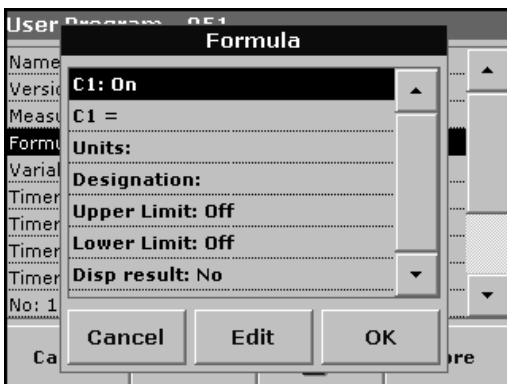

2. Highlight the line **C1: Off** and press **Edit**.
3. Select **C1: Off** again and press **Edit**. The display changes to **C1: On**.
4. Highlight the next line **C1 =** to define the formula and press **Edit**.

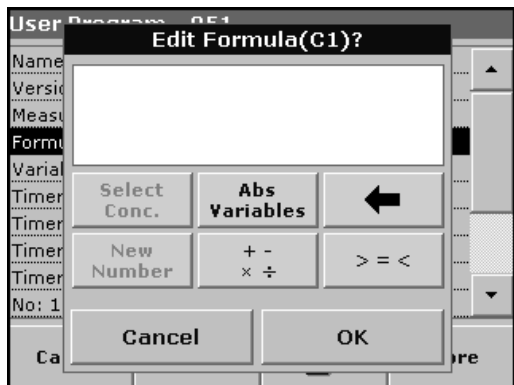

Refer to Table 11 for detailed information on the Edit formula keys.


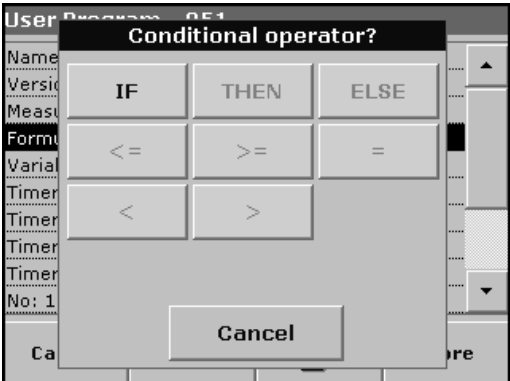
**Note:** The evaluation formula is built up successively in the display in accordance with the input.


**Note:** The **arrow** key deletes the most recently entered element of the formula.

**Table 11 Edit formula key descriptions**

Screen	Key	Description
	<p><b>Abs/Variables</b></p>	<p>Press the <b>Abs/Variables</b> key to select, from the displayed list, the required element of the defined measuring sequence and therefore the corresponding measurement wavelength, so that this can be taken into account in the formula.</p>
	<p><b>New Number</b></p>	<p>Press <b>New Number</b> to enter a new factor or constant.</p>

Table 11 Edit formula key descriptions (continued)


Screen	Key	Description
	<p><b>+ - ÷ x</b></p>	<p>Press <b>+ - ÷ x</b>, to enter a mathematical operation. Select the operation and press <b>OK</b> to confirm. The available choice of mathematical operations depends on the defined formula. This means that functions such as "( )" or "ln"/"log" etc. are only active if a term in parentheses or the calculation of a logarithm is mathematically permissible in the defined formula (this also applies to the basic mathematical operations). The following mathematical operations are available:</p> <ul style="list-style-type: none"> <li>• + (Addition)</li> <li>• - (Subtraction)</li> <li>• ÷ (Division)</li> <li>• x (Multiplication)</li> <li>• ^ (Exponent)</li> <li>• Ln (Natural logarithm)</li> <li>• Log (Common logarithm)</li> </ul>
	<p><b>&gt;=&lt;</b></p>	<p>Press <b>&gt;=&lt;</b> to include logic statements/links/conditions in the formula. The following functions are available: = (Equal to), &lt; (Less than), &gt; (Greater than), &lt;= (Less than or equal to), &gt;= (Greater than or equal to), IF, THEN, ELSE</p> <p>When the evaluation formula C1 has been entered completely, press <b>OK</b> to confirm. Press <b>OK</b> again to return to the Formula display.</p> <p>When the formula C1 has been entered and confirmed, the parameter name, upper and lower limits of the measuring range and display result (yes, no) can be entered.</p>

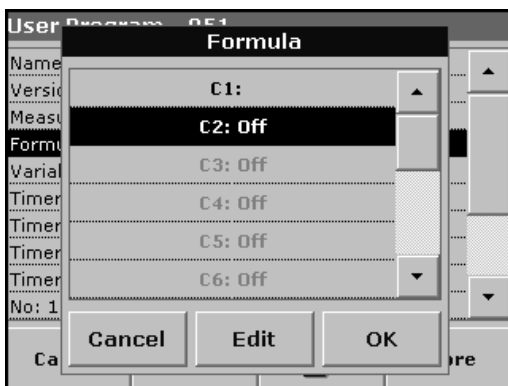

**Enter the next calibration formula (C2 or Cn)**

5. Highlight **C2: Off** and press **Edit**.
6. Select **C2: Off** again and press **Edit**. The display switches to **C2: On**.
7. Highlight the next line **C2 =** to define the formula and press **Edit**.

In addition to the keys described in Table 11, only one other function is available:

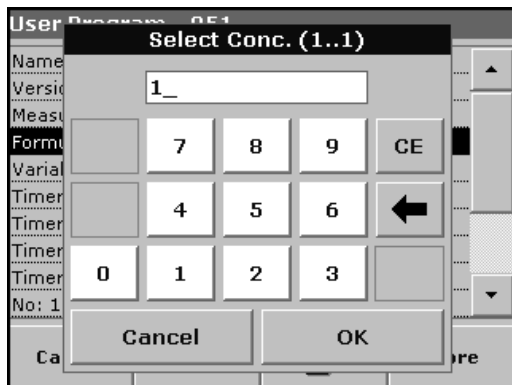

### Select Conc. key

If an already defined formula, in this case C1, is to be taken into account in the formula for C2, press **Select conc.**


Enter the number of the formula (e.g. 1 for C1) and press **OK** to confirm.

Cn can now be linked with a mathematical operation.

**Note:** The Cn concentrations that are to be calculated are numbered in sequence: C1, C2, C3, etc.

**Note:** When the first Cn formula has been defined, the Formula list is automatically extended by Cn+1.

### 6.1.2.4 Enter variables

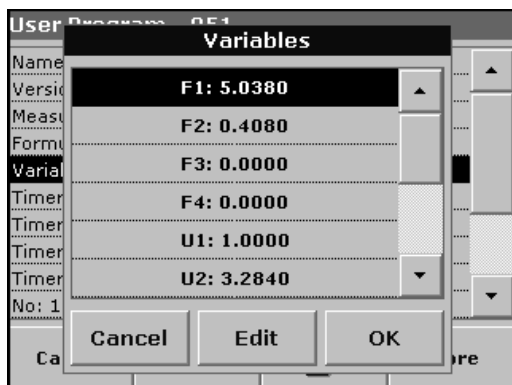

1. Highlight the **Variables** line in the data overview and press **Edit**.
2. Select the variable to be edited, press **Edit** and use the alphanumeric keypad to enter the data specified in the working procedure (for F1, F2, I1, U1 etc.). Press **OK** to confirm each entry.

Abbreviation of Variables:

F1: Factor 1

F2: Factor 2

$\lambda$ 1: Wavelength 1

U1: Conversion Factor 1 for the first chemical form

U2: Conversion Factor 2 for the further chemical form etc.


### 6.1.2.5 Save a free programming user program

1. Press **Store** to save the entered data. The data can be stored under any data point (Measurement sequence, Formula, Timer, etc.).
2. Press the **PC & Printer** icon to send the program data to a printer or to a USB memory stick (connect the USB memory stick to the USB interface first).

**Note:** The test data will be formatted in the preinstalled PrgData folder as a **.csv** and as a **.lst** file.

3. To transfer the program data from one instrument to another create a new folder on the USB memory stick named **dbhl** and copy the **.lst** files into this folder (see section 6.8.2 on page 106).

### 6.1.3 Select a user program

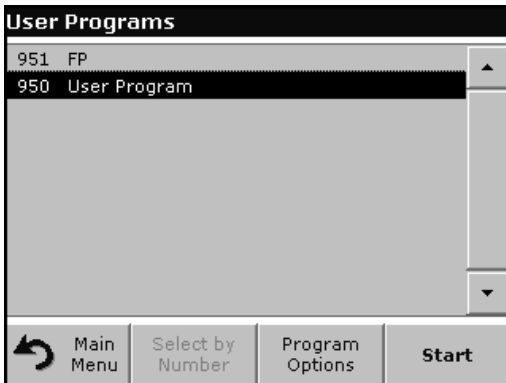

1. Press **User Programs** in the Main Menu to view an alphabetical list of user programs with program numbers.

The User Programs list will appear.


**Note:** Use the scroll bar to scroll through the list quickly.

**Note:** Highlight the selection by pressing it or press **Select by Number** to search for the program by number.

2. Press **Start** to run the program.

### 6.1.4 Add, edit and delete user programs from the favorites list

The most frequently used tests/methods in the User Program menu can also be added to the list of favorites to simplify their selection.

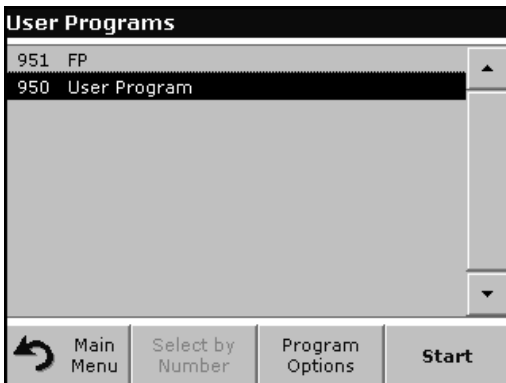

1. Press **User Programs** in the Main Menu.

The User Programs list will appear.

2. Highlight the selection by pressing it or press **Select by Number** to search for the program by number.

**Note:** Use the scroll bar to scroll through the list quickly. Use the alphanumeric keypad to enter the test number (program number) and press OK to confirm.



3. Press **Program Options**.

4. Press **Add to Favorites**, **Edit** or **Delete** and press **OK** to confirm.

**Note:** If the stored program is deleted in User Programs, it will also be deleted in Favorites Programs.

## Advanced Operations


### 6.1.4.1 Add to Favorites



1. Press **Add to Favorites** and press **OK** to confirm.

The program is added to the Favorites.

### 6.1.4.2 Edit

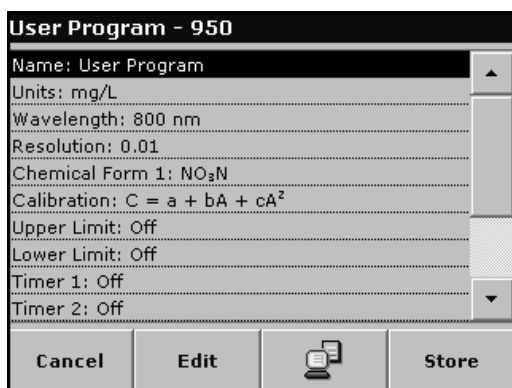

1. Press **Edit** and press **OK** to confirm.

An overview of the specifications of the programmed test is displayed. More information about the input options is provided in the section 6.1.2 on page 72.

### 6.1.4.3 Delete



1. Press **Delete** and press **OK** to confirm.

The program is deleted from the list of User Programs.

**Note:** If the stored program is deleted in User Programs, it will also be deleted in Favorites Programs.




## 6.2 Favorite Programs

The most frequently used tests/methods in the **Stored Programs** menu and the **User Programs** menu can also be added to the list of favorites to simplify their selection.

To add **Stored Programs** and/or **User Programs** to the favorites list or the favorite programs, see section 6.1.4 on page 79.

### 6.2.1 Recall a favorite program

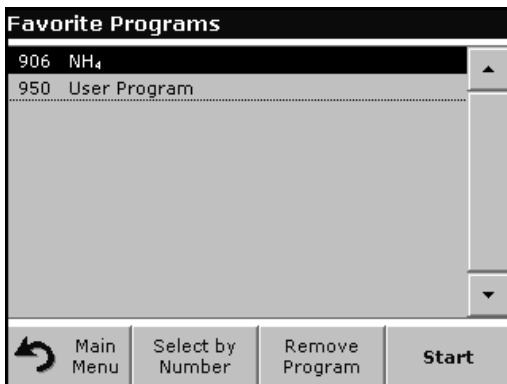

1. Press **Favorite Programs** in the Main Menu.

The Favorite Programs list will appear.

2. Highlight the selection by pressing it.

*Note:* Use the scroll bar to scroll through the list quickly.

*Note:* Press **Select by Number** to search for the program by number. Use the alphanumeric keypad to enter the test number (program number) and confirm the input by pressing **OK**.

3. Press **Start**.

### 6.2.2 Delete a favorite program



1. Press **Favorite Programs** in the Main Menu.

The Favorite Programs list will appear.

2. Highlight the selection by pressing it.

*Note:* Use the scroll bar to scroll through the list quickly.

*Note:* Press **Select by Number** to search for the program by number. Use the alphanumeric keypad to enter the test number (program number) and confirm the input by pressing **OK**.

3. Press **Remove Program** and press **OK** to confirm.

*Note:* If a **Favorite Programs** is deleted, it will stay in the **User Programs** or **Stored Programs**.

*Note:* If the stored program is deleted in **User Programs**, it will also be deleted in **Favorites Programs**.

### 6.3 Standard Addition – monitoring/checking results

The accuracy of measured values (their correspondence with the actual concentration of the analyte in the sample) and their precision (correspondence of the measurement results obtained from several samples containing the same concentration of the test analyte) can be determined or improved using the standard addition method.

This method (also referred to as spiking) serves to identify sample-specific interference factors, e.g. substances in the sample that falsify the analysis (sample matrix effect), a defective measuring instrument or contaminated reagents.

**Method:**

A defined amount (concentration) of a standard solution of the test substance is added to the sample. The detection rate should be close to 100%.

$$\text{Detection rate} = \frac{\text{Measured value after a standard addition}}{\text{Expected value after a standard addition}}$$

Detection rate	Conclusion	
100%	Probability that the measurement results are correct is high.	
< 100%	Assumption: The analysis was falsified by substances in the sample (sample matrix effect) <b>Test to determine whether a sample matrix effect is present:</b> Use distilled water instead of the sample. Add standard solution as described in the procedure.	
	Detection rate	Conclusion
	100%	Ions in the sample are interfering with the analysis, causing false results to be obtained.
≠ 100%	No interfering ions - consider other interference factors.	

**Measures to identify other interference factors:**

**Checklist:**

1. Check if the procedure is completed correctly:
  - a. Are the reagents added in the correct order?
  - b. Is enough time allowed for color development?
  - c. Is the correct glassware in use?
  - d. Is the glassware clean?

- e. Does the test require the sample to be at a certain temperature?
  - f. Was the pH of the sample in the correct range?
  - g. Is the pipette volume correct?
2. Check the used reagents by repeating the standard addition procedure with freshly prepared reagents.

Detection rate	Conclusion	
100%	The originally used reagents were defective. <b>Check the standard solution:</b> Repeat the standard addition procedure with a freshly prepared standard solution.	
	Detection rate	Conclusion
	100%	The originally used standard solution was defective.

If none of these measures resolves the problem, please contact the manufacturer or a sales representative.

### 6.3.1 Complete a standard addition

Carry out the standard addition in accordance with the corresponding procedure.

There are two different methods:


**Peak volume (Standard addition):**

Defined volumes of a standard solution are added step by step to an already analyzed sample. The sample is measured again after each addition.

**Sample volume:**


A defined volume of a standard solution of known concentration is added to the volume of sample specified in the procedure and the sample is measured after each addition. In most cases three different standard solutions are prepared and the procedure is repeated for each of them.

**Note:** The units and chemical forms used for the sample are used for the standard solutions. Be careful to ensure that to use the correct units for subsequent entries.

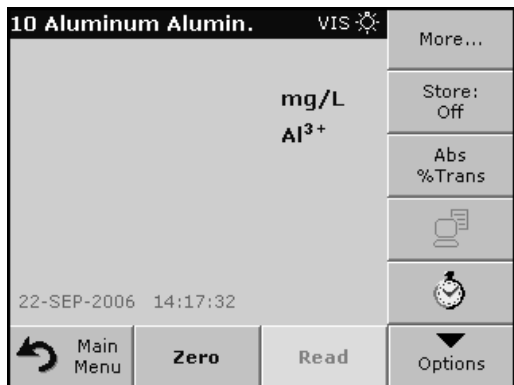


## Peak volume/sample volume methods

1. Select **Stored Programs** in the main menu. Select the required program.
2. Press **Start**.
3. Analyze a sample without added standard solution in accordance with the instructions in the Procedures Manual. When the measurement is complete, leave the sample cuvette/cell in the cuvette/sample cell holder.

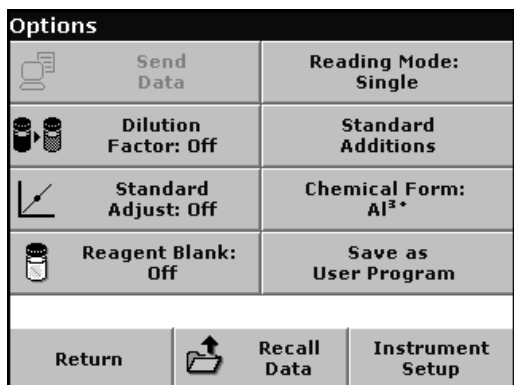


4. Press **Options>More>Standard Addition**.

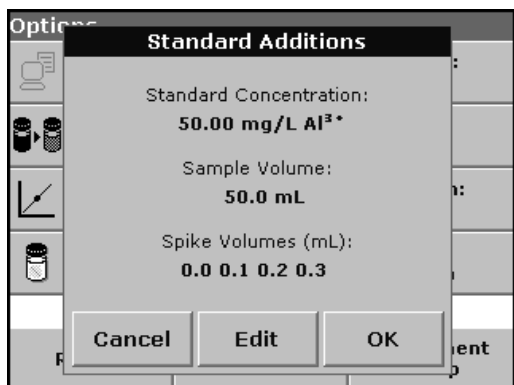


An overview of the data of the standard addition procedure is displayed.

5. Press **OK** to accept the standard values for standard concentration, sample volume (total) and standard addition volume. Press **Edit** to change any of these values.

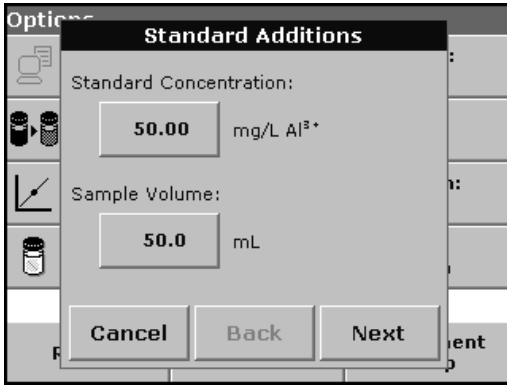

6. Press the key of the value to change it. Use the alphanumeric keypad to change the value. Press **OK** to confirm.

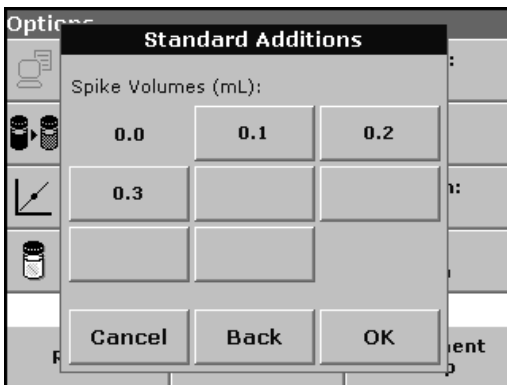

7. Press the keys to enter the standard addition volumes. Use the alphanumeric keypad to enter the new data and press **OK**.

Standard Additions		
mL	mg/L	%
Standard	Al <sup>3+</sup>	Recovery
0.0	0.188	100
0.1	0.265	92
0.2	0.345	89.3
0.3	0.431	88.8
Exit	Zero	Read
		Graph


**Description of the table of readings**

- The **first** column shows the standard addition volume. 0 mL stands for a sample containing no added standard.
- The **second** column shows the reading of the samples with and without added standard.
- The **third** column shows the detection rate of the added standard.

*Note: The black highlighted line is active.*

8. The reading of the sample in the instrument, without added standard, is automatically shown under 0 mL.
9. Prepare the standard addition solution as described in the procedure.
10. Use the arrow keys to select the first standard addition volume in the table and insert the cuvette/cell with the corresponding volume of added standard into the cell compartment. Press **Read**.

Repeat the procedure from point 8 with all the other standard addition solutions.

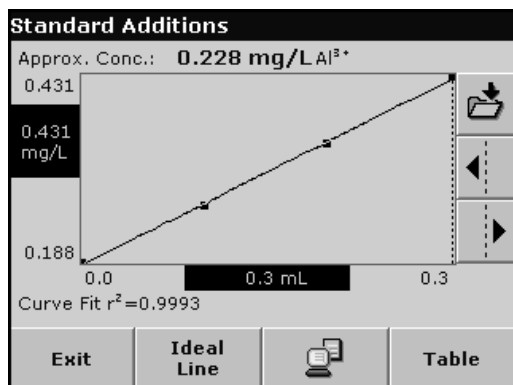

11. After all the standard addition solutions have been measured, press **Graph**.

The regression line through the standard addition data points is displayed.

The correlation coefficient  $r^2$  indicates how close the data points are to the line.

If the correlation coefficient = 1, the curve is linear.

The concentration shown above the curve is the estimated concentration of the sample without the added standard.

**Note:** In the curve menu, the name on the **Curve** key switches to **Table**. Press **Table** to display all the data in the table again.

12. Press **Ideal line** to display the relationship between the added standard solutions and the ideal line (detection rate 100%).

## 6.4 Single Wavelength (absorbance, concentration and transmittance measurements)

The Single Wavelength mode can be used in three ways. For sample measurements at a single wavelength, the instrument can be programmed to measure the absorbance, % transmittance or concentration of the analyte.

Absorbance measures the amount of light absorbed by the sample, in units of absorbance.

% transmittance measures the percent of the original light that passes through the sample and reaches the detector.

Turn the concentration factor on to select a specific multiplier for converting absorbance readings to concentration. In a graph of concentration versus the absorbance, the concentration factor is the slope of the line.

### 6.4.1 Set up single wavelength mode

Press **Single Wavelength** in the Main Menu. Press **Options** for Parameter Setup.



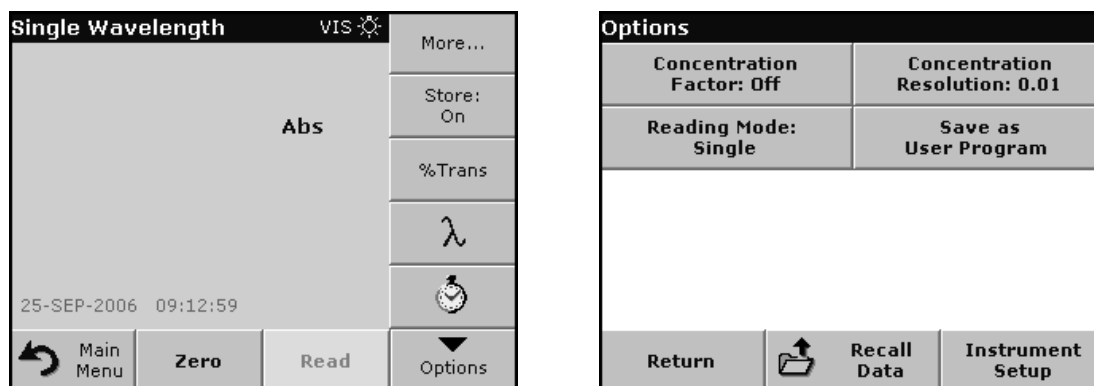



Table 12 Single wavelength setup options

Options	Description
<b>More</b>	For further Options
<b>Store Off/On</b>	With the <b>Store On</b> setting, all measurement data are stored automatically. With the <b>Store Off</b> setting, no measurement data are stored.
<b>% Trans/Abs</b>	To switch to % transmittance, concentration or absorbance readings
<b>λ Wavelength</b>	To enter the measurement wavelength. Use the alphanumeric keypad to enter the measurement wavelength. The entered wavelength must be in the range from 190–1100 nm.
<b>Timer icon</b>	This functions as a stopwatch. It helps to ensure that the steps of an analysis are correctly timed (e.g. reaction times, wait times, etc., can be exactly specified). When the specified time has elapsed, an acoustic signal is emitted. The use of the timer has no influence on the measurement program.
<b>Concentration Factor</b>	Multiplication factor for converting absorbance values into concentration values.
<b>Concentration Resolution</b>	To select the position of the decimal point in the calculated concentration readings.

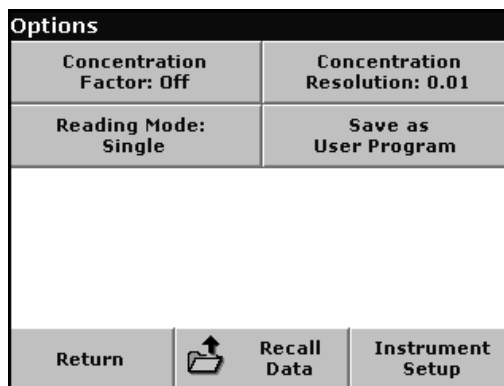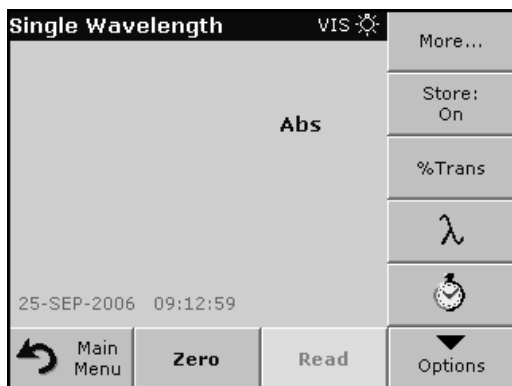


**Table 12 Single wavelength setup options (continued)**

Options	Description
<b>Reading Mode</b>	<b>Single Reading Mode:</b> A reading is only displayed after a measurement has been carried out (press Read; standard setting) (see section 6.4.2 on page 89). <b>Continuous Reading Mode:</b> After the zero measurement, all readings are displayed automatically and continuously (see section 6.4.3 on page 89).
<b>Save as User Program</b>	To store the selected parameters as a User Program, see section 6.1 on page 61
<b>Recall Data</b>	Call up saved measurement data, wavelength scans or time courses, see section 5.3 on page 35.
<b>Instrument Setup</b>	Basic data of the instrument, see section 5.2 on page 23.

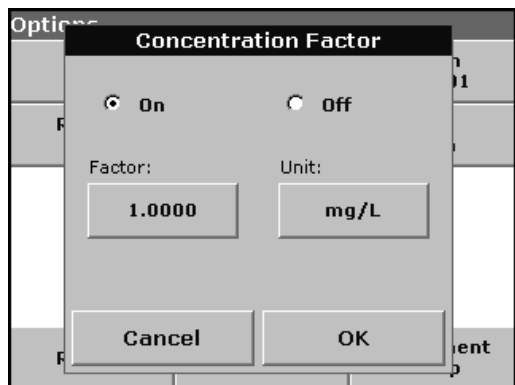

### Concentration factor:


1. Press **Concentration Factor: Off** in the Options menu. Press **On** to highlight this feature.
2. Press the "Factor" key and use the alphanumeric keypad to enter the factor by which absorbance readings are to be multiplied. Press the "Unit" key to select the units for concentration measurements or to create a new unit.
3. Press **OK** to confirm.

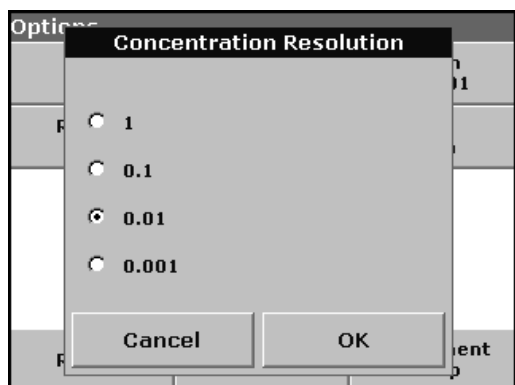

### Concentration resolution:

1. Press **Concentration Resolution** in the Options menu.
2. Select the resolution and press **OK** to confirm.




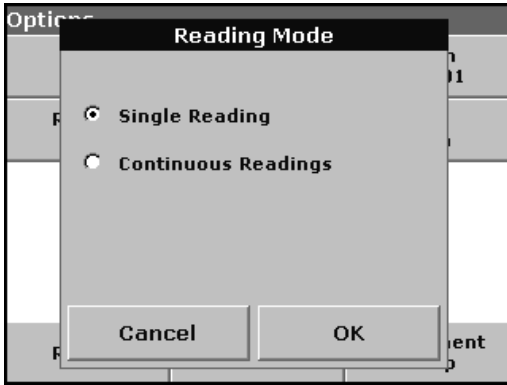

**Reading mode:**

1. To highlight the required mode, start by pressing **Reading mode**.
2. Select the required mode, then press **OK**, then **Return** to return to the result display.

**6.4.2 Take single wavelength measurements (single reading)**

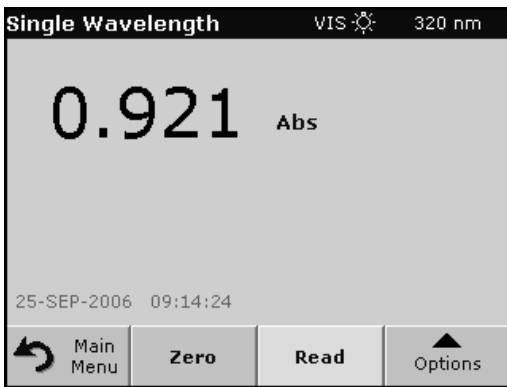

1. Insert the blank cuvette/cell into the cuvette/sample cell holder. Press **Zero**.  
*Note: The **Read** key is only active after the zero measurement has been completed.*
2. Insert the sample cuvette/cell into the cuvette/sample cell holder. Press **Read**.
3. For data storage, see section 5.3.2 on page 38.

**6.4.3 Take single wavelength measurements (continuous readings)**



1. Insert the blank cuvette/cell into the cuvette/sample cell holder. Press **Zero**.  
*Note: In the reading mode "Continuous" only the Zero key is shown to start the reading. The reading sequence is started automatically.*
2. Insert the sample cuvette/cell into the cuvette/sample cell holder.
3. Press **Options** and then the **Store** icon to store the displayed data in the Data Log.

*Note: For data storage, see section 5.3.2 on page 38.*

## 6.5 Multi-Wavelength mode – measurements at more than one wavelength

In the multi-wavelength mode, absorbance values can be measured at up to four wavelengths and the results can be mathematically processed to obtain sums, differences and relationships.

Absorbance measures the amount of light absorbed by the sample, in units of absorbance.

% Transmittance measures the percent of the original light that passes through the sample and reaches the detector.

Turning on the concentration factor allows selection of a specific multiplier for converting absorbance readings to concentration. In a graph of concentration versus the absorbance, the concentration factor is the slope of the line. Concentration is calculated using a single factor for each wavelength, which is input by the user.

### 6.5.1 Set the reading mode at different wavelengths

Press **Multi Wavelength** in the Main Menu. Press **Options** for Parameter Setup.

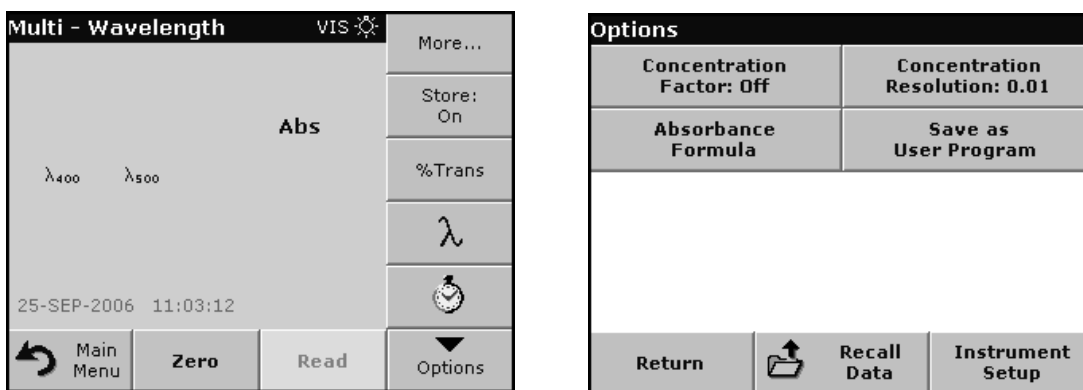


Table 13 Multi-wavelength setup options

Options	Description
More	For further Options
Store Off/On	With the <b>Store On</b> setting, all measurement data are stored automatically. With the <b>Store Off</b> setting, no measurement data are stored.
% Trans/Abs	To switch to % transmittance, concentration or absorbance readings
λ Wavelength	To enter the measurement wavelength. Use the alphanumeric keypad to enter the measurement wavelength. The entered wavelength must be in the range from 190–1100 nm.
Timer icon	This functions as a stopwatch. It helps to ensure that the steps of an analysis are correctly timed (e.g. reaction times, wait times, etc., can be exactly specified). When the specified time has elapsed, an acoustic signal is emitted. The use of the timer has no influence on the measurement program.
Concentration Factor	Multiplication factor for converting absorbance values into concentration values.
Concentration Resolution	To select the position of the decimal point in the calculated concentration readings.
Absorbance Formula	Calculation basis for evaluating samples
Save as User Program	To store the selected parameters as a User Program, see section 6.1 on page 61


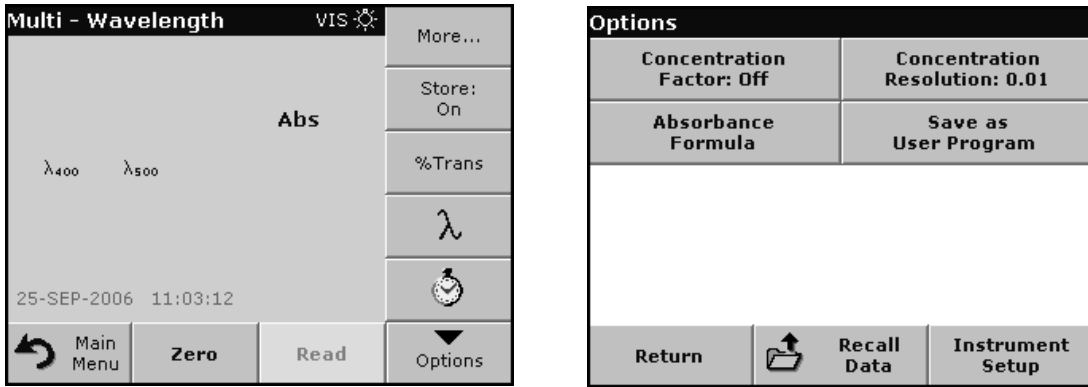


Table 13 Multi-wavelength setup options (continued)

Options	Description
Recall Data	Call up saved measurement data, wavelength scans or time courses, see section 5.3 on page 35.
Instrument Setup	Basic data of the instrument, see section 5.2 on page 23.

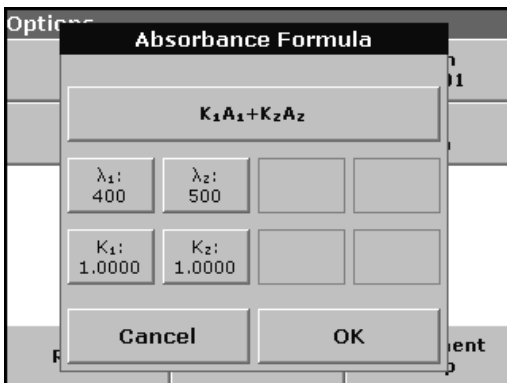

$\lambda$  / Absorbance formula:

1. Press **Absorbance Formula**.
2. The formula selected in the top key determines the number of wavelength and coefficient keys that will appear below. To change the absorbance formula, press the top key, select a formula from the displayed list and press **OK**. When a new formula is selected, the number of variables below changes to match.

The following formulas are available:

$$K_1 A_1 + K_2 A_2$$

$$K_1 A_1 + K_2 A_2 + K_3 A_3$$

$$K_1 A_1 + K_2 A_2 + K_3 A_3 + K_4 A_4$$

$$K_1 A_1 / K_2 A_2$$

$$(K_1 A_1 + K_2 A_2) / K_3 A_3$$

$$(K_1 A_1 + K_2 A_2) / (K_3 A_3 + K_4 A_4)$$


$A_1$  refers to the absorbance at wavelength 1

$A_2$  refers to the absorbance at wavelength 2, etc.

$K_1$  refers to the coefficient at wavelength 1


$K_2$  refers to the coefficient at wavelength 2, etc.

Coefficients can be set negative where subtraction is required.

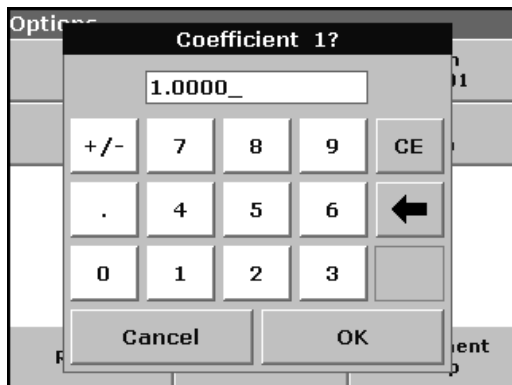


3. To change a wavelength, press one of the "λx:" keys. Enter the desired wavelength coefficient into the numeric keypad. Press **OK** to confirm.
4. To change a coefficient, press one of the "Kx:" keys. Enter the desired coefficient into the numeric keypad. Press **OK** to confirm.

**Note:** The instrument allows entry of up to 5 significant digits, with a maximum of 4 significant digits after the decimal point.

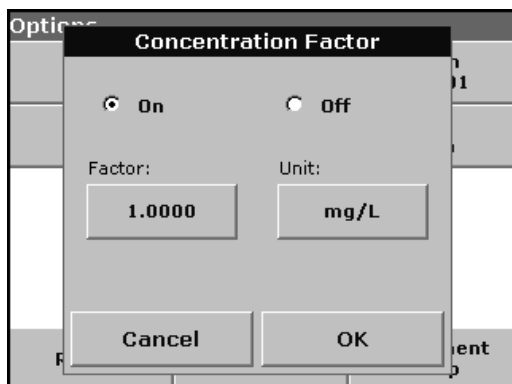

### Concentration factor:

1. Press **Concentration Factor: Off** in the Options menu. Press **On** to highlight this feature.
2. Press the "Factor" key to enter the factor by which absorbance readings are to be multiplied. Press the "Unit" key to select the units for concentration measurements or to create a new unit.
3. Press **OK** to confirm.

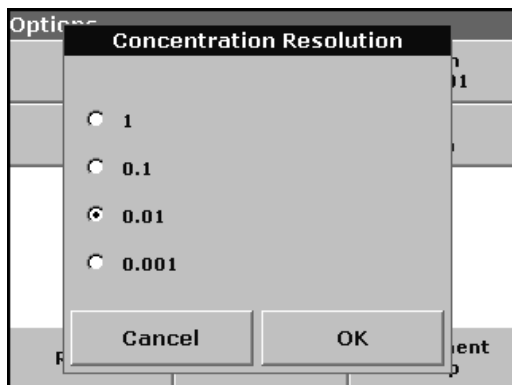

### Concentration resolution:

1. Press **Concentration Resolution** in the Options menu.
2. Select the resolution and press **OK** to confirm.

## 6.5.2 Complete a measurement in the multi wavelength mode



1. Insert the blank cuvette/cell into the cuvette/sample cell holder. Press **Zero**.

*Note: The **Read** key does not become active until the zero measurement has been completed.*

2. Insert the sample cuvette/cell into the cuvette/sample cell holder. Press **Read**.

*Note: The UV lamp is not switched on until a wavelength in the UV spectrum has been entered and the blank reading has been started. During the warm-up phase of the UV lamp, the message "Warming up..." is displayed and the "UV lamp" symbol flashes. As soon as the UV lamp is ready, the blank reading is carried out.*

*Note: **Zero** and **Read** are disabled until the cell compartment is closed.*

3. For data storage, see section 5.3.1.1 on page 35.

## 6.6 Wavelength Scan mode – recording of absorbance and transmittance spectrums

In the wavelength scan mode, the absorbance of the light in a solution over a defined wavelength spectrum is measured.

The measurement results can be displayed as a curve, as percent transmittance (%T) or as absorbance (Abs). The collected data can be printed as a table or a curve.

The data are available for formatting changes. These include automatic scaling and zoom functions. Maximum and minimum values are determined and shown as a table.

The cursor can be moved to any point on the curve for the purpose of reading off the absorbance or transmittance value and the wavelength. The data associated with each data point can also be shown as a table.

### 6.6.1 Set up the wavelength scan

Press **Wavelength Scan** in the Main Menu. Press **Options** for Parameter Setup.

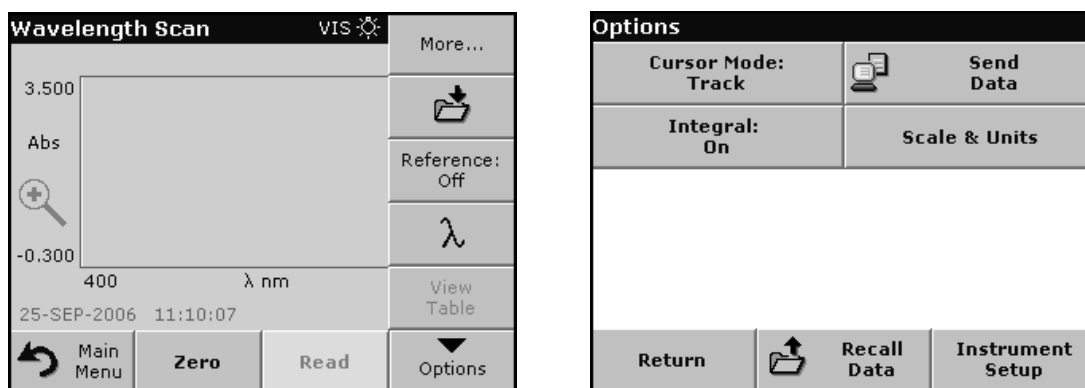


Table 14 Wavelength scan setup options

Option	Description
More	For further Options
Store icon	To store the scan data
Reference Off/On	From the displayed list of stored scans, a record is selected for use as a reference scan/superimposed scan. This can be highlighted or shown in the background in comparison with the actual measured scan. <i>Note: This option is only available when there are stored scans with the same wavelength range and step.</i>
λ	To enter the wavelength spectrum and the scan interval
Select View	Enables the user to switch the display back and forth between the scan data tables (wavelength/absorbance) and the graph of the curve. <i>Note: Select View will be activated after the first reading.</i>
Cursor Mode	To select <b>Track</b> or <b>Peak/Valley</b> . The selection for this menu item determines to which points on the graph the cursor moves.
Send Data	To send Data to a printer, computer or USB memory stick (Type A)
Integral: On/Off	The integral gives the area and the derivative of the integral gives the original function


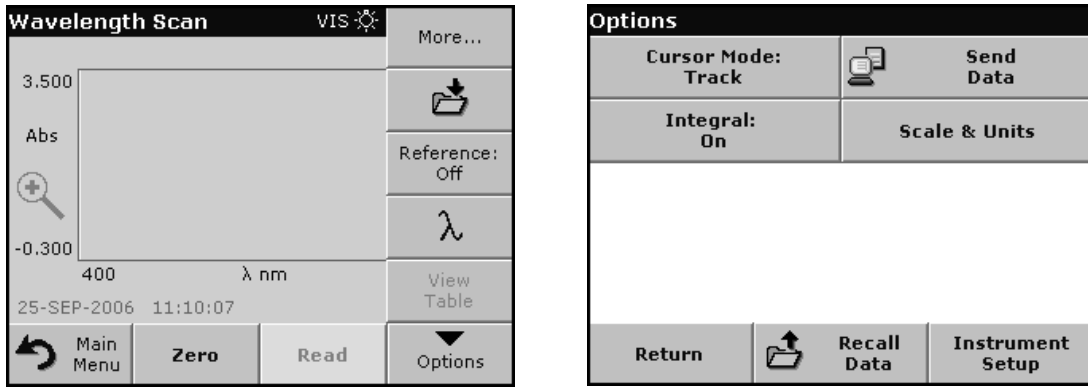


Table 14 Wavelength scan setup options (continued)

Option	Description
Scale & Units	<b>Scale:</b> In the automatic scaling mode, the y-axis is automatically adjusted so that the total scan is displayed. The manual scaling mode allows sections of the scan to be displayed. <b>Units:</b> Choice of absorbance or transmittance.
Recall Data	Call up saved measurement data, wavelength scans or time courses, see section 5.3 on page 35.
Instrument Setup	Basic data of the instrument, see section 5.2 on page 23.

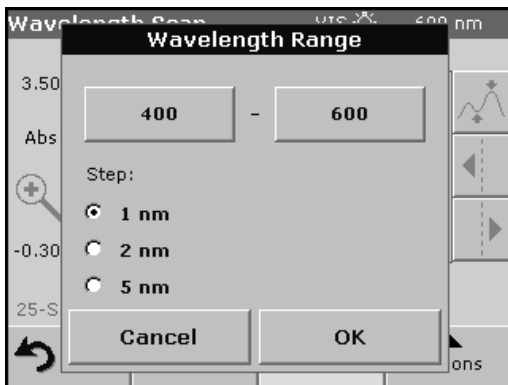

**λ Setting wavelength**

1. Press the λ key in the Options menu to select the wavelength range and the wavelength step.
2. Press the upper left key to open the numeric keypad and select the minimum wavelength. Press **OK** to confirm.
3. Press the upper right key to open the numeric keypad and select the maximum wavelength. Press **OK** to confirm.

**Note:** Do not select the same wavelength for minimum and maximum.

4. Highlight the required wavelength step.

**Note:** Scan recordings of high resolution data take a longer time than recordings of low resolution data. Selecting a larger step allows the instrument to scan faster, but decreases the resolution of the collected data.

- According to the selected wavelength range, select the interval from the actively displayed wavelength steps. In summary maximally 780 measuring steps can be accomplished during a Scan.

**Note:** The maximum wavelength adjusts automatically if the difference between the maximum and minimum wavelength is not a multiple of the interval.

- Press **OK** to return to the scan mode. Selected parameters are displayed along the graph's x-axis.

**Wavelength Scan** VIS 400 nm


nm	Abs	Min/Max	nm	Abs	Min/Max
400	0.515		401	0.509	
402	0.538		403	0.523	Valley
404	0.530		405	0.557	Peak
406	0.514		407	0.481	Valley
408	0.564		409	0.616	
410	0.634	Peak	411	0.580	Valley

Integral: 323.54

Main Menu Zero Read Options


### Select view (displaying table)

- Press **Select View** in the Options menu after a reading is taken.
- A table with the results is displayed.
- To return to the graph press **Options** and then **View Graph**.

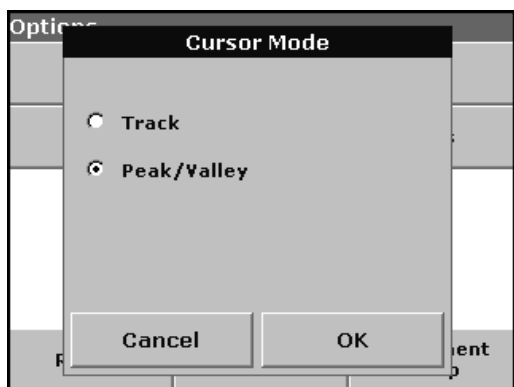

### Cursor mode

- Press **Cursor Mode: Track** in the Options menu.
- The selection for this menu item determines what data are displayed in the table. Highlight **Track** or **Peak/Valley**.
- Press **OK** to confirm.
- Press **Return** to return to the scan mode.

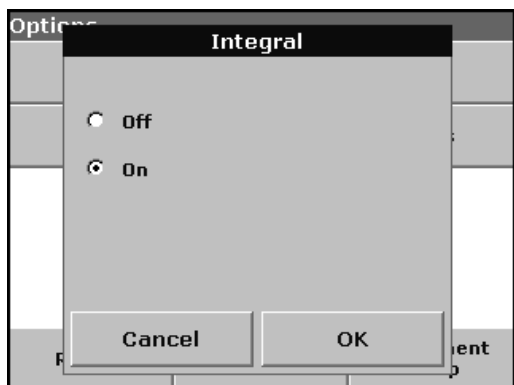

### Integral

The Integral applies to the whole wavelength range of the scan.


- Press **Integral: Off** in the Options menu.
- Highlight **On** to show the Integral. To find the integral of other wavelength ranges, change the wavelength range and scan again.
- Press **OK** to confirm.



4. Press **Return** to return to the scan mode.

*Note: The Integral is shown instead of the date on the display.*

*Note: For the next scan measurement the setting for the Integral will be On.*

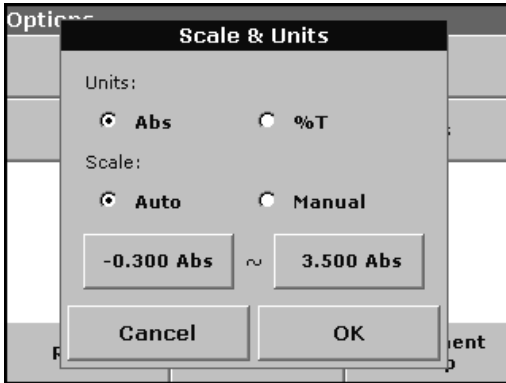

### Scale & units


1. Press **Scale & Units**.
2. Highlight the required units (**Abs** or **%T**).
3. Highlight **Auto** or **Manual** scaling on the graph's y-axis.

*Note: If manual scaling is selected, use the alphanumeric keypad to set the limits  $y_{min}$  and  $y_{max}$ . The graph is adjusted to display only the values in the selected range. If automatic scaling is selected, the instrument sets the limits automatically so that the total range can be displayed.*


4. Press **OK**.
5. Press **Return** to return to the scan mode.

## 6.6.2 Wavelength scan reading


After the scanning parameters have been selected, the baseline must be scanned. Changing any of the scanning parameters requires a new baseline scan. When the baseline has been scanned, the instrument is ready to scan one or more samples.

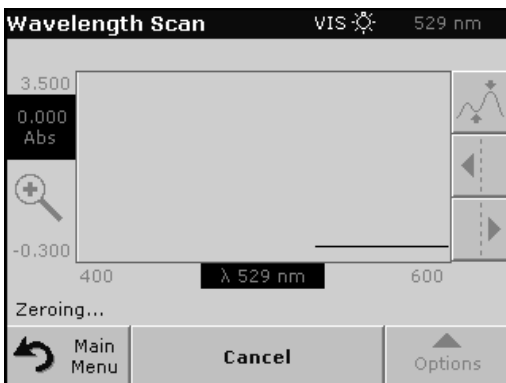

1. Press **Wavelength Scan** in the Main Menu.
2. Insert the blank cuvette/cell into the cuvette/sample cell holder.

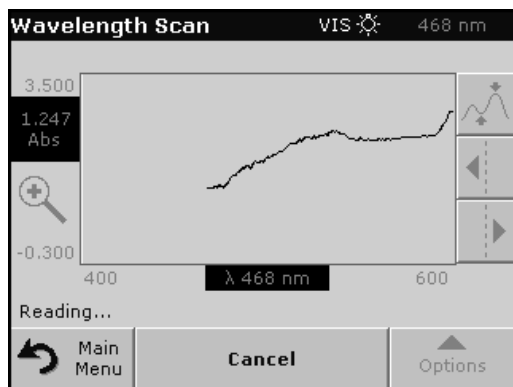

3. Press **Zero**. "Zeroing...." appears below the graph as the baseline scan begins.
4. Insert the prepared sample cuvette/cell into the cuvette/sample cell holder.
5. Press **Read**. Below the graph "Reading..." appears and a graph of the absorbance or transmittance values at the scanned wavelengths is displayed continuously.

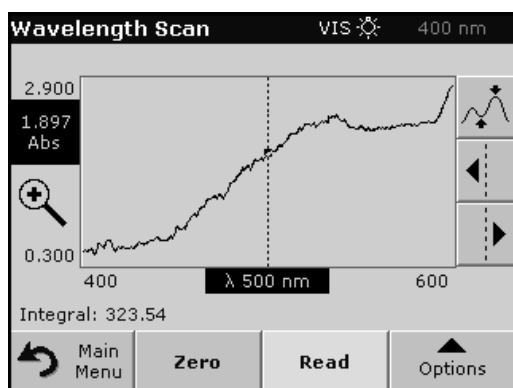

The Wavelength Scan is complete, if

- the graph is shown full size,
- the scaling of the x-axis fits automatically,
- the Cursor functions in the vertical navigation bar are highlighted.


## 6.6.2.1 Navigation of the wavelength scan graph or a wavelength scan analysis

**Table 15 Navigating the wavelength scan**

Cursor Function/ Zoom Function	Description
<b>Curve Icon (Choice of Cursor Mode)</b>	Choice of Cursor Mode <b>Peak/Valley</b> (cursor moves between minimum/maximum absorbance values) or Cursor Mode <b>Tracking</b> (cursor moves over each data point of the scan).
<b>Arrow keys</b>	The arrow keys are used (right/left) to move the cursor (depending on the selected mode) to the next data point. The data of the data point (wavelength/absorbance or transmittance value) are highlighted on the x and y axes. <b>Note:</b> Press any point on the curve to display the associated data.
<b>Zoom Icon</b>	This function is used to magnify the section of the curve in the vicinity of the cursor. The original curve size can be restored by pressing the zoom icon again.

### 6.6.3 Work with reference scans

There are two options to work with **Reference Scan**:

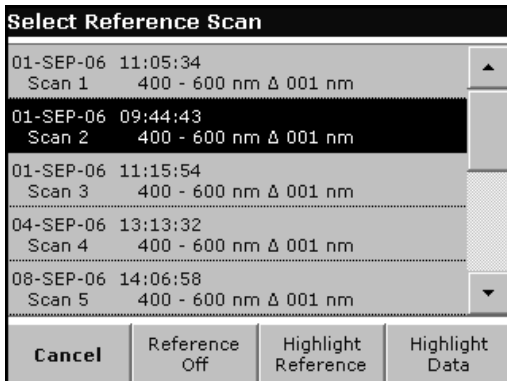

#### Option 1:


1. Press **Reference: Off** in the Options menu to select another scan to display on the same screen with the current scan. Highlight the required scan number and press **Highlight Reference**.

**Note:** After selecting a reference scan the **Reference: Off** key in the Options menu turns into **Reference: On**.


**Note:** Only scans that have the same wavelength range and step can be displayed using the overlay option. A reference curve can only be displayed with another active curve. Two previously stored scans cannot be overlaid.

2. The reference curve is shown in gray. The absorbance or transmittance value and the associated wavelength are highlighted in gray.

**Note:** A black and a gray box are shown in the left corner of the display. The gray box relates to the reference scan and the black one relates to the current wavelength scan.

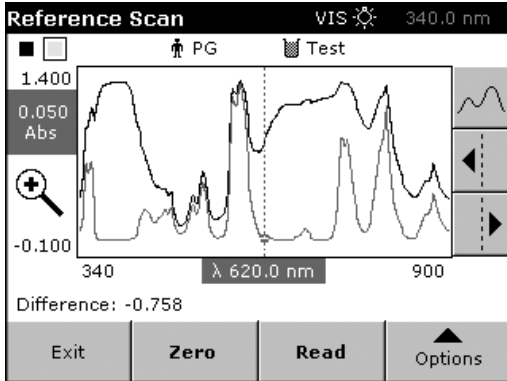

3. To complete the wavelength scan reading, see section 6.6.2 on page 97.
  - The newly plotted wavelength scan curves are shown in black.
  - The absorbance or transmittance value and the associated wavelength are highlighted in black.
  - In addition, the display shows the difference between the wavelength scan curve and the reference curve against the wavelength.
4. Press the black or gray small box in the left upper corner on the screen to switch between the actual wavelength scan and reference scan.

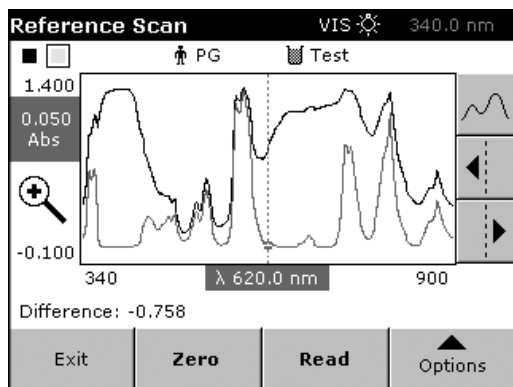

### Option 2:

1. Insert the blank cuvette/cell into the cuvette/sample cell holder. Press **Zero**.
2. Insert the sample cuvette/cell into the cuvette/sample cell holder. Press **Read**.

- The newly plotted wavelength scan curves are shown in black.
- The absorbance or transmittance value and the associated wavelength are highlighted in black.

3. Press **Options** and then **Reference: Off** in the Options menu to select another scan to display on the same screen with the current scan. Highlight the required scan number and press **Highlight Reference**.

**Note:** After selecting a reference scan the **Reference: Off** key in the Options menu turns into **Reference: On**.

**Note:** Only scans that have the same wavelength range and step can be displayed using the overlay option. A reference curve can only be displayed with another active curve. Two previously stored scans cannot be overlaid.

4. The reference curve is shown in gray. The absorbance or transmittance value and the associated wavelength are highlighted in gray.
  - Additionally the difference of the extinction and/or transmission value between the two Scans (measured Scan and reference Scan) is indicated/highlighted at each position of the cursor.

**Note:** A black and a gray box are shown in the left corner of the display. The gray box relates to the reference scan and the black one relates to the current wavelength scan.


5. Press the black or gray small box in the left upper corner on the screen to switch between the actual wavelength scan and reference scan.

## 6.7 Time course of absorbance/transmittance

The Time Course Mode is used to collect data in either absorbance or transmittance for a user-specified length of time. After the data are collected, they can be displayed in either graphic or tabular format.

### 6.7.1 Time course setup parameters

Press **Time Course** mode in the Main Menu. Press **Options** to configure parameters.

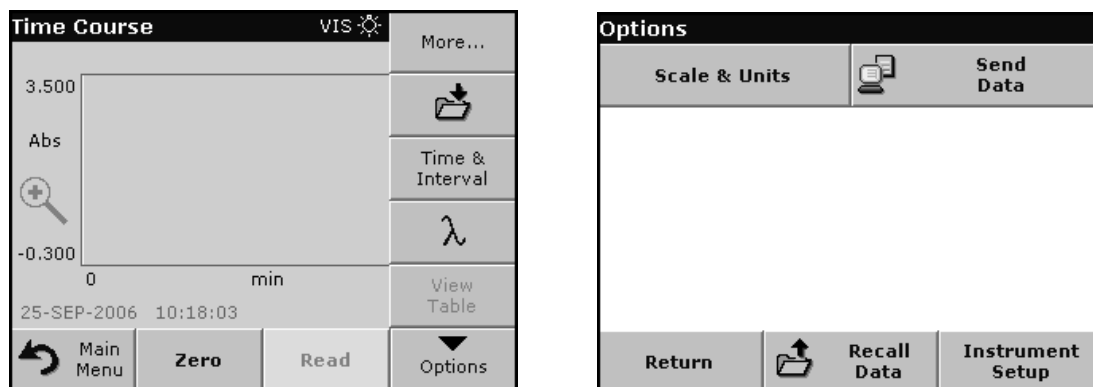

**Table 16 Time course setup options**


Option	Description
<b>More</b>	For further Options
<b>Store icon</b>	To store the scan data
<b>Time &amp; Interval</b>	To input the total time for data collection and the time interval between the collection of data points
$\lambda$	To input the wavelength setting
<b>View Table/Graph</b>	To display readings in absorbance, transmittance or concentration. This can be changed after sample data are collected
<b>Scale &amp; Units</b>	<b>Scale:</b> In the automatic scaling mode, the y-axis is automatically adjusted so that the total scan is displayed. The manual scaling mode allows sections of the scan to be displayed. <b>Units:</b> Choice of absorbance or transmittance.
<b>Send Data</b>	To send Data to a printer, computer or USB memory stick (Type A)
<b>Recall Data</b>	Call up saved measurement data, wavelength scans or time courses, see section 5.3 on page 35.
<b>Instrument Setup</b>	Basic data of the instrument, see section 5.2 on page 23.



### Time & interval:

1. Press **Time & Interval** in the Options menu.
2. Input the total time and the reading time and press **OK** to confirm.

**Note:** In total 500 measuring steps are possible. To select a total time and a time interval that would cause this number of measurements to be exceeded, the time interval is defined automatically and the **OK** key is inactivated.

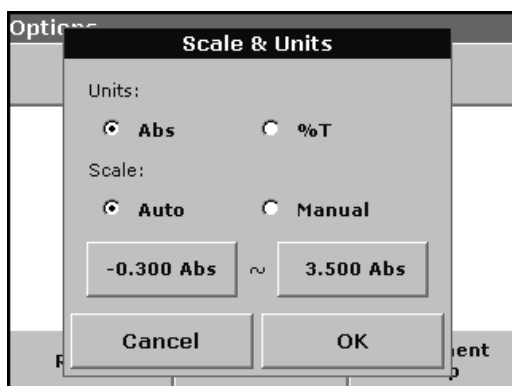


### Scale & units:

1. Press **Scale & Units** in the Options menu.
2. Highlight **Abs** or **%T** as the required units.
3. Highlight **Auto** or **Manual** scaling on the graph's y-axis.

**Note:** If manual scaling is selected, use the alphanumeric keypad to set the limits  $y_{min.}$  and  $y_{max.}$ . The graph is adjusted to display only the values in the selected range. If automatic scaling is selected, the instrument sets the limits automatically so that the total range can be displayed.

4. Press **OK** to confirm.
5. Press **Return** to return to the scan mode.


## 6.7.2 Time course scan reading

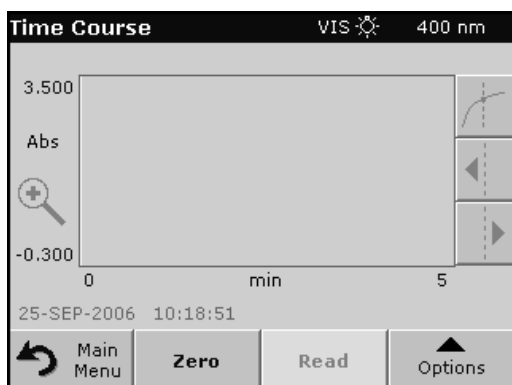

After the parameters have been selected, the instrument must be blanked, then the sample can be analyzed.

1. Insert the blank cuvette/cell into the cuvette/sample cell holder. Press **Zero**. The blank reading is shown on the display.
2. Insert the sample cuvette/cell into the cuvette/sample cell holder. Press **Read**. Start collecting time course (kinetic) data.


**Note:** During the measurement the **Zero** and **Read** keys change to **Mark** and **Stop**.

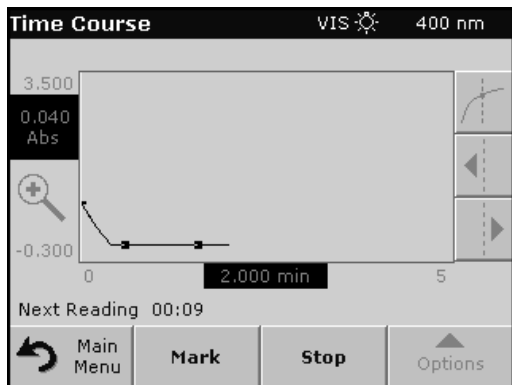

- Select **Mark** to mark the next data point collected. This mark is not used by the instrument, but is available for the user and may indicate a significant event, such as the addition of a sample or other reagent. The mark is also shown in the table.
- Select **Stop** to stop taking sample readings.

### 6.7.3 Analysis of time course data

After the data are collected, the following manipulations can be done on the graphic data:

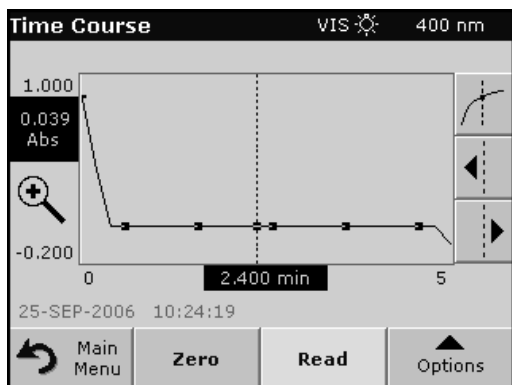

The **Time Course** Program is complete, if

- the sound is turned on, the instrument beeps when the readings are done
- the graph is shown fullsize,
- the scaling of the x-axis occurs automatically,
- the Cursor functions in the vertical navigation bar are highlighted.

### 6.7.3.1 Navigation of a time scan or a time scan analysis

After a time scan has been completed, the time and the absorbance/transmittance data are displayed as a curve.


Where the cursor is positioned on the curve, the elapsed time up to this point and the corresponding absorbance are highlighted.

**Table 17 Navigating the time scan**


Cursor Function/ Zoom Function	Description
<b>Curve Icon (Choice of Cursor Mode)</b>	<b>Delta mode:</b> A second cursor is highlighted. The position of the fixed cursor was previously defined in Cursor Mode Single. Use the active cursor to select any point on the measurement curve. The difference to the fixed cursor is shown on the curve. The delta values are correspondingly highlighted and displayed on the x and y axes. The gradient of the curve and the correlation coefficient ( $r^2$ ) between the cursor points in the Delta mode are shown under the curve.
	<b>Cursor Mode Single:</b> The cursor moves to each selected measurement point of the scan.
<b>Arrow keys</b>	The arrow keys (right/left) are used to move the cursor (depending on the selected mode) to the next data point. The data of the data point (wavelength/absorbance or transmittance value) are highlighted on the x and y axes. <i>Note: Press any point on the curve to display the associated data.</i>
<b>Zoom Icon</b>	This function is used to magnify the section of the curve in the vicinity of the cursor. The original curve size can be restored by pressing the zoom icon again.



## 6.8 System checks

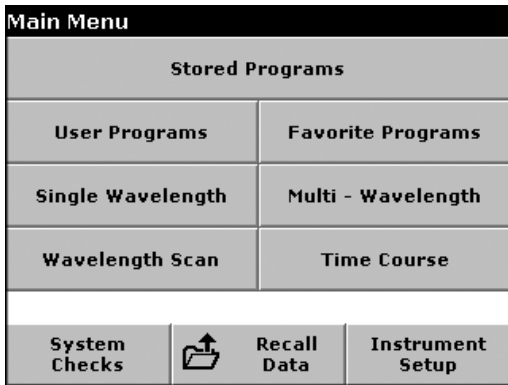


1. Press **System Checks** in the Main Menu.

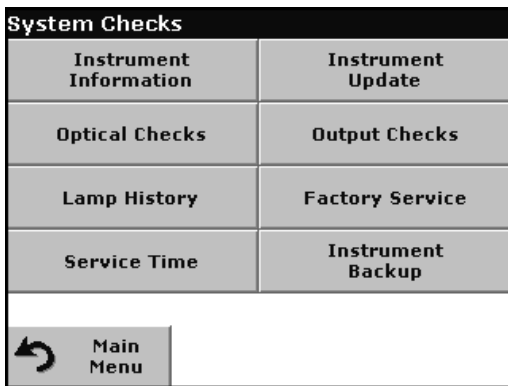

The System Checks menu contains instrument information and various performance tests.

### 6.8.1 Instrument information

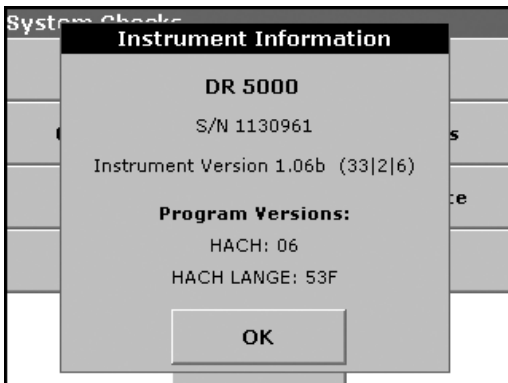

1. Press **Instrument Information** in the System Checks menu.
2. The model, serial number and software version are displayed.

### 6.8.2 Upgrade of the instrument software

To obtain the software for the update from the Internet at [www.hach.com](http://www.hach.com):




1. Go to <http://www.hach.com>.
2. On the DR 5000 product page, click **Lab System Software/Software Update Downloads** under the heading **Downloads**.
3. Locate the appropriate download and follow the prompts for saving the file(s) to the USB memory stick.
4. On the DR 5000, press **Instrument Update** in the System Checks menu.
5. Connect the USB stick to the USB interface on the DR 5000 (section 3.3 on page 14). Press **OK**. The link is established automatically and the software is updated.
6. Press **OK** to return to the System Checks menu.

**Note:** When the instrument software has been updated, a prompt to restart the instrument is displayed.

### 6.8.3 Optical checks

For each optical check, the measured results must be evaluated against user-specific requirements. The check options do not define tolerances.

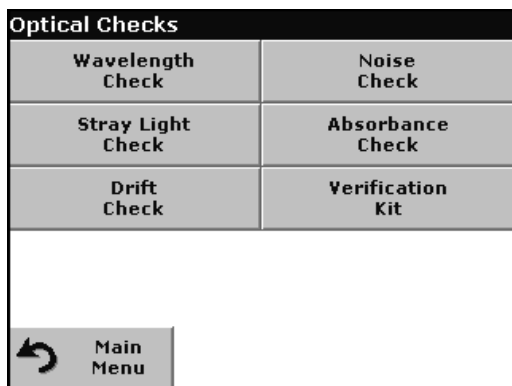


1. Press **Optical Checks** in the System Checks menu.

The Optical Checks menu contains programs for checking the wavelength accuracy, stray light and photometric accuracy.


An optional test filter set (Verification Kit) (Section 9 on page 129) containing six precision glass filters, target values, tolerances and instructions is available and is recommended as an aid for carrying out comprehensive in-house instrument checks.

### 6.8.3.1 Wavelength check

The Wavelength Check test is used to check wavelength accuracy at 656.1 and 486.0 nm and to check bandwidth at 656.1 nm, only.

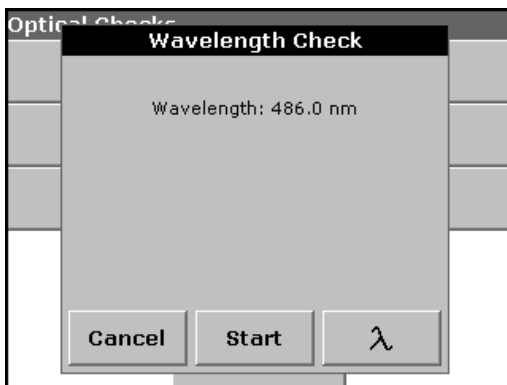


1. Press **Wavelength Check** in the Optical Checks menu.
2. Press  $\lambda$  to select the wavelength.

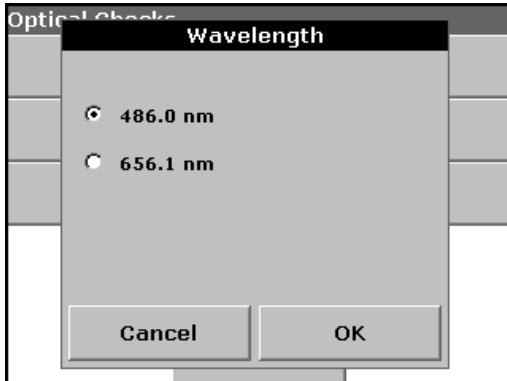

3. Make a choice and confirm with **OK**.
4. Press **Start**.



**Note:** During the warm-up phase of the UV lamp, the message "Lamp Warmup..." is displayed.

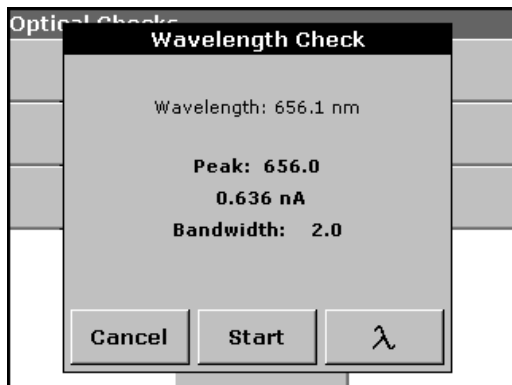


5. The result is displayed.
6. Press **Cancel** to return to Optical Checks.

### 6.8.3.2 Noise check

The Noise Check test is used to test the photometric noise in the instrument.


This test can be used to test noise at any wavelength and at an absorbance level determined by a specific sample.

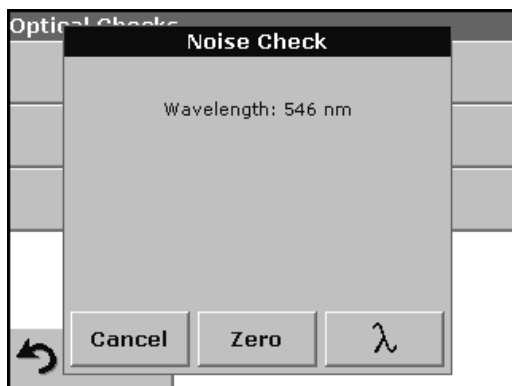


1. Press **Noise Check** in the Optical Checks menu.
2. Press  $\lambda$  to input the wavelength.
3. Input the wavelength and press **OK** to confirm.



4. Press **Zero**.

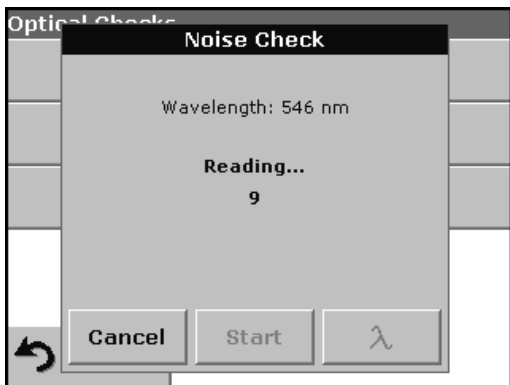

5. Insert sample cuvette/cell into the cell compartment. Close the cell compartment.
6. Press **Start**.

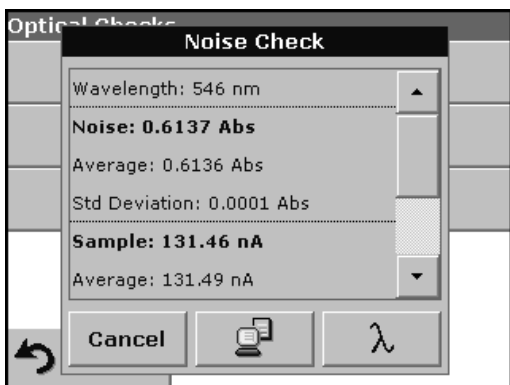


7. The result will be displayed. Thirty readings are averaged for the blank. The "Average" and "Std Deviation" are calculated from 100 consecutive absorbance readings (Noise, Sample, Reference).
 

*Note: Failings and passings will be defined by the user.*
8. Press **PC & Printer** icon to send data to a PC or Printer.
9. Press **Cancel** to return to Optical Checks.

### 6.8.3.3 Stray light check

The Stray Light test is used to measure the stray light in the instrument at 220.0 nm and 340.0 nm.

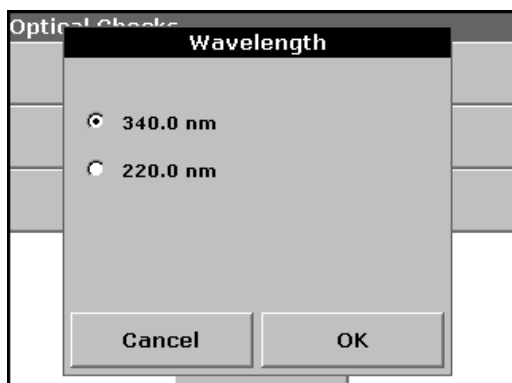


1. Press **Stray Light Check** in the Optical Checks menu.
2. Press  $\lambda$  to select the wavelength.
3. Make a choice and confirm with **OK**.



4. Press **Zero**.



5. Insert the sample cuvette/cell or reference filter into the cell compartment. Close the cell compartment. Press **Start**.

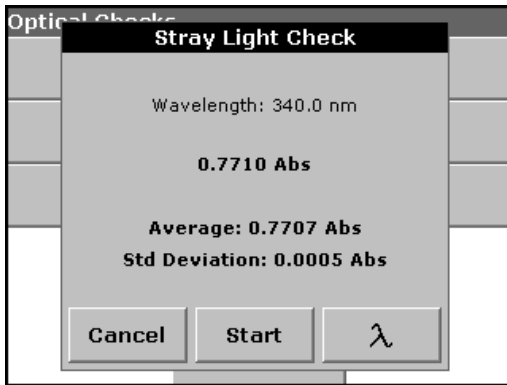


- The "Average" and "Std Deviation" are calculated from 100 successive absorbance measurements. The result is displayed and is to be compared with the nominal/standard data (given in the quality control certificate) of the sample cell/test filter.

*Note: Failings and passings will be defined by the user.*

- Press **Cancel** to return to Optical Checks.


#### 6.8.3.4 Absorbance check

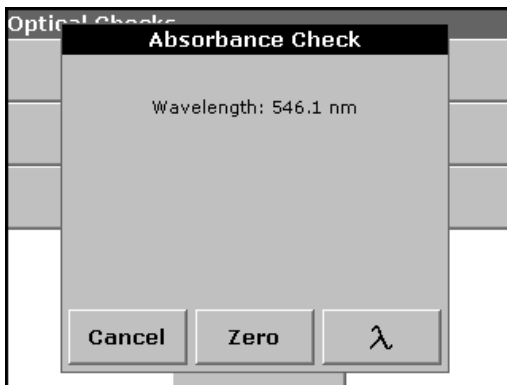


The Absorbance Check test is used to test the photometric accuracy and repeatability of the instrument.

This test can be used to test absorbance at any wavelength by a specific sample or test filter set (Refer to section 6.8.3.6 on page 113).

- Press **Absorbance Check** in the Optical Checks menu.
- Press  $\lambda$  to input the wavelength.
- Input the wavelength and press **OK** to confirm.
- Remove any cuvette/sample cell from the cell compartment, close the cell compartment, and press **Zero**.

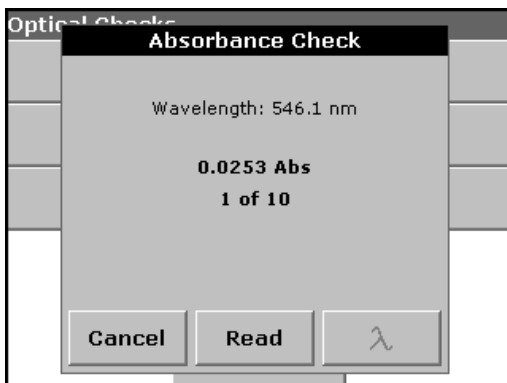

- Insert sample cuvette/cell or test filter into the cell compartment, close the cell compartment, and press **Read**.
- Repeat steps 4 and 5 for a total of 10 replicates to obtain final results.

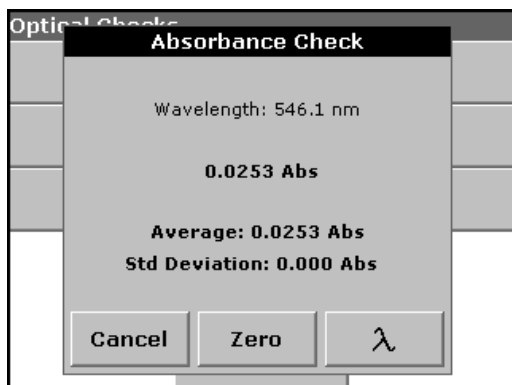


7. The result is displayed and is to be compared with the nominal/standard data of the test filter. If a specific sample is used, the measured results must be evaluated against user-defined requirements.
8. Press **Cancel** to return to Optical Checks.


### 6.8.3.5 Drift check

The Drift Check test is used to test the stability of the instrument.

**Note:** *The Drift Check runs 1 hour.*

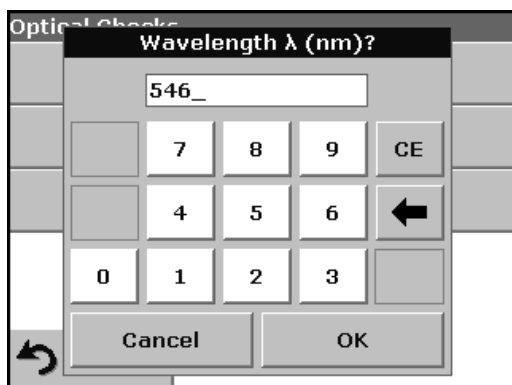


1. Press **Drift Check** in the Optical Checks menu.
2. Press  $\lambda$  to input the wavelength.
3. Input the wavelength and press **OK** to confirm.

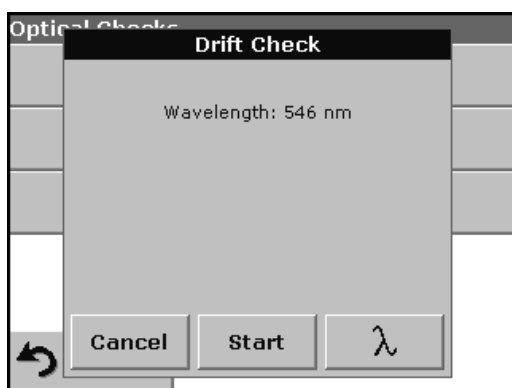


4. Press **Start**.



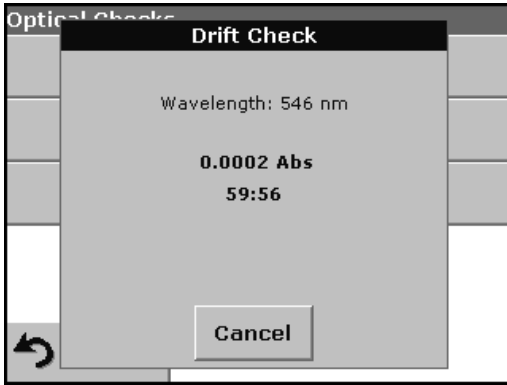

5. The Drift Check runs 1 hour.

This test takes a reading every minute for one hour. Every 15 minutes, linear regression is used to calculate the slope (rate of change) for the previous 15 minute interval.

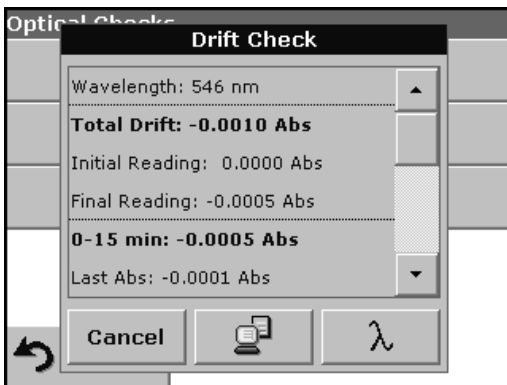


6. The last reading and slope for each 15 minute interval are displayed. At the end of the hour, the overall values are calculated and displayed. The test proceeds to completion.

7. Press **Cancel** to return to Optical Checks.

### 6.8.3.6 Verification kit


The Verification Kit, LZV 537, (see Section 9 on page 129) is designed for periodic monitoring of scattered light, photometric accuracy and the wavelength accuracy of the spectrophotometers.

When results exceed allowable tolerances (given in the quality control certificate), contact the manufacturer.

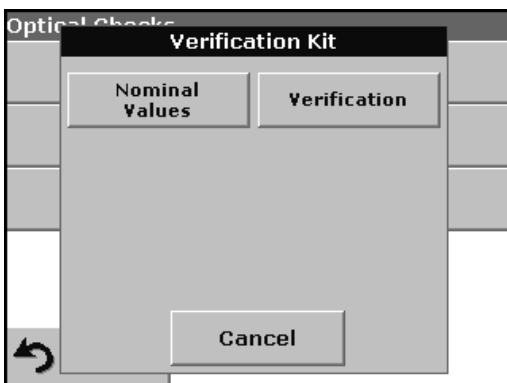

1. Press **Verification Kit** in the Optical Checks menu.

2. Press **Nominal Values**.

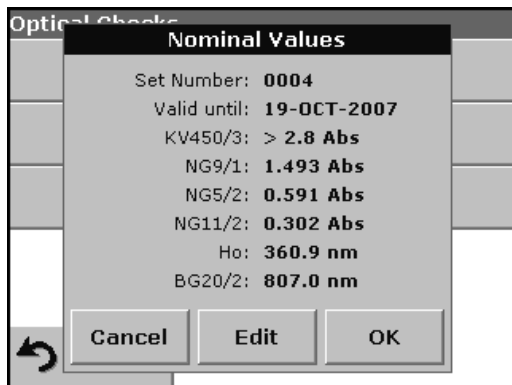


3. Press **Edit**.

An automatic menu guidance queries values (filters, wavelength, nominal values and tolerances) given in the quality control certificate, to the following specifications:

- Stray Light
- Photometrical accuracy
- Wavelength accuracy

4. Press **OK** when all values are entered and the overview is displayed.

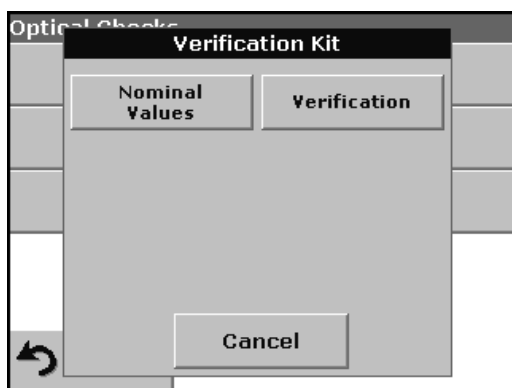

5. Press **Verification**.


6. Insert the Multi Cell Holder in the cell compartment in such a way that the cell opening for 10 mm square cuvettes/cells is at the front. Secure the holder with two locking screws.



7. Remove any cuvettes/cells from the cell compartment, close the cell compartment, and press **Start**.



8. Insert the different filter in the given order one after the other with the filter glass facing the front of the instrument. Press **Next** after inserting a filter and closing the cell compartment

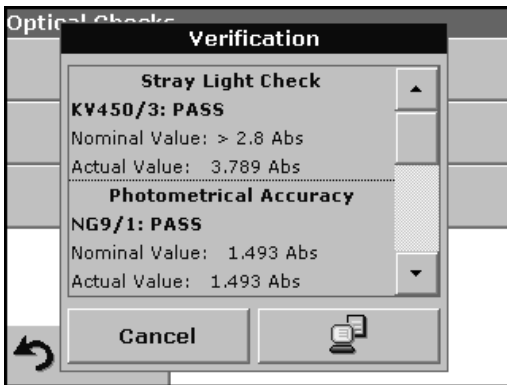

After the last measurement the results are displayed.

9. Press **PC & Printer** icon to send the data to a USB memory stick, PC or to a printer.

The files will be stored automatically as CSV file (Comma Separated Value). The file name will be formatted as "Verification.csv".


### 6.8.4 Output checks

If a printer is connected a test printing of the current screen will be printed.

### 6.8.5 Lamp history

The Lamp History menu provides the following information:


- The amount of time that the lamp has been on (Hours)
- The number of times that the lamp has been turned on (Cycles)
- The number of times that the instrument tried to turn the UV lamp on (triggers)

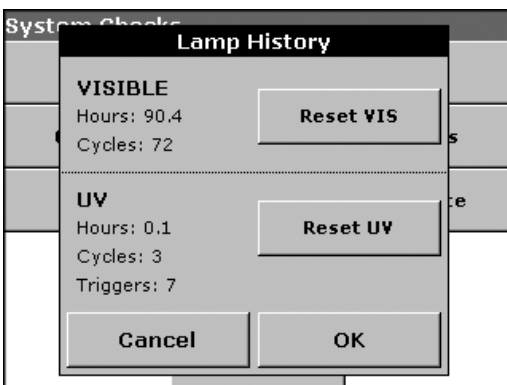

After a lamp is replaced and the Lamp History is reset, the display of the total operating time is reset to 0.

1. Press **Lamp History** in the System Checks menu.
2. Press **Reset VIS** and the Visible Lamp will be reset.
3. Press **Reset UV** and the UV Lamp will be reset.
4. Press **OK** to return to System Checks.

### 6.8.6 Factory service

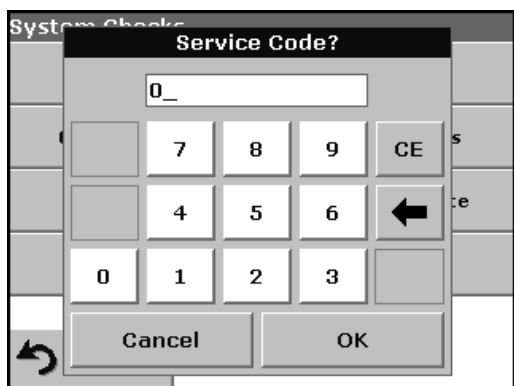

The Factory Service menu is password protected. This menu is not intended for customer use.

### 6.8.7 Service time

In order to ensure a regular inspection, an automatic memory reference for the service times can be entered. After switching the instrument on this memory reference will be activated and indicated at the appropriate time.



1. Press **Service Time** in the System Checks menu.
2. Select **On** and then **Last Service** to enter the date of the last inspection.
3. Press **OK** to confirm.



4. Select **Next Service** to determine a specific period of time up to the next inspection.
5. Press **OK** to confirm.




If the next service is due, the message "**Next service is due!**" is displayed after switching on the instrument.

6. Press **OK** to return to the Main Menu.

Contact the manufacturer or distributor to arrange an appointment for the next service.

### 6.8.8 Instrument Backup

Before the next service date the Instrument Backup menu offers the possibility to store all programs, measuring data, Operator ID, Sample ID, passwords and all adjustable data on a USB stick.

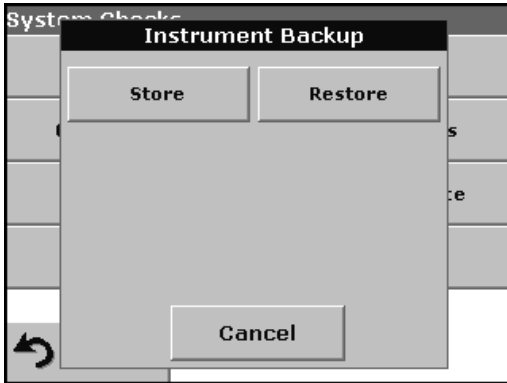

1. Press **Instrument Backup** in the System Checks menu.
2. Connect the USB memory stick (section 3.3 on page 14).
3. Press **Store** to start a Backup.



**Note:** If the USB stick is not connected, the message "Please insert USB Memory" is displayed. Connect a USB stick, in order to store the data. Press **OK** to confirm and press **Store** again.




**Note:** If the Backup was already stored before, the message "Data already exists. Overwrite?" is displayed. Press **OK** to overwrite the data.



If the file was stored the message Instrument Backup is stored to USB stick will be displayed.

4. Press **OK** to return to the System Checks menu.

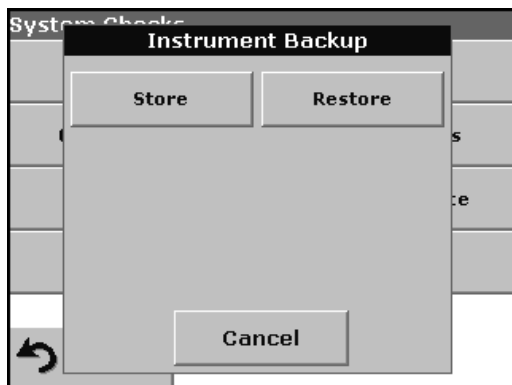

### Restore backup data:

**Important Note:** All current data will be overwritten when restoring the Backup file!


1. Press **Instrument Backup** in the System Checks menu.
2. Connect the USB memory stick containing the Backup (section 3.3 on page 14).
3. Press **Restore** to pass back the data.



**Note:** If the USB stick is not connected, the message "Please insert USB Memory" is displayed. Connect a USB stick, in order to store the data. Press **OK** to confirm and press **Restore** again.



4. Press **OK** to confirm after the message "Instrument Backup from S/N XXXXXXXX. Restore?" is displayed.

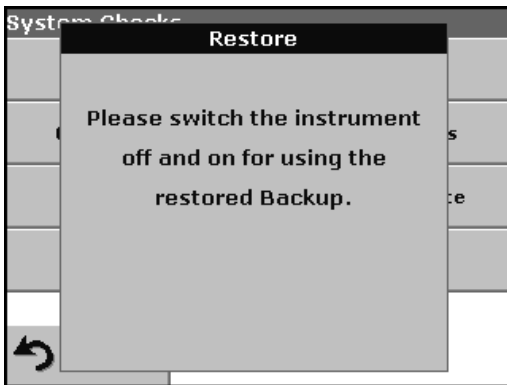

5. After the backup start the instrument again.





### **CAUTION**

**Potential Chemical, Biological Eye and Skin Hazards.**  
**Only qualified personnel should conduct the tasks described in this section of the manual.**

**Important Note:** Remove any cuvettes/cells that are still in the instrument and dispose of them or their contents using an approved disposal method.

## 7.1 Cleaning requirements

### **CAUTION**

**Potential Pinch, Eye, Burn and Chemical Hazards.**  
**Always disconnect power from the instrument before attempting any cleaning operations.**

**Important Note:** Under no circumstances should the instrument, display or the accessories be cleaned with solvents such as white spirit, acetone, etc.

### 7.1.1 Spectrophotometer

- Clean the enclosure, cuvette/sample cell compartments and all accessories with a soft damp cloth. A mild soap solution can also be used. Do not get excess water in the cuvette/sample cell compartments. Do not insert a brush or sharp object into Cell Compartment #1 to avoid damaging the mechanical components.
- Dry the cleaned parts carefully with a soft cotton cloth.

### 7.1.2 Display

- Take care not to scratch the display. Do not touch the screen with ball pens, pencils or similar pointed objects.
- Clean the display with a soft, lint-free and oil-free cotton cloth. Diluted window cleaner liquid can also be used.

### 7.1.3 Cuvettes/sample cells

#### **CAUTION**

**Potential Chemical/ Biological Exposure Hazards.**  
**Use proper laboratory practices whenever there is a risk of chemical exposure.**

1. After performing a procedure, clean glass cuvettes/sample cells with cleaning agents.
2. Afterwards, rinse the cuvettes/sample cells several times with tap water and then thoroughly with deionized water.

**Important Note:** Glass cuvettes/sample cells that have been used for organic solvents (such as chloroform, benzene, toluene, etc.) must be rinsed with acetone before being treated with cleaning agents. In addition, another rinse with acetone is necessary as a final treatment step before the cuvettes/sample cells are dried.

### 7.2 Lamp replacement

#### **CAUTION**

***The UV lamp generates UV light. Do not look directly at an operating lamp without wearing UV protective eye glasses.***

The Lamp compartment is on the left side behind the display and is provided with ventilation on the back side. The tungsten and deuterium (UV) lamp are installed in the lamp compartment. On the back side a fan is installed for cooling of electric components. The ventilation system operates automatically.

Refer to section 7.2.1 on page 122, section 7.2.2 on page 123 and Figure 8 on page 124 for instructions on replacing the tungsten lamp and deuterium lamp (UV).

#### 7.2.1 Tungsten lamp replacement

1. Switch the instrument off.
2. Unplug the power cord.

#### **DANGER**

***The instrument must remain switched off until the change of lamp has been completed.***

#### **WARNING**

***Burn Hazard. Wait until the lamp cools down. Contact with the hot lamp can cause burns.***

3. Use a screwdriver to remove the cover from the back of the instrument (the screws may be slotted or cross-headed).
4. Place the cover and the attached fan carefully beside the instrument (take special care with the fan cable).
5. Push the spring (Figure 8 on page 124, item 2) up and remove the tungsten lamp (Figure 8, item 1) from the lamp compartment.
6. Unplug plug contact (Figure 8, item 7) from the tungsten lamp.

***Important Note: The lamp should only be held by the lamp socket. Avoid touching the glass.***

7. Push plug contact firmly onto the new tungsten lamp.
8. Insert the tungsten lamp into the lamp compartment. Push the spring down.
9. Check that the spring and the lamp socket are positioned correctly.
10. Use a screwdriver to screw the back cover onto the instrument again.
11. Plug in power supply again.

12. Switch the instrument on.
13. Reset the Lamp History, see section 6.8.5 on page 115.

## 7.2.2 Deuterium lamp (UV) replacement

1. Switch the instrument off.
2. Unplug the power cord.

### **DANGER**

***The instrument must remain switched off until the change of lamp has been completed.***

### **WARNING**

***Burn Hazard. Wait until the lamp cools down. Contact with the hot lamp can cause burns.***

3. Use a screwdriver to remove the cover from the back of the instrument (the screws may be slotted or cross-headed).
4. Place the cover and the attached fan carefully beside the instrument (take special care with the fan cable).
5. Unplug the deuterium lamp (Figure 8 on page 124, item 4) from the socket by pushing down on the safety contact.
6. Use a screwdriver to unscrew the two fastening screws (Figure 8, item 3) (the screws may be slotted or cross-headed) out of the socket.
7. Holding the lamp socket, lift the deuterium lamp up out of the lamp compartment (remove the complete unit, including the cable).
8. Carefully insert the new deuterium lamp into the lamp compartment.

***Important Note: Do not touch the glass envelope on the new lamp. If it is touched, clean with alcohol.***

9. Screw in both fastening screws again until they are finger-tight.
10. Insert the deuterium lamp in the socket so that the safety contact clicks into place.
11. Use a screwdriver to screw the back cover onto the instrument again.
12. Plug in power supply again.
13. Switch the instrument on.
14. Reset the Lamp History, see section 6.8.5 on page 115.

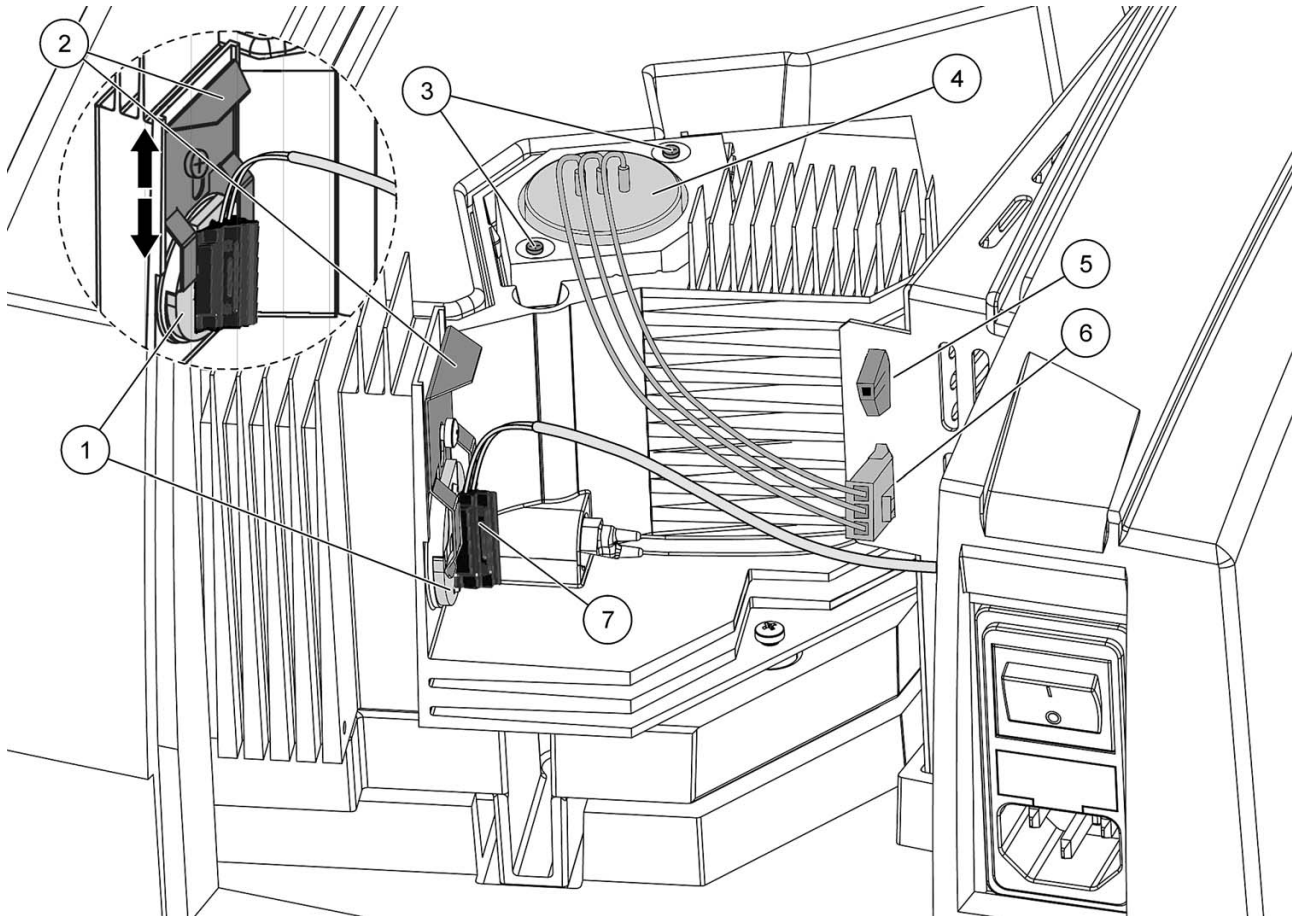


Figure 8 Lamp compartment

1 Tungsten lamp	5 Plug contacts for fan
2 Spring	6 Plug contact for deuterium lamp
3 Screws	7 Plug contact for tungsten lamp (with grip and cable)
4 Deuterium lamp (lamp socket)	

### 7.3 Fuse replacement

**DANGER**

*Certain electrical circuits within this equipment are protected by fuses against over-current conditions. For continued protection against a risk of fire, replace fuses only with the same type and rating specified.*

**DANGER**

*Failed fuses are generally an indication of a problem with the equipment. If fuse failure continues, contact a service representative for instructions on how to return the equipment for repair. Do not attempt to repair the equipment yourself.*

1. Switch the instrument off.
2. Unplug the power cord to remove all power from the instrument.

3. Pull out the plastic base over the power cable socket.
4. Remove the defective fuse, see Figure 1 on page 14.
5. Insert new fuse (T 2 A H; 250 V)
6. Replace the plastic base.
7. Plug in the power cord.
8. Switch the instrument on.

## **7.4 Filter pad maintenance**

To determine when the filter mat needs to be replaced, inspect the filter mat every 3–6 months (in a relatively dust-free environment, this interval can be longer).

1. Remove any cuvettes and cuvette/sample cell adapter from the cell compartment.
2. Turn the instrument off.
3. Unplug the power cord.
4. Lift the instrument and check the color of the filter mat. Replace the filter mat if is dark gray or black.

### **7.4.1 Filter pad replacement**

5. Carefully turn the instrument over and place it on a soft surface.
6. Use a screwdriver (standard or cross-head) to open the filter grid (Figure 9 on page 126, item 1).
7. Lift the filter grid (Figure 10 on page 126, item 1).
8. Remove the old air filter pad and replace it with a new air filter pad (Figure 10, item 2).
9. Screw the grid back in place.
10. Carefully stand the instrument upright.
11. Plug the instrument in.

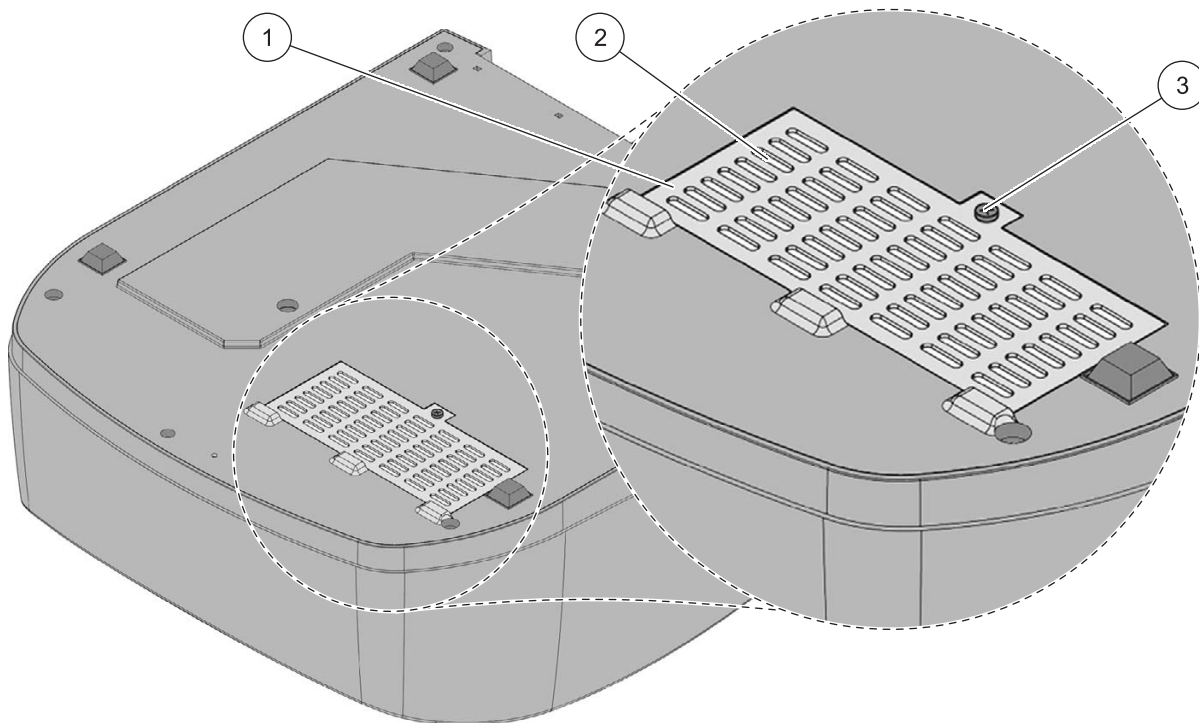


Figure 9 Base of DR 5000 with filter grid

1 Filter grid	3 Phillips screw
2 Filter pad	

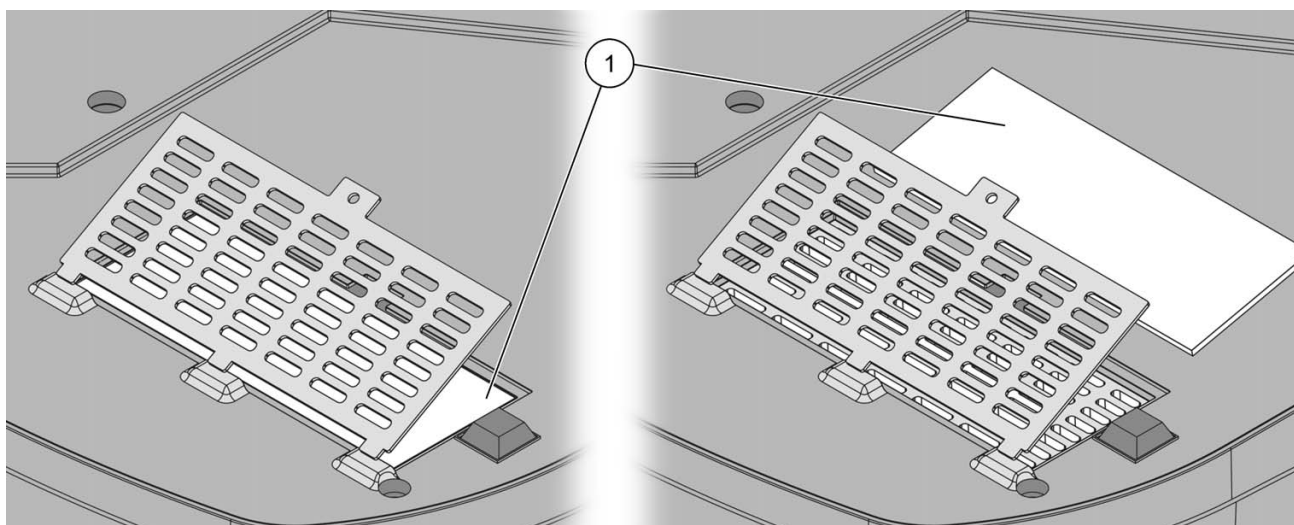


Figure 10 Change of filter pad

1 Filter pad
--------------

## Section 8 Troubleshooting

Problem/Display screen	Likely Cause	Action
<b>Absorbance &gt; 3.5!</b>	The measured absorbance exceeds 3.5	Dilute the sample and repeat the measurement
<b>Concentration too high!</b>	Calculated concentration is higher than 999999	Dilute the sample and repeat the measurement
<b>Error Barcode controlnumber? Please update program data!</b>	Deviation to the stored data	Data updating
<b>Error Program not available. Please update program data!</b>	Barcode cuvette test missing	Data updating
<b>Error Clean Cuvette</b>	The cuvette is soiled or there are undissolved particles in the cuvette	Clean the cuvette; allow the particles to settle
<b>Error Please check the UV Lamp.</b>	The lamp output is too low	Check the lamp and replace it if necessary
<b>Error Selfcheck stopped. Please check the lamp. Please close the lid. Error [xx]</b>	Self Check Test stops while starting the instrument	Check the lamp and replace, if necessary. Close the lid. Press <b>Start Again</b> .
<b>Error Selfcheck stopped. Please remove the cuvette Please close the lid.</b>	Self Check Test stops while starting the instrument	Remove the cuvette/sample cell from the cell compartment. Press <b>OK</b> .
<b>Error Sipper?</b>	Kink in tube Peristaltic pump failure	Check the tube connections
<b>Error Selfcheck stopped. Hardware error. Error [x]</b>	Electronic defect	Contact the manufacturer or a sales representative and indicate the error number
<b>Negative result!</b>	The calculated result is negative	Check the concentration of the sample
<b>No evaluation!</b>	Error in the test database / user database	Check the programming Contact the manufacturer or a sales representative
<b>Over measuring range</b>	The measured absorbance is above the calibration range of the test	Dilute the sample and repeat the measurement
<b>Under measuring range</b>	The measured absorbance is below the calibration range of the test	If possible, select a test with a lower measurement range or use a cuvette with a longer path length





## Section 9 Replacement Parts

Description	Cat.-No.
Sipper Module DR 5000	LZV485
Peltier temperature controlled Module DR 5000	LZV513
Pour-Thru Module DR 5000	LZV479
Multi-Cell Holder DR 5000	A23618
Carousel Holder (Sample changer)	A23620
USB-Keyboard (keyboard layout: US)	LZV582
USB-Barcode Scanner (handset scanner for entering sample IDs)	LZV566
Hach Data Trans (PC software for data transfer)	LZY274
Brewery Analysis Software	LZV659
Sipper Tubing, Inlet and Waste	A23800
Sipper Tubing, PHARMED quality	LZV580
Sipper Tubing, comes standard with Sipper Module	LZV780
Pour Thru Tube Kit	LZV569
USB-Memory Stick	2946900
USB-Interface Cable, (1 m)	LZV567
Tungsten (VIS) Lamp	A23778
Deuterium (UV) Lamp	A23792
Fuse	A23772
Filter Pad	A23766
Certified test filter set for self-checks (Verification Kit) (6 precision glass filters with target values)	LZV537
Semi-Micro Cell, 50 mm, Glass	LZP269
Macro Cell, 20 mm, Glass	LZP331
Sample Cell, 1 cm Square Glass, (10 mm path length) matched pair	2095100
Sample Cell, 1 cm Square Quartz, (10 mm path length) matched pair	4822800
Sample Cell, 50 mm Rectangular Glass	2629250
Sample Cell, 50 mm Rectangular Quartz	2624450
Sample Cell, 100 mm Rectangular Glass	2629201
Sample Cell, 100 mm Rectangular Quartz	2624401
Sample Cell, 1" Round Plastic, (1 cm path length) 10 mL, w/cap	4864302
Sample Cell, 1" Square Glass, 10 mL, matched sets	2495402
Sample Cell, 1" Square Glass, 10&25 mL, matched sets	2665902
Sample Cell, 1" Square Plastic	2410212
Pour Thru Cell, 1x1 cm (Quartz/160 µL/ Glass center height 10 mm)	A24209
Pour Thru Cell, 1x1 cm (Quartz/450 µL/ Glass center height 10 mm)	LZV510
Pour Thru Cell, 5x1 cm (Quartz/370 µL/ Glass center height 10 mm)	LZV649
Pour Thru Cell, 1" Square (1" path length)	5913700
Adapter for 100 mm Sample Cell	LZY421
Dust Cover	LZV728



# Section 10 Contact Information

## **HACH Company World Headquarters**

P.O. Box 389  
Loveland, Colorado  
80539-0389 U.S.A.  
Tel (800) 227-HACH  
(800) -227-4224  
(U.S.A. only)  
Fax (970) 669-2932  
orders@hach.com  
www.hach.com

## **HACH LANGE GMBH**

Willstätterstraße 11  
D-40549 Düsseldorf  
Tel. +49 (0)2 11 52 88-320  
Fax +49 (0)2 11 52 88-210  
info@hach-lange.de  
www.hach-lange.de

## **DR. BRUNO LANGE AG**

Juchstrasse 1  
CH-8604 Hegnau  
Tel. +41(0)44 9 45 66 10  
Fax +41(0)44 9 45 66 76  
info@hach-lange.ch  
www.hach-lange.ch

## **HACH LANGE APS**

Åkandevej 21  
DK-2700 Brønshøj  
Tel. +45 36 77 29 11  
Fax +45 36 77 49 11  
info@hach-lange.dk  
www.hach-lange.dk

## **HACH LANGE LDA**

Av. do Forte nº8  
Fracção M  
P-2790-072 Carnaxide  
Tel. +351 214 253 420  
Fax +351 214 253 429  
info@hach-lange.pt  
www.hach-lange.pt

## **HACH LANGE KFT.**

Hegyalja út 7-13.  
H-1016 Budapest  
Tel. +36 (06)1 225 7783  
Fax +36 (06)1 225 7784  
info@hach-lange.hu  
www.hach-lange.hu

## **HACH LANGE D.O.O.**

Fajfarjeva 15  
SI-1230 Domžale  
Tel. +386 (0)59 051 000  
Fax +386 (0)59 051 010  
info@hach-lange.si  
www.hach-lange.si

## **Repair Service in the United States:**

HACH Company  
Ames Service  
100 Dayton Avenue  
Ames, Iowa 50010  
Tel (800) 227-4224  
(U.S.A. only)  
Fax (515) 232-3835

## **HACH LANGE LTD**

Pacific Way  
Salford  
GB-Manchester, M50 1DL  
Tel. +44 (0)161 872 14 87  
Fax +44 (0)161 848 73 24  
info@hach-lange.co.uk  
www.hach-lange.co.uk

## **HACH LANGE FRANCE S.A.S.**

33, Rue du Ballon  
F-93165 Noisy Le Grand  
Tél. +33 (0)1 48 15 68 70  
Fax +33 (0)1 48 15 80 00  
info@hach-lange.fr  
www.hach-lange.fr

## **HACH LANGE AB**

Vinthusdsvägen 159A  
SE-128 62 Sköndal  
Tel. +46 (0)8 7 98 05 00  
Fax +46 (0)8 7 98 05 30  
info@hach-lange.se  
www.hach-lange.se

## **HACH LANGE SP.ZO.O.**

ul. Opolska 143 a  
PL-52-013 Wrocław  
Tel. +48 (0)71 342 10-83  
Fax +48 (0)71 342 10-79  
info@hach-lange.pl  
www.hach-lange.pl

## **HACH LANGE S.R.L.**

Str. Leonida, nr. 13  
Sector 2  
RO-020555 Bucuresti  
Tel. +40 (0) 21 201 92 43  
Fax +40 (0) 21 201 92 43  
info@hach-lange.ro  
www.hach-lange.ro

## **HACH LANGE E.Π.E.**

Αυλιδος 27  
GR-115 27 Αθήνα  
Τηλ. +30 210 7777038  
Fax +30 210 7777976  
info@hach-lange.gr  
www.hach-lange.gr

## **Repair Service in Canada:**

Hach Sales & Service  
Canada Ltd.  
1313 Border Street, Unit 34  
Winnipeg, Manitoba  
R3H 0X4  
Tel (800) 665-7635  
(Canada only)  
Tel (204) 632-5598  
Fax (204) 694-5134  
canada@hach.com

## **HACH LANGE LTD**

Unit 1, Chestnut Road  
Western Industrial Estate  
IRL-Dublin 12  
Tel. +353(0)1 46 02 5 22  
Fax +353(0)1 4 50 93 37  
info@hach-lange.ie  
www.hach-lange.ie

## **HACH LANGE SA**

Motstraat 54  
B-2800 Mechelen  
Tél. +32 (0)15 42 35 00  
Fax +32 (0)15 41 61 20  
info@hach-lange.be  
www.hach-lange.be

## **HACH LANGE S.R.L.**

Via Riccione, 14  
I-20156 Milano  
Tel. +39 02 39 23 14-1  
Fax +39 02 39 23 14-39  
info@hach-lange.it  
www.hach-lange.it

## **HACH LANGE S.R.O.**

Lešanská 2a/1176  
CZ-141 00 Praha 4  
Tel. +420 272 12 45 45  
Fax +420 272 12 45 46  
info@hach-lange.cz  
www.hach-lange.cz

## **HACH LANGE**

8, Kr. Sarafov str.  
BG-1164 Sofia  
Tel. +359 (0)2 963 44 54  
Fax +359 (0)2 866 04 47  
info@hach-lange.bg  
www.hach-lange.bg

## **HACH LANGE E.P.E.**

27, Avlidos str  
GR-115 27 Athens  
Tel. +30 210 7777038  
Fax +30 210 7777976  
info@hach-lange.gr  
www.hach-lange.gr

## **Repair Service in Latin America, the Caribbean, the Far East, Indian Subcontinent, Africa, Europe, or the Middle East:**

Hach Company World  
Headquarters,  
P.O. Box 389  
Loveland, Colorado,  
80539-0389 U.S.A.  
Tel +001 (970) 669-3050  
Fax +001 (970) 669-2932  
intl@hach.com

## **HACH LANGE GMBH**

Hütteldorferstr. 299/Top 6  
A-1140 Wien  
Tel. +43 (0)1 9 12 16 92  
Fax +43 (0)1 9 12 16 92-99  
info@hach-lange.at  
www.hach-lange.at

## **DR. LANGE NEDERLAND B.V.**

Laan van Westroijen 2a  
NL-4003 AZ Tiel  
Tel. +31(0)344 63 11 30  
Fax +31(0)344 63 11 50  
info@hach-lange.nl  
www.hach-lange.nl

## **HACH LANGE S.L.U.**

Edif. Arteaga Centrum  
C/Larrauri, 1C- 2ª Pl.  
E-48160 Derio/Vizcaya  
Tel. +34 94 657 33 88  
Fax +34 94 657 33 97  
info@hach-lange.es  
www.hach-lange.es

## **HACH LANGE S.R.O.**

Roľnícka 21  
SK-831 07 Bratislava –  
Vajnory  
Tel. +421 (0)2 4820 9091  
Fax +421 (0)2 4820 9093  
info@hach-lange.sk  
www.hach-lange.sk

## **HACH LANGE SU ANALİZ SİSTEMLERİ LTD.ŞTİ.**

Hilal Mah. 75. Sokak  
Arman Plaza No: 9/A  
TR-06550 Çankaya/ANKARA  
Tel. +90 (0)312 440 98 98  
Fax +90 (0)312 442 11 01  
bilgi@hach-lange.com.tr  
www.hach-lange.com.tr



## Section 11 Limited Warranty

---

Hach Company warrants its products to the original purchaser against any defects that are due to faulty material or workmanship for a period of one year from date of shipment unless otherwise noted in the product manual.

In the event that a defect is discovered during the warranty period, Hach Company agrees that, at its option, it will repair or replace the defective product or refund the purchase price excluding original shipping and handling charges. Any product repaired or replaced under this warranty will be warranted only for the remainder of the original product warranty period.

This warranty does not apply to consumable products such as chemical reagents; or consumable components of a product, such as, but not limited to, lamps and tubing.

Contact Hach Company or your distributor to initiate warranty support. Products may not be returned without authorization from Hach Company.

### Limitations

This warranty does not cover:

- Damage caused by acts of God, natural disaster, labor unrest, acts of war (declared or undeclared), terrorism, civil strife or acts of any governmental jurisdiction
- Damage caused by misuse, neglect, accident or improper application or installation
- Damage caused by any repair or attempted repair not authorized by Hach Company
- Any product not used in accordance with the instructions furnished by Hach Company
- Freight charges to return merchandise to Hach Company
- Freight charges on expedited or express shipment of warranted parts or product
- Travel fees associated with on-site warranty repair

This warranty contains the sole express warranty made by Hach Company in connection with its products. All implied warranties, including without limitation, the warranties of merchantability and fitness for a particular purpose, are expressly disclaimed.

Some states within the United States do not allow the disclaimer of implied warranties and if this is true in your state the above limitation may not apply to you. This warranty gives you specific rights, and you may also have other rights that vary from state to state.

This warranty constitutes the final, complete, and exclusive statement of warranty terms and no person is authorized to make any other warranties or representations on behalf of Hach Company.

### Limitation of Remedies

The remedies of repair, replacement or refund of purchase price as stated above are the exclusive remedies for the breach of this warranty. On the basis of strict liability or under any other legal theory, in no event shall Hach Company be liable for any incidental or consequential damages of any kind for breach of warranty or negligence.



# Index

---

## A

Absorbance Check ..... 111  
Absorbance Formula ..... 90, 91  
Alphanumeric keypad ..... 21  
Audio signals ..... 25

## B

Barcode Programs ..... 51, 52, 53  
    edit ..... 53  
Basic data ..... 42

## C

Calibration formula ..... 75  
Calibration Settings ..... 66  
Chemical Form ..... 43, 47, 53  
Chemical form ..... 51, 52  
Cleaning  
    Display ..... 121  
    Spectrophotometer ..... 121  
Concentration Factor ..... 87, 88, 90, 92  
Concentration Resolution ..... 87, 88, 90, 92  
Continuous Reading ..... 43, 45  
Cursor Mode ..... 94, 96  
Cuvettes/cells ..... 121

## D

Data  
    delete ..... 35, 37  
    recall ..... 35  
    send ..... 35  
    store ..... 35  
Data Storage ..... 38  
Date ..... 25  
Date and Time ..... 25  
Dilution Factor ..... 43, 45, 53  
Display ..... 21, 121  
Display and Sound ..... 25  
Drift Check ..... 112

## F

Factory Service ..... 116  
Favorites ..... 22, 49, 80, 81  
    delete ..... 80, 81  
    edit ..... 80  
    recall ..... 81  
Filter Settings ..... 36  
Formula ..... 54, 55  
Free programming ..... 72, 73, 75, 78, 80

## H

HACH Data Trans ..... 31, 36  
Hazard Information ..... 9

## I

Instrument Backup ..... 117, 118  
Instrument Information ..... 105  
Instrument Setup ..... 22, 23, 43, 53, 88, 91, 95, 101  
Integral ..... 94, 96  
Interfaces ..... 29  
Internet ..... 58, 106

## L

Language ..... 19

## M

Main Menu ..... 22  
Maintenance ..... 121  
Mark ..... 103  
Measurement Process ..... 54, 55, 72  
Measuring sequence ..... 73  
Multi Wavelength ..... 22, 64, 66, 90, 91, 92, 93

## N

Name ..... 54  
Noise Check ..... 108

## O

Operating environment ..... 13  
Operator ID ..... 23  
    creating ..... 23  
    delete ..... 23  
Optical Checks ..... 106, 107, 108, 110, 111, 112, 113  
Output Checks ..... 115

## P

Password ..... 32  
    activate ..... 32  
PC and Printer ..... 29, 31  
Print data ..... 31  
Printer Setup ..... 29  
Printing data ..... 31  
Program ..... 61  
Programm overview ..... 22

## R

Reading Mode ..... 43, 44, 88, 89  
Reagent Blank ..... 43, 48  
Recall Data ..... 22, 38, 43, 53, 88, 91, 95, 101  
Reference Scan ..... 94, 99

## S

Safety Information ..... 9  
Sample blank ..... 53  
Sample ID ..... 24  
    create ..... 24  
    delete ..... 24  
Save as User Program ..... 43, 53, 88, 90  
Scale & Units ..... 95, 97, 101, 102  
Security List ..... 32  
Select View ..... 96  
Send Data ..... 36, 43, 53, 94, 101  
Service Time ..... 116  
Single Reading ..... 43  
Single Wavelength ..... 22, 43, 45, 63, 66, 87, 88, 89  
Software ..... 58, 106  
Standard Addition ..... 43, 82  
Standard Adjust ..... 43, 46  
Stored Data ..... 35  
Stored Programs ..... 22, 42, 43, 45, 46, 47, 48, 49  
    send Data ..... 43  
Stray Light Check ..... 110  
System Check ..... 116  
System Checks 22, 105, 107, 108, 110, 111, 112, 113,  
    ..... 115, 116, 117

---

**T**

Time & Interval .....	101, 102
Time Course .....	22, 38, 39, 40, 41, 101, 103, 104
Time Course Scan Reading .....	102
Timer .....	25, 43, 53, 54, 56, 87, 90
Touch Screen .....	21

**U**

Unpack the instrument .....	13
Update .....	54, 58, 106
User Programs .....	22, 61

**V**

Variables .....	54, 56, 78
Verification Kit .....	113
Version .....	54, 55
View Table .....	101

**W**

Warranty .....	133
Wavelength .....	87, 90
Wavelength Check .....	107
Wavelength Range .....	95
Wavelength Scan .....	22, 38, 39, 40, 41, 94, 95, 98
Wavelength setting .....	101
Wavelength spectrum .....	94



## Free Manuals Download Website

<http://myh66.com>

<http://usermanuals.us>

<http://www.somanuals.com>

<http://www.4manuals.cc>

<http://www.manual-lib.com>

<http://www.404manual.com>

<http://www.luxmanual.com>

<http://aubethermostatmanual.com>

Golf course search by state

<http://golfingnear.com>

Email search by domain

<http://emailbydomain.com>

Auto manuals search

<http://auto.somanuals.com>

TV manuals search

<http://tv.somanuals.com>